

# BM series

## INSTRUCTION MANUAL

---

### Analytical Balance

BM-5

BM-5D

BM-20

BM-22

BM-252

BM-200

BM-300

BM-500



A&D Company, Ltd.

© 2019 A&D Company, Limited. All rights reserved.

No part of this publication may be reproduced, transmitted, transcribed, or translated into any language in any form by any means without the written permission of A&D Company, Limited.

The contents of this manual and the specifications of the instrument covered by this manual are subject to change for improvement without notice.

Windows, Word and Excel are registered trademarks of Microsoft Corporation.

# Contents

## Basic Operation

|       |  |    |
|-------|--|----|
| 1.    | Introduction.....                                    | 5  |
| 1.1.  | Features.....  | 6  |
| 1.2.  | Compliance.....                                      | 7  |
| 2.    | Unpacking the Balance.....                           | 8  |
| 3.    | Installing the Balance and Precautions.....          | 11 |
| 3.1.  | Installing the Balance.....                          | 11 |
| 3.2.  | Precautions Before Use.....                          | 14 |
| 3.3.  | Precautions During Use.....                          | 16 |
| 3.4.  | Precautions After Use.....                           | 17 |
| 3.5.  | Precautions for Power Supply.....                    | 17 |
| 4.    | Display Symbols and Key Operation.....               | 18 |
| 5.    | Weighing Units.....                                  | 19 |
| 5.1.  | Units.....   | 19 |
| 5.2.  | Storing Active Units.....                            | 23 |
| 5.3.  | Selecting Units and Weighing Modes.....              | 24 |
| 6.    | Weighing.....  | 26 |
| 6.1.  | Basic Operation (Gram Mode).....                     | 26 |
| 6.2.  | Smart Range.....                                     | 28 |
| 6.3.  | Counting Mode (PC).....                              | 29 |
| 6.4.  | Percent Mode (%).....                                | 31 |
| 6.5.  | Ionizer (Neutralization Device).....                 | 32 |
| 6.6.  | Monitoring the Clock and Environmental Sensors.....  | 34 |
| 6.7.  | Repeatability Test.....                              | 35 |
| 7.    | Response Adjustment.....                             | 36 |
| 7.1.  | Automatic Response Adjustment.....                   | 36 |
| 7.2.  | Manual Response Adjustment.....                      | 37 |
| 8.    | Sensitivity Adjustment / Calibration Test.....       | 38 |
| 8.1.  | Automatic Sensitivity Adjustment.....                | 39 |
| 8.2.  | Sensitivity Adjustment Using the Internal Mass.....  | 40 |
| 8.3.  | Calibration Test Using the Internal Mass.....        | 41 |
| 8.4.  | Sensitivity adjustment Using an External Weight..... | 42 |
| 8.5.  | Calibration Test Using an External Weight.....       | 43 |
| 8.6.  | Correcting the Internal Mass Value.....              | 44 |
| 9.    | Function Switch and Initialization.....              | 45 |
| 9.1.  | Permit or Inhibit.....                               | 45 |
| 9.2.  | Initializing the Balance.....                        | 46 |
| 10.   | Function Table.....                                  | 47 |
| 10.1. | Setting the Function Table.....                      | 47 |
| 10.2. | Details of the Function Table.....                   | 49 |
| 10.3. | Description of the Class "Environment, Display"..... | 51 |
| 10.4. | Description of the Item "Data Output Mode".....      | 53 |
| 10.5. | Description of the Item "Data Format".....           | 54 |
| 10.6. | Data Format Examples.....                            | 57 |
| 10.7. | Clock and Calendar Function.....                     | 59 |

|       |  |     |
|-------|--|-----|
| 11.   | ID Number and GLP Report.....  | 60  |
| 11.1. | Setting the ID Number .....  | 60  |
| 11.2. | GLP Report.....  | 61  |
| 12.   | Data Memory.....   | 67  |
| 12.1. | Notes on Using Data Memory .....                                       | 67  |
| 12.2. | Data Memory for Weighing Data .....                                    | 67  |
| 12.3. | Data Memory for Sensitivity adjustment and Calibration Test Data ..... | 71  |
| 13.   | Underhook.....   | 73  |
| 14.   | Density Measurement (Specific gravity) .....                           | 74  |
| 15.   | Password Function.....   | 77  |
| 15.1. | Balance Software Version 1.50 to 1.64 .....                            | 77  |
| 15.2. | Balance Software Version 1.660 or later .....                          | 78  |
| 15.3. | Using the Password Function.....                                       | 79  |
| 15.4. | Entering the Password Before Weighing .....                            | 80  |
| 15.5. | Logging Out .....  | 82  |
| 15.6. | Storing or Changing the Password.....                                  | 83  |
| 15.7. | Changing the Password .....  | 84  |
| 15.8. | Deleting the Password (USER 01~10) .....                               | 85  |
| 15.9. | Forgotten Password.....  | 85  |
| 16.   | Standard Input and Output Interface .....                              | 86  |
| 16.1. | RS-232C Interface .....  | 86  |
| 16.2. | Connecting to Peripheral Equipment.....                                | 87  |
| 16.3. | Commands .....   | 91  |
| 17.   | Key Lock Function.....   | 96  |
| 17.1. | Locking All the Keys.....  | 96  |
| 17.2. | Locking the Specified Keys .....                                       | 97  |
| 18.   | Checking the Balance Software Version .....                            | 97  |
| 19.   | Maintenance.....   | 98  |
| 19.1. | Cleaning and Handling the Balance .....                                | 98  |
| 19.2. | Error Codes.....   | 98  |
| 19.3. | Other Display .....  | 100 |
| 19.4. | Checking the Balance Performance and Environment.....                  | 101 |
| 19.5. | Asking for Repair .....  | 101 |
| 20.   | Specifications .....   | 102 |
| 20.1. | External Dimensions.....   | 104 |
| 20.2. | Options and Peripheral Equipment .....                                 | 105 |
| 21.   | Terms.....   | 110 |

# 1. Introduction

This manual describes how the **BM** series balances work and how to get the most out of them in terms of performance. Read this manual thoroughly before using the balance and keep it at hand for future reference.

## About this manual

This manual consists of the following five parts:

Basic operation ..... Describes precautions, the balance's construction and basic operation.

Adapting to the environment..... Describes response (and stability) adjustment to adapt to the environment where there is vibration or drafts, the way to maintain weighing precision in a variation of ambient temperature, sensitivity adjustment and calibration test.

Selecting functions ..... Describes functions of the balance.

Interface and communication..... Describes the serial interface used for communicating with a computer that requests weighing data and controls the balance, and for use with a printer

Maintenance ..... Describes maintenance, error codes, troubleshooting, specifications and options.

## Balance software version

Some balance operations vary depending on the balance software version. For details on checking the software version, refer to “[18. Checking the Balance Software Version](#)”.

## 1.1. Features

- A built-in ionizer can eliminate static electricity from the weighing sample before weighing, reducing weighing errors.
- Each electrode unit of the ionizer is designed to be removed, cleaned and replaced.
- Automatic sensitivity adjustment, using the internal mass, adapting to changes in temperature.
- Response adjustment adapting to drafts and/or vibration in the environment automatically.
- Memory function to store weighing data, sensitivity adjustment data and calibration test data. When only weighing data is stored, a maximum of 200 data can be stored. Interval mode to store the weighing data periodically.
- Good laboratory practice (GLP) / Good manufacturing practice (GMP) data can be output using the RS-232C serial interface.
- A built-in clock and calendar that can add the time and date to the output data. (Changing the clock setting can be limited to the administrator. Refer to “[15. Password Function](#)”.)
- The password function can be used to restrict balance users or changes to the function settings.
- The key lock function disables the balance key operation and enables the balance to be operated only by commands from external devices.
- Underhook, for measuring density and weighing magnetic materials.
- Multiple weighing units with most of the common units used around the world. Gram, Milligram, Counting mode, Percent mode, Ounce (Avoir), Troy Ounce, Metric carat, Momme, Pennyweight, Grain (UK), Tael (preset at the factory), Tola (India), Messghal and Density mode
- The **BM-5D** and **BM-22** are equipped with a smart range function to weigh in the precision range (readability: 1 µg) after zeroing a tare weight, if within the weighing capacity.
- The micro tube holder is included in the accessories of the **BM-5**, **BM-5D**, **BM-20** and **BM-22**.
- Density mode for calculating the density of a solid.
- The **BM** series balances are equipped with an RS-232C serial interface and a USB interface to communicate with a computer. While outputting data to a printer using the RS-232C interface, data can be transmitted to a computer at the same time using the USB interface. Using the RS-232C interface and the Windows communication tools software (WinCT) enables bi-directional communication between a Windows computer and the balance. The current version of the WinCT can be downloaded from the A&D website.
- Connecting the USB interface of the balance by a cable to a Windows computer allows transmission of the weighing data to Excel or Word.
- When multiple balances have the BM-08 Ethernet interface installed in place of the USB interface and connected to a LAN, data can be acquired from each of them using the WinCT-Plus software.
- When the accessory data logger (**AD-1688**) is connected, weighing data can be stored without using a computer.

## **1.2. Compliance**

### **1.2.1. Compliance with FCC Rules**

---


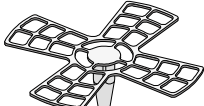




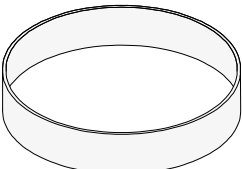
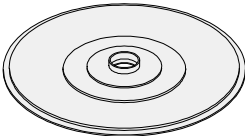



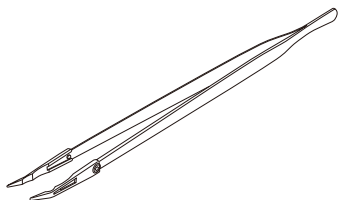
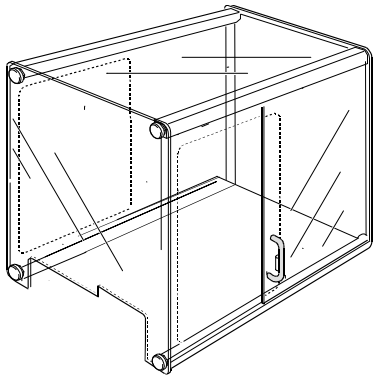

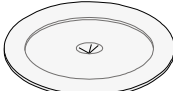
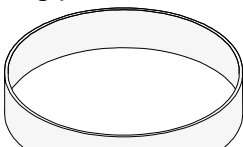
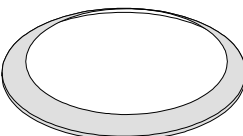
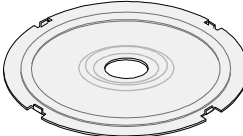
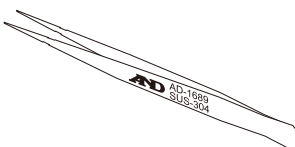
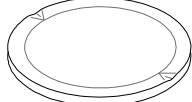
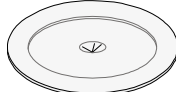
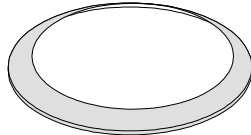
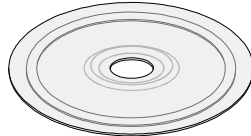

Please note that this equipment generates, uses and can radiate radio frequency energy. This equipment has been tested and has been found to comply with the limits of a Class A computing device pursuant to Subpart J of Part 15 of FCC rules. These rules are designed to provide reasonable protection against interference when equipment is operated in a commercial environment. If this unit is operated in a residential area, it may cause some interference and under these circumstances the user would be required to take, at his own expense, whatever measures are necessary to eliminate the interference.

(FCC = Federal Communications Commission in the U.S.A.)

## 2. Unpacking the Balance


Unpack the balance carefully. Keep the packing material to be used for transporting the balance in the future. See the illustrations to confirm that everything is included.
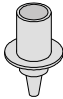
### Accessories for each model

| BM-5, BM-5D, BM-20, BM-22   | BM-252   | BM-200, BM-300<br>BM-500   |
|---|--|--|
|  Small weighing pan for filter<br> Large weighing pan for filter<br> Weighing pan<br> Pan support<br> Micro tube holder<br> Small fine range ring<br> Fine range ring<br> Dust plate<br> $\phi 15$ , 0.8 mL Large<br> $\phi 12$ , 0.3 mL Medium<br> $\phi 8$ , 0.05 mL Small<br>Round aluminum analytical pans (10 pieces each)<br> Tweezers<br>AX-TWEEZERS-25<br> Tabletop Breeze break (M) |  Weighing pan<br> Pan support<br> Fine range ring<br> Breeze break ring<br> Dust plate<br> Tweezers<br>AD-1689<br>SUS-304 |  Weighing pan<br> Pan support<br> Breeze break ring<br> Dust plate<br> Antistatic balance tray<br>AX-TRAY-10ML |





## Common accessories

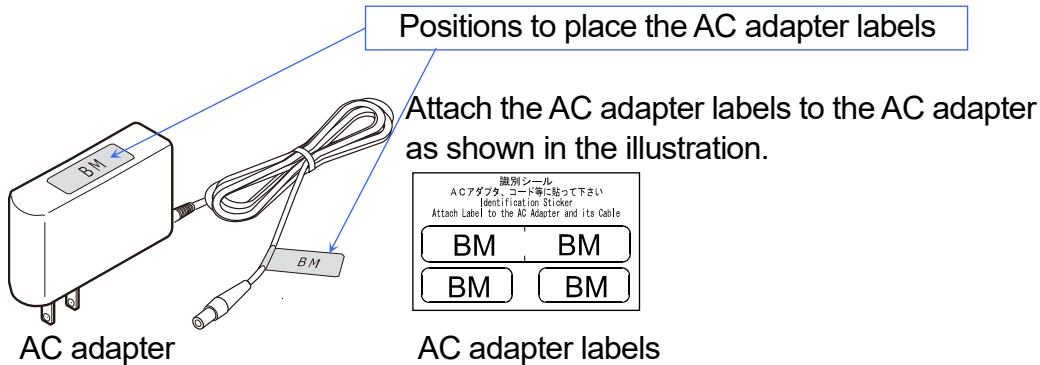


### Notes

- Please confirm that the AC adapter type is correct for your local voltage and receptacle type.
- Please use the dedicated AC adapter specified for the balance.
- Do not use the AC adapter provided with the balance for other models or equipment with which the AC adapter may not be compatible.
- If you use the wrong AC adapter, the balance and other equipment may not operate properly.

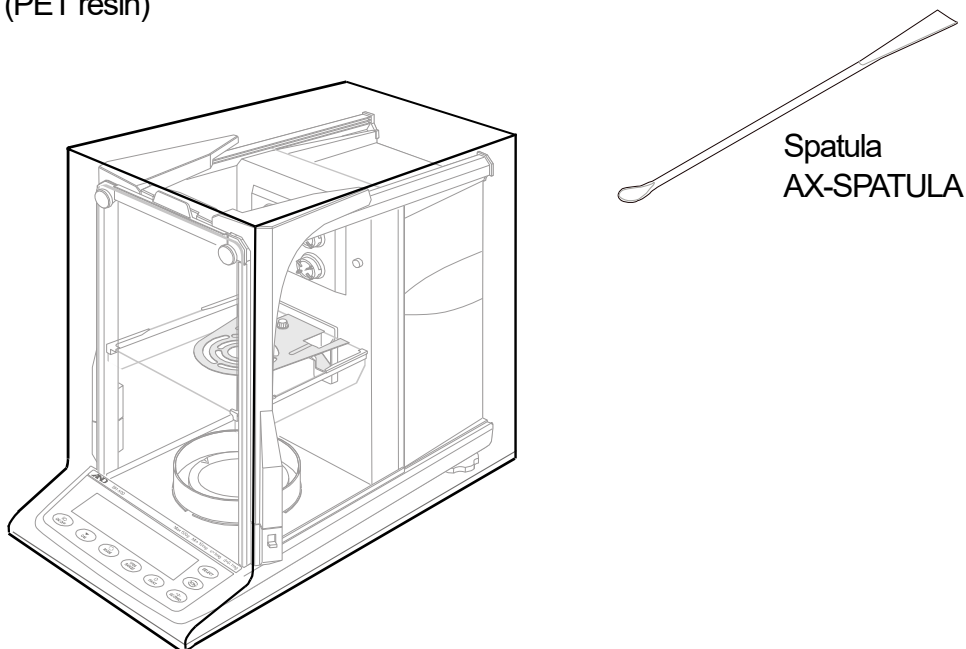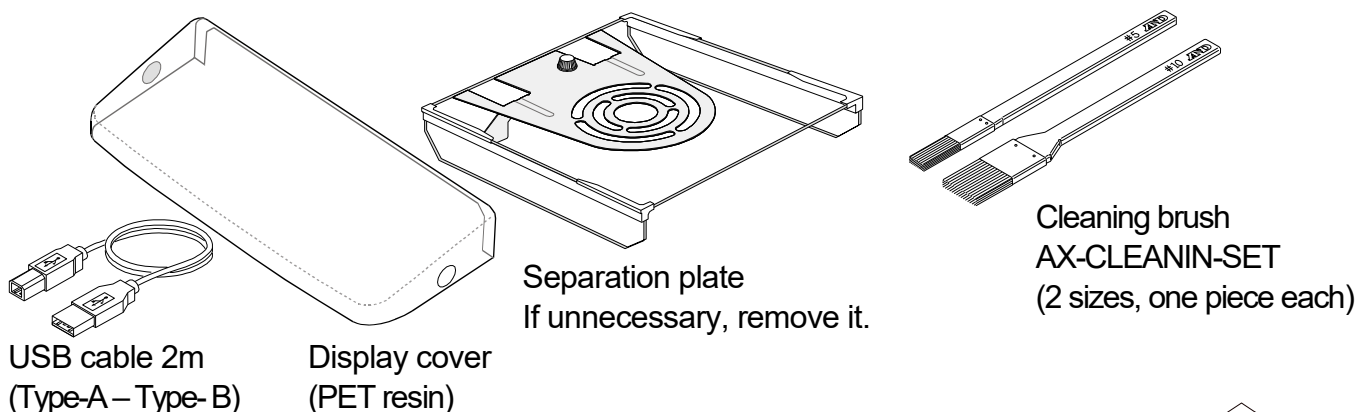

Cover during storage to prevent dust and dirt contamination.  
The cover can be used for the balance with the AC adapter and other cables connected.

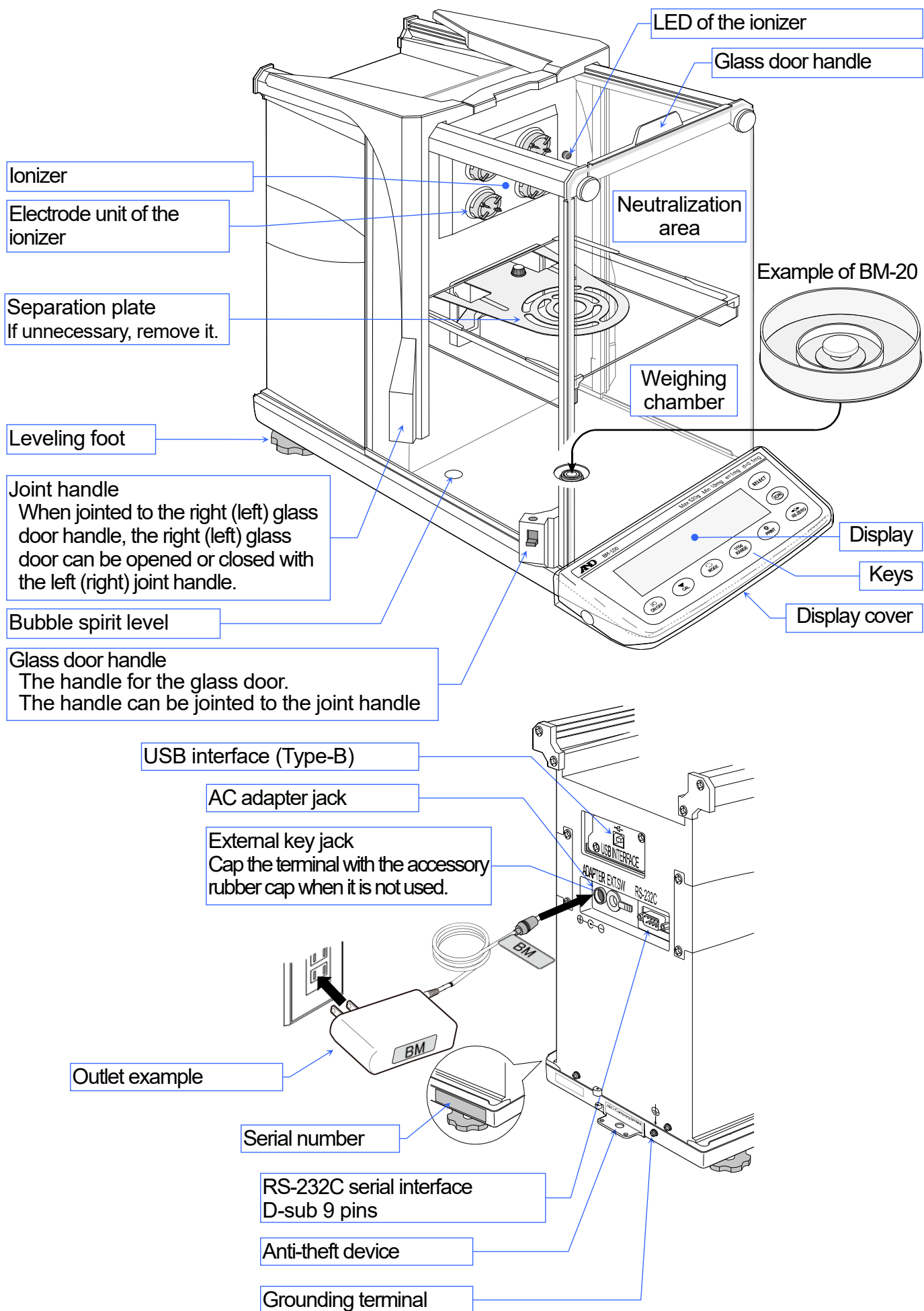


## 3. Installing the Balance and Precautions


### 3.1. Installing the Balance

- 1 Place the balance on a solid weighing table. Refer to "3.2. Precautions Before Use" for details on a place to install the balance.
- 2 Assemble the weighing pan and other parts in accordance with the model and use as shown below.

#### BM-200, BM-300, BM-500

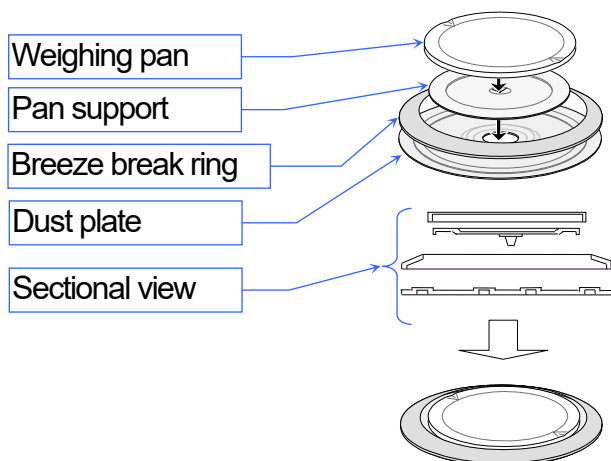


#### BM-252

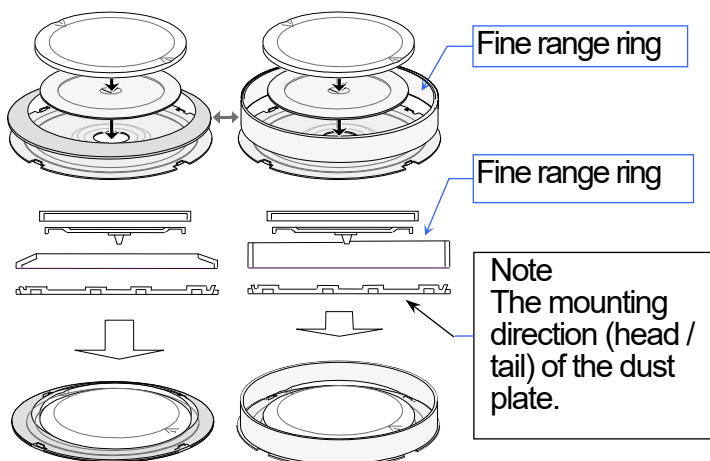

Use the fine range ring in place of the breeze break ring to avoid errors caused by drafts when weighing with a readability of 0.01 mg.

#### BM-5, BM-5D, BM-20, BM-22


**Note** When using the micro tube holder for the BM-5 and BM-5D:  
The micro tube holder is approximately 2 g heavier than the standard pan (weighing pan and pan support). When a micro tube holder is used for the BM-5 and BM-5D, the available weighing range will be up to approximately 3 g including the tare such as a micro tube and test tube.

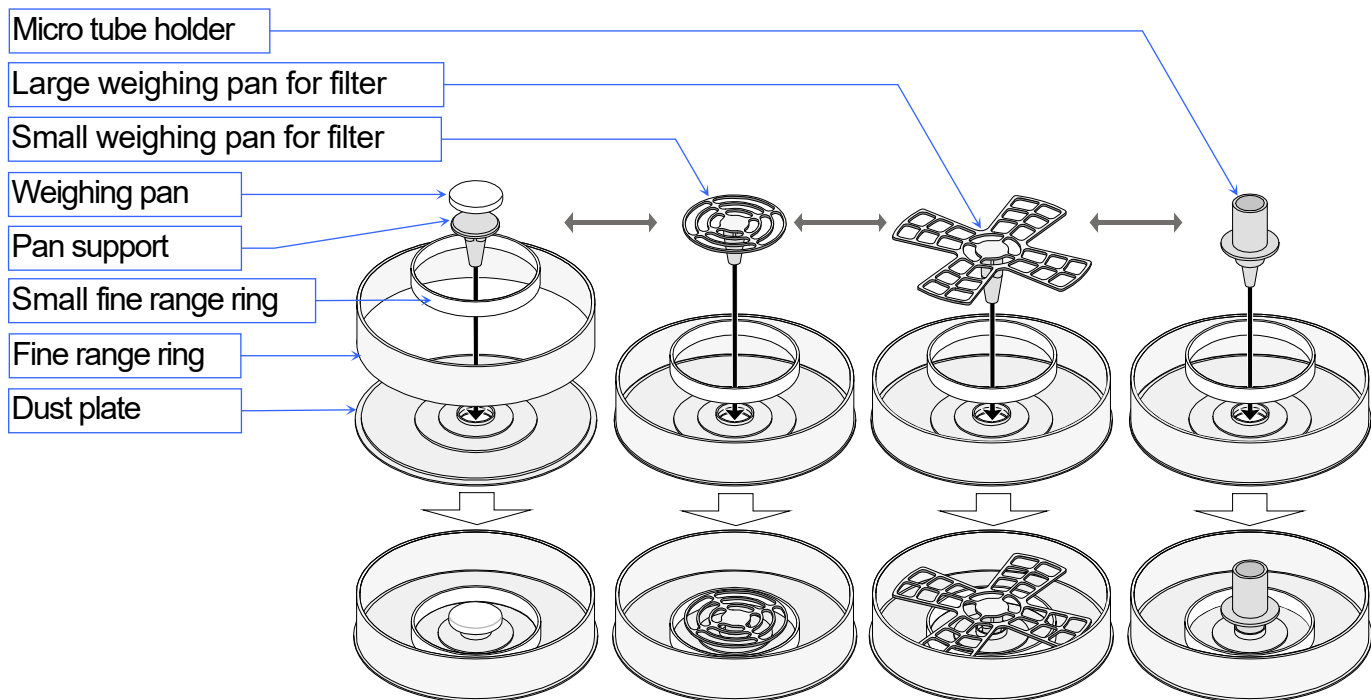


Choose the weighing pan appropriate to the filter size.

Use the holder for micro tube or test tube of  $\phi 12$  mm or less.

## Assembling the parts in the weighing chamber

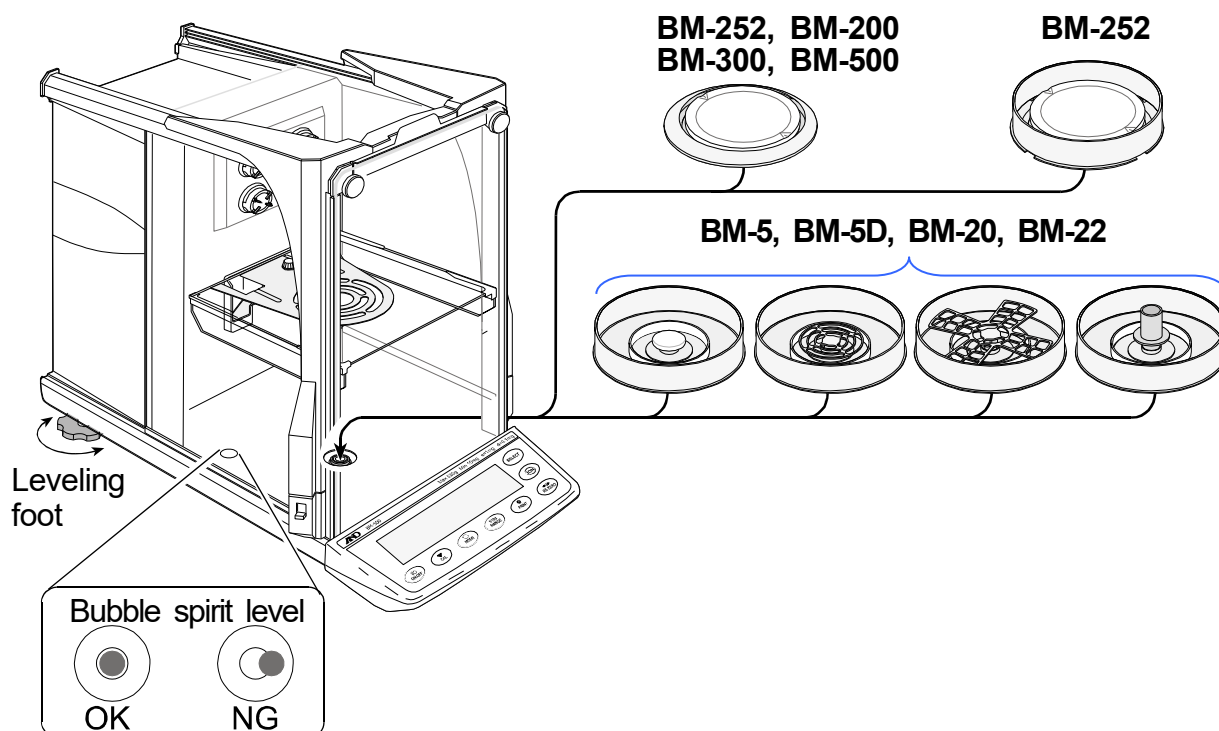

- 3 If the separation plate is not necessary, remove the screw and take the plate off. Refer to "[3.1.2. Removing and attaching the separation plate](#)".
- 4 Adjust the level of the balance using the leveling feet. (Refer to "[Adjusting the level](#)" on page 15.)
- 5 Confirm that the AC adapter type is correct for your local voltage and power receptacle type.
- 6 Connect the specified AC adapter to the balance. Warm up the balance for at least one hour with nothing on the weighing pan.

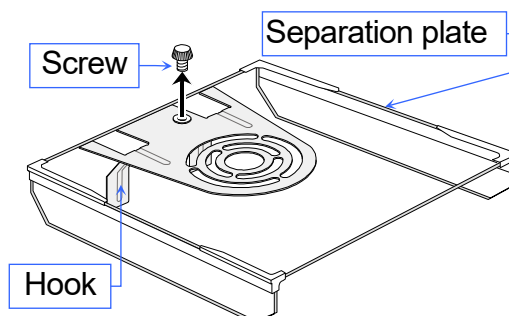


### 3.1.1. How to use the Tabletop breeze break (For BM-5, BM5D, BM-20, BM-22)

---

#### Features

- This Tabletop breeze break is used to prevent from ambient drafts due to air conditioning and body motion so as to reduce weighing error.
- Be sure to install this breeze break on the balance, as it is affected by even the slightest wind.
- This breeze break is designed for the BM series with most suitable use.
- The transparent panel assembly consists of antistatic plastic material, which protects the balance from static electricity.
- The rear panel is equipped the hole that is used to pass the RS-232C and AC adapter cables through.

#### How to install

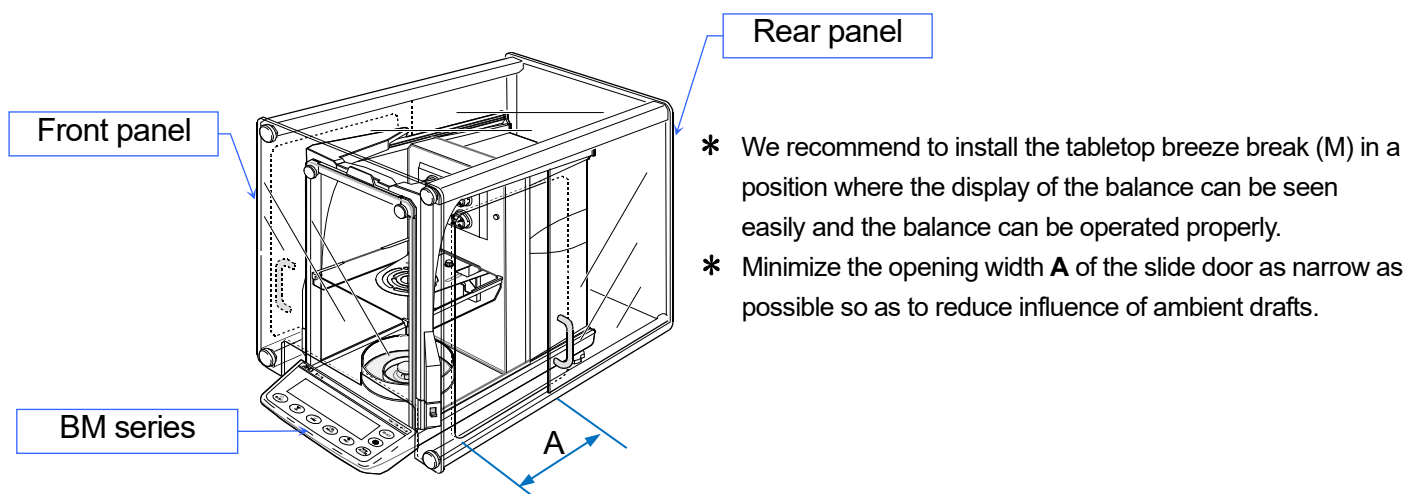

#### Caution

- Do not use organic solvents to clean the breeze break because it may damage the transparency and anti-static properties of the panels.
- Use a soft lint free cloth dampened with warm water and a mild detergent when cleaning the breeze break.

### 3.1.2. Removing and attaching the separation plate

**Note** Take care not to damage the separation plate when handling it.


#### Removing the separation plate

- 1 Support the separation plate and remove the screw from the plate.
- 2 Lift the plate and unhook the hooks.
- 3 Rotate the plate on the axis of the front edge.  
Rotate the plate on the axis of the back edge.
- 4 Remove the plate from the chamber.

#### Attaching the separation plate

Use the arrows in reverse to attach the plate.

- 5 Insert the plate into the chamber.
- 6 Rotate the plate on the axis of the back edge.  
Rotate the plate on the axis of the front edge.  
Level the plate.
- 7 Insert the hooks and hook the plate.
- 8 Support the plate and secure the screw.

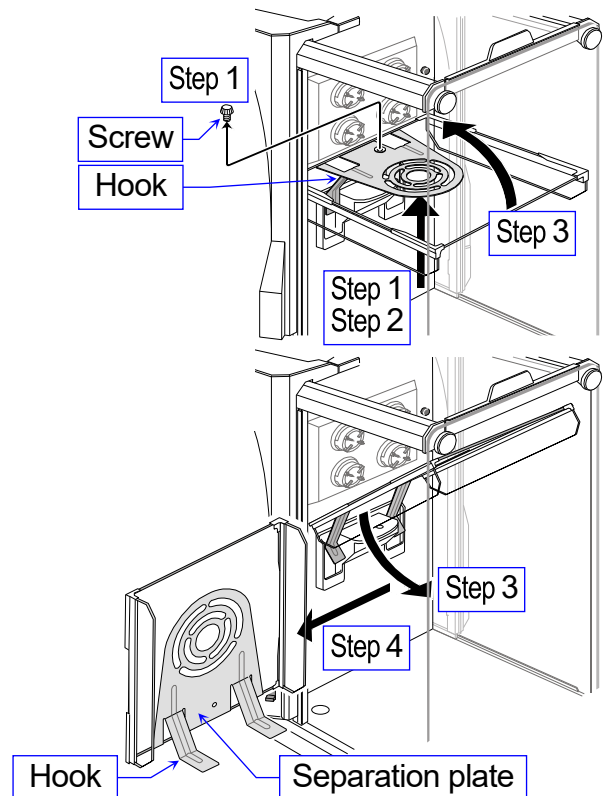



## 3.2. Precautions Before Use

To ensure that you get the most from your balance, please try to follow these conditions as closely as possible, especially for the **BM-5**, **BM-5D**, **BM-20** and **BM-22** that are sensitive.

- The best operating temperature is about 20°C / 68°F at about 50% Relative Humidity.
- The weighing room should be free of dust.
- Clean the weighing chamber and its surroundings to prevent foreign matter from invading, especially when weighing a very small amount of sample.
- The weighing table should be solid and free from vibration, drafts (such as frequently opening doors or windows) and as level as possible. We recommend that you use the anti-vibration table (**AD-1671**) and the remote controller (**AD-8922A**) for the **BM-5**, **BM-5D**, **BM-20** and **BM-22**.
- Do not install the balance where it will be subject to vibration. Corners of rooms on the first floor are best.
- Do not install the balance near a heater, air conditioner, or in a breeze. The influence of breezes can be reduced by using the tabletop breeze break (L) (**AD-1672/AD-1672A**) or the tabletop breeze break (M) (**AD-1676**).
- Do not install the balance in direct sunlight. Avoid excessive temperature changes.
- Do not use the balance near other equipment which produces magnetic fields.
- Adjust the level of the balance using the leveling feet.
- Please warm up the balance for at least one hour. Plug in the AC adapter as usual.
- Adjust the sensitivity of the balance if it is installed for the first time or relocated, or when a weighing pan is replaced.
- Ensure a stable power source when using the AC adapter.

**⚠ Do not place or use the balance where there is flammable or corrosive gas present.**

## Adjusting the level

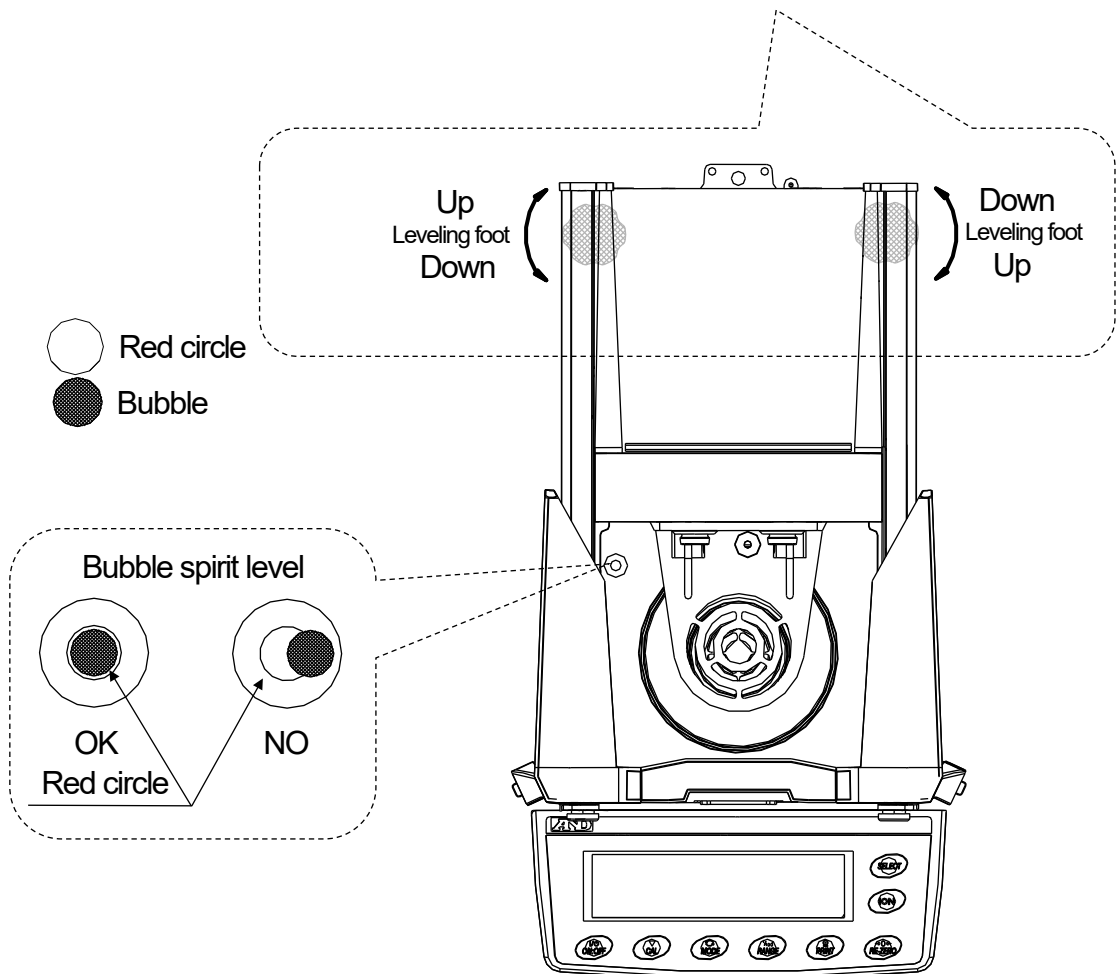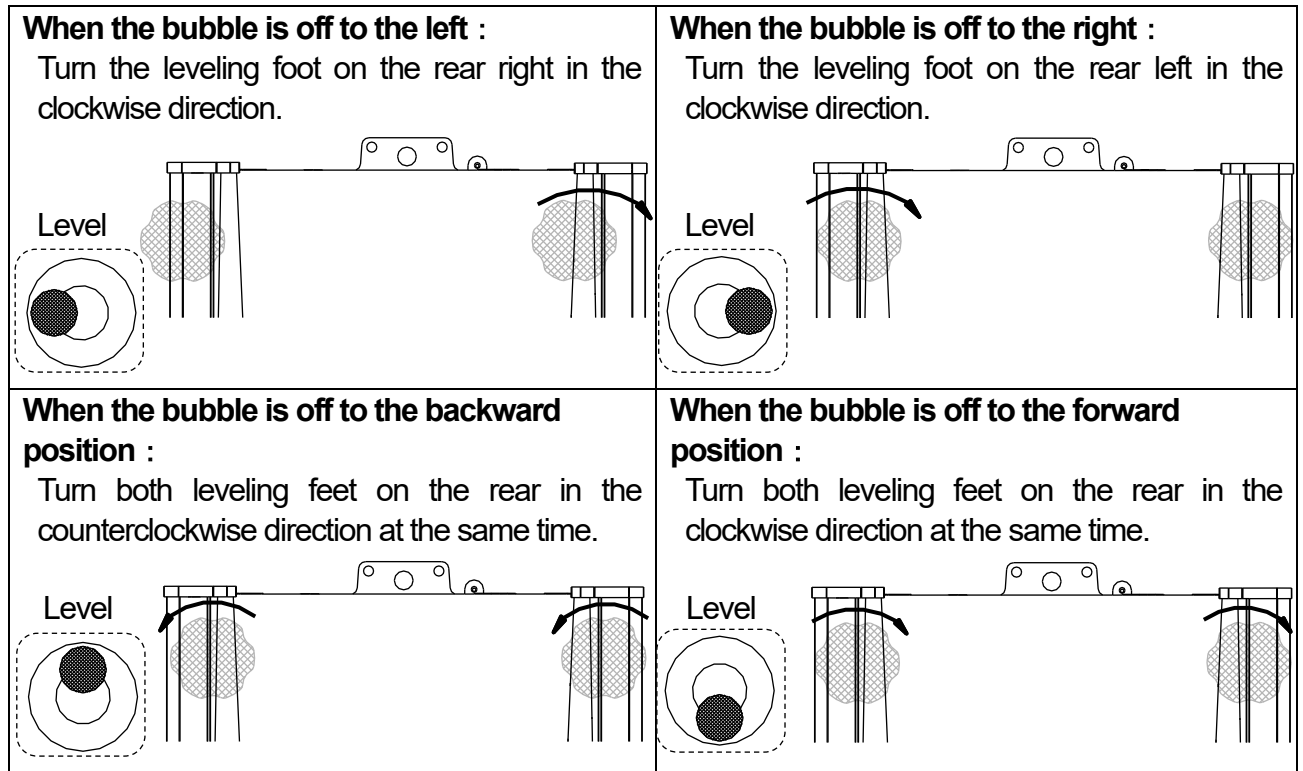




### 3.3. Precautions During Use

Note the following items to get accurate weighing data.

- Discharge static electricity from the weighing sample. When a weighing sample (plastics, insulator, etc.) could have a static charge, the weight value is influenced. Ground the balance, and
  - Eliminate the static electricity using the built-in ionizer.
  - Try to keep the ambient humidity at or above 45% RH in the room.
  - Use a metal shield case.
  - Wipe a charged material (plastic sample, etc.) with a damp cloth.
  - Set up the balance using a ground terminal.
- This balance uses a strong magnet as part of the balance assembly, so please use caution when weighing magnetic materials. If there is a problem, use the underhook (refer to "[13. Underhook](#)") to suspend the material away from the influence of the magnet.
- Eliminate temperature differences between the weighing sample and the environment. When a sample is warmer (cooler) than the ambient temperature, the sample will appear lighter (heavier) than the true mass. This error is due to the rising (falling) draft around the sample. If you touch the sample, the same type of error will occur. Do not touch the sample directly with your hand. Use tweezers or other tools.
- Do not drop things upon the weighing pan, or place a weight beyond the range of the balance on the weighing pan.
- Make each weighing gently and quickly to avoid errors due to changes in the environmental conditions.
- Use the **fine range ring** and attach the **separation plate** (in condition at shipping from factory) in order to avoid weighing errors caused by drafts when weighing using the **BM-252** (readability: 0.01 mg).
- Use the **small fine range ring and fine range ring** and attach the **separation plate** (in condition at shipping from factory) in order to avoid weighing errors caused by drafts when weighing using the **BM-5**, **BM-5D**, **BM-20** or **BM-22** (readability: 0.01 mg and 0.001 mg). To reduce the influence of breezes, minimize operation of doors of the weighing chamber. In addition, in order to weigh more accurately while reducing the effects of the surrounding environment, be sure to install the tabletop breeze break (M) to the weighing chamber.
- The weighing value of the **BM-5**, **BM-5D**, **BM-20** and **BM-22** may change after the stabilization indicator is displayed. We recommend that you set a certain amount of wait time (approx. 10 seconds) before reading the value after the stabilization indicator is displayed.
- If contamination may be a problem, perform sampling operations outside the weighing chamber in order to prevent the sample from scattering in the weighing chamber.
- Take into consideration the affect of air buoyancy on a sample when more accuracy is required.

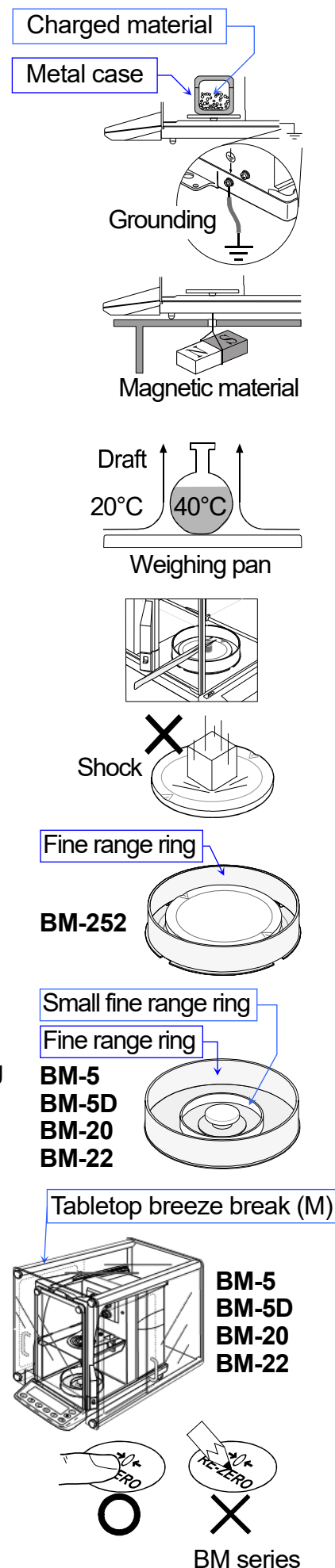



- Do not use a sharp instrument (such as a pencil or ball point pen) to press the keys, use your finger only.
- Press the **RE-ZERO** key before each weighing to prevent possible errors.
- Keep the balance interior free of foreign matter (dust, liquid or metal fragments).

### 3.4. Precautions After Use


- Avoid mechanical shock to your balance.
- Do not disassemble the balance. Contact your local A&D dealer if your balance needs service or repair.
- Do not use solvents to clean the balance. For best cleaning, wipe with a dry lint free cloth or a lint free cloth that is moistened with warm water and a mild detergent.
- Keep the balance interior free of foreign matter (dust, liquid or metal fragments).

### 3.5. Precautions for Power Supply

- Do not remove the AC adapter while the internal mass is in motion, for example, right after the AC adapter is connected, or during sensitivity adjustment using the internal mass.  
If the AC adapter is removed under the conditions described above, the internal mass will be left unsecured, that may cause mechanical damage when the balance is moved.  
Before removing the AC adapter, press the **ON:OFF** key and confirm that zero is displayed in the weighing mode.
- After connecting the AC adapter to the balance, warm up the balance for at least one hour before weighing. ( Four hours or more with the BM-5, BM-5D, and BM-20, BM-22)
- When the AC adapter is connected, the balance is in the standby mode if the standby indicator is on. This is a normal state and does not harm the balance. For accurate weighing, we recommend that the AC adapter be kept connected, especially for the **BM-5, BM-5D, BM-20** and **BM-22**.

## 4. Display Symbols and Key Operation


### Display

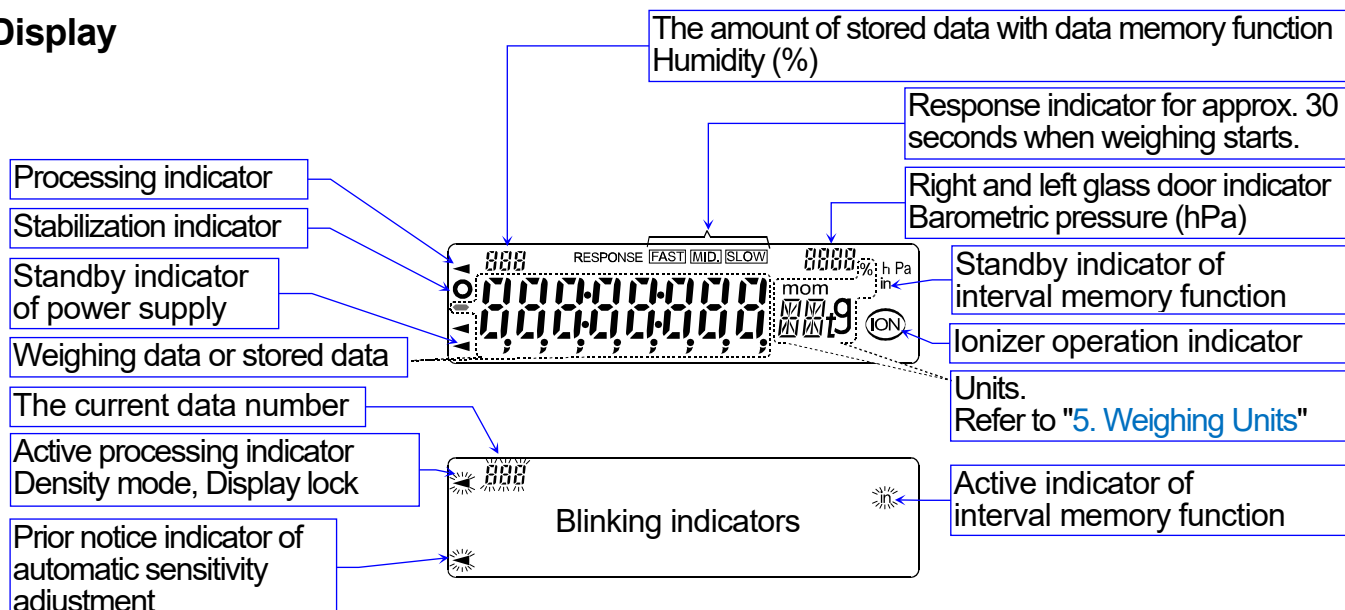

The interval memory function is used to store the weighing data periodically. Refer to "12. Data Memory".

### Key operation

- Press and release the key immediately or "Click the key"
- Press and hold the key (for approx. 2 seconds)

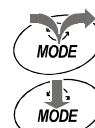


| Key | When pressed and released   | When pressed and held (approx. 2 seconds)  |
|-----|---|--|
|     | Turns the display on and off. The standby indicator is displayed when the display is turned off. When the display is turned on, the weighing mode is enabled. If the password function is used, <b>PASS</b> is displayed to prompt you to enter a password. Refer to "15.4. Entering the Password Before Weighing". This key is available anytime. Pressing the key during operation will interrupt the operation and turn the display off. |  |
|     | Cancels the operation when performing function settings. Enters the sensitivity adjustment mode using the internal mass.  | Displays other items of the sensitivity adjustment menu.   |
|     | Switches the preset weighing units stored in the function table. Refer to "5. Weighing Units".  | Performs automatic response adjustment.  |
|     | <ul style="list-style-type: none"> <li>□ In the weighing mode, turns the minimum weighing value on and off.</li> <li>□ In the counting or percent mode, enters the unit mass storing mode or reference mass storing mode.</li> </ul>  | <ul style="list-style-type: none"> <li>□ Displays the function table menu. Refer to "10. Function Table".</li> <li>□ Performs a repeatability test when pressed and held for another 2 seconds after the function table menu is displayed. Refer to "6.7 Repeatability Test".</li> </ul>             |
|     | The key to output the weighing data to a printer or personal computer (or store it in memory) depending on the function table settings. (Factory setting = output)  | <ul style="list-style-type: none"> <li>□ No function at the factory setting.</li> <li>□ By changing the function table: <ul style="list-style-type: none"> <li>■ "Title block" and "End block" for GLP / GMP report are output.</li> <li>■ The data memory menu is displayed.</li> </ul> </li> </ul> |
|     | The key to set the display to zero.   |  |
|     | The key to turn the ionizer on and off.   |  |
|     | The key to monitor the time and date, and the environmental sensors (temperature, humidity, barometric pressure), for several seconds. Refer to "6.6 Monitoring the Clock and Environmental Sensors".   |  |

## 5. Weighing Units

### 5.1. Units

- The units and weighing modes can be selected and stored in the function table as described in "5.2. Storing Active Units".  
The sequence of displaying them can be arranged to fit the frequency of use.  
They are stored and are maintained in non-volatile memory, even if the AC adapter is removed.
- If the law in your area permits, you may use all of the units. You can disable the units that you don't regularly use. And you are able to turn them back on.
- If a weighing mode (or unit of mass) has been turned off, that mode or unit will be missing in the sequence. Tael has four varieties, one of which can be selected and installed at the factory.
- Press the **[MODE]** key to select a unit or mode for weighing.
- For details about the units and modes, see the table below:

| Name (unit, mode)            | Abbreviation | Display unit   | Conversion factor |
|------------------------------|--------------|--|-------------------|
| Gram                         | g            | g  | 1 g               |
| Milligram                    | mg           | mg   | 0.001 g           |
| Counting mode                | PC           | PC   | -                 |
| Percent mode                 | %            | %  | -                 |
| Ounce (Avoir)                | oz           | oz   | 28.349523125 g    |
| Troy Ounce                   | ozt          | ozt  | 31.1034768 g      |
| Metric Carat                 | ct           | ct   | 0.2 g             |
| Momme                        | mom          | mom  | 3.75 g            |
| Pennyweight                  | dwt          | dwt  | 1.55517384 g      |
| Grain (UK)                   | GN           | GN   | 0.06479891 g      |
| Tael (HK general, Singapore) | tl           | tL   | 37.7994 g         |
| Tael (HK jewelry)            |              |  | 37.429 g          |
| Tael (Taiwan)                |              |  | 37.5 g            |
| Tael (China)                 |              |  | 31.25 g           |
| Tola (India)                 | t            | t  | 11.6638038 g      |
| Messghal                     | m            | m  | 4.6875 g          |
| Density mode                 | DS           | Refer to "14. Density Measurement (Specific gravity) " |                   |

- Density mode
  - To use the density mode, it must be stored in the function table as described on page 23.  
For details about this mode, refer to "14. Density Measurement (Specific gravity) ".
  - To select the density mode, press the **[MODE]** key until the processing indicator  blinks with the unit "g" displayed.

## Capacity and readability for the BM-252, BM-200, BM-300, BM-500

- The tables below indicate the weighing capacity and the readability for each balance.

| Unit                         | BM-200   | BM-300 | BM-500 |             |
|------------------------------|----------|--------|--------|-------------|
|                              | Capacity |        |        | Readability |
| Gram                         | 220      | 320    | 520    | 0.0001      |
| Milligram                    | 220000   | 320000 | 520000 | 0.1         |
| Ounce (Avoir)                | 7.76     | 11.2   | 18.3   | 0.00001     |
| Troy Ounce                   | 7.07     | 10.2   | 16.7   | 0.00001     |
| Metric Carat                 | 1100     | 1600   | 2600   | 0.001       |
| Momme                        | 58.6     | 85.3   | 138    | 0.0001      |
| Pennyweight                  | 141      | 205    | 334    | 0.0001      |
| Grain (UK)                   | 3395     | 4938   | 8024   | 0.002       |
| Tael (HK general, Singapore) | 5.82     | 8.46   | 13.7   | 0.00001     |
| Tael (HK jewelry)            | 5.87     | 8.54   | 13.8   | 0.00001     |
| Tael (Taiwan)                | 5.86     | 8.53   | 13.8   | 0.00001     |
| Tael (China)                 | 7.04     | 10.2   | 16.6   | 0.00001     |
| Tola (India)                 | 18.8     | 27.4   | 44.5   | 0.00001     |
| Messghal                     | 46.9     | 68.2   | 110    | 0.0001      |

| Unit                         | BM-252   |             |
|------------------------------|----------|-------------|
|                              | Capacity | Readability |
| Gram                         | 250      | 0.00001     |
| Milligram                    | 250000   | 0.01        |
| Ounce (Avoir)                | 8.81     | 0.000001    |
| Troy Ounce                   | 8.03     | 0.000001    |
| Metric Carat                 | 1250     | 0.0001      |
| Momme                        | 66.6     | 0.00001     |
| Pennyweight                  | 160      | 0.00001     |
| Grain (UK)                   | 3858     | 0.0002      |
| Tael (HK general, Singapore) | 6.61     | 0.000001    |
| Tael (HK jewelry)            | 6.67     | 0.000001    |
| Tael (Taiwan)                | 6.66     | 0.000001    |
| Tael (China)                 | 8.00     | 0.000001    |
| Tola (India)                 | 21.4     | 0.000001    |
| Messghal                     | 53.3     | 0.00001     |

## Capacity and readability for the BM-5 and BM-5D

- The **BM-5D** is equipped with the precision range and standard range of the smart range function.

| Unit                         | <b>BM-5</b>            |             |
|------------------------------|------------------------|-------------|
|                              | <b>Precision range</b> |             |
|                              | Capacity               | Readability |
| Gram                         | 5.20                   | 0.000001    |
| Milligram                    | 5200                   | 0.001       |
| Ounce (Avoir)                | 0.183                  | 0.0000001   |
| Troy Ounce                   | 0.167                  | 0.0000001   |
| Metric Carat                 | 26.0                   | 0.00001     |
| Momme                        | 1.38                   | 0.000001    |
| Pennyweight                  | 3.33                   | 0.000001    |
| Grain (UK)                   | 80.1                   | 0.00002     |
| Tael (HK general, Singapore) | 0.137                  | 0.0000001   |
| Tael (HK jewelry)            | 0.138                  | 0.0000001   |
| Tael (Taiwan)                | 0.138                  | 0.0000001   |
| Tael (China)                 | 0.166                  | 0.0000001   |
| Tola (India)                 | 0.444                  | 0.0000001   |
| Messghal                     | 1.10                   | 0.000001    |

| Unit                         | <b>BM-5D</b>           |             |                       |             |
|------------------------------|------------------------|-------------|-----------------------|-------------|
|                              | <b>Precision range</b> |             | <b>Standard range</b> |             |
|                              | Capacity               | Readability | Capacity              | Readability |
| Gram                         | 2.10                   | 0.000001    | 5.20                  | 0.00001     |
| Milligram                    | 2100                   | 0.001       | 5200                  | 0.01        |
| Ounce (Avoir)                | 0.0737                 | 0.0000001   | 0.183                 | 0.000001    |
| Troy Ounce                   | 0.0671                 | 0.0000001   | 0.167                 | 0.000001    |
| Metric Carat                 | 10.5                   | 0.00001     | 26.0                  | 0.0001      |
| Momme                        | 0.560                  | 0.000001    | 1.38                  | 0.00001     |
| Pennyweight                  | 1.34                   | 0.000001    | 3.33                  | 0.00001     |
| Grain (UK)                   | 32.4                   | 0.00002     | 80.1                  | 0.0001      |
| Tael (HK general, Singapore) | 0.0551                 | 0.0000001   | 0.137                 | 0.000001    |
| Tael (HK jewelry)            | 0.0560                 | 0.0000001   | 0.138                 | 0.000001    |
| Tael (Taiwan)                | 0.0560                 | 0.0000001   | 0.138                 | 0.000001    |
| Tael (China)                 | 0.0671                 | 0.0000001   | 0.166                 | 0.000001    |
| Tola (India)                 | 0.179                  | 0.0000001   | 0.444                 | 0.000001    |
| Messghal                     | 0.444                  | 0.000001    | 1.10                  | 0.00001     |

## Capacity and readability for the BM-20 and BM-22

- The **BM-22** is equipped with the precision range and standard range of the smart range function.

| Unit                         | <b>BM-20</b>           |             |
|------------------------------|------------------------|-------------|
|                              | <b>Precision range</b> |             |
|                              | Capacity               | Readability |
| Gram                         | 22.0                   | 0.000001    |
| Milligram                    | 22000                  | 0.001       |
| Ounce (Avoir)                | 0.776                  | 0.0000001   |
| Troy Ounce                   | 0.707                  | 0.0000001   |
| Metric Carat                 | 110                    | 0.00001     |
| Momme                        | 5.86                   | 0.000001    |
| Pennyweight                  | 14.1                   | 0.000001    |
| Grain (UK)                   | 339                    | 0.00002     |
| Tael (HK general, Singapore) | 0.582                  | 0.0000001   |
| Tael (HK jewelry)            | 0.587                  | 0.0000001   |
| Tael (Taiwan)                | 0.586                  | 0.0000001   |
| Tael (China)                 | 0.704                  | 0.0000001   |
| Tola (India)                 | 1.88                   | 0.0000001   |
| Messghal                     | 4.69                   | 0.000001    |

| Unit                         | <b>BM-22</b>           |             |                       |             |
|------------------------------|------------------------|-------------|-----------------------|-------------|
|                              | <b>Precision range</b> |             | <b>Standard range</b> |             |
|                              | Capacity               | Readability | Capacity              | Readability |
| Gram                         | 5.10                   | 0.000001    | 22.0                  | 0.00001     |
| Milligram                    | 5100                   | 0.001       | 22000                 | 0.01        |
| Ounce (Avoir)                | 0.179                  | 0.0000001   | 0.776                 | 0.000001    |
| Troy Ounce                   | 0.163                  | 0.0000001   | 0.707                 | 0.000001    |
| Metric Carat                 | 25.5                   | 0.00001     | 110                   | 0.0001      |
| Momme                        | 1.36                   | 0.000001    | 5.86                  | 0.00001     |
| Pennyweight                  | 3.27                   | 0.000001    | 14.1                  | 0.00001     |
| Grain (UK)                   | 78.7                   | 0.00002     | 339                   | 0.0001      |
| Tael (HK general, Singapore) | 0.134                  | 0.0000001   | 0.582                 | 0.000001    |
| Tael (HK jewelry)            | 0.136                  | 0.0000001   | 0.587                 | 0.000001    |
| Tael (Taiwan)                | 0.136                  | 0.0000001   | 0.586                 | 0.000001    |
| Tael (China)                 | 0.163                  | 0.0000001   | 0.704                 | 0.000001    |
| Tola (India)                 | 0.437                  | 0.0000001   | 1.88                  | 0.000001    |
| Messghal                     | 1.08                   | 0.000001    | 4.69                  | 0.00001     |

## 5.2. Storing Active Units

- The units and modes can be selected and stored in the function table.  
The sequence of displaying them can be arranged to fit the frequency of use.  
The units stored are maintained in non-volatile memory, even if the AC adapter is removed.


- 1 In the weighing mode, press and hold the **RANGE** key (for approx. 2 seconds) until **bASFnC** is displayed, and then release the key.
- 2 Press the **RANGE** key several times to display **Unit**.
- 3 Press the **PRINT** key to enter the unit selection mode.
- 4 Specify a unit or mode in the order to be displayed using the following keys.  
**RANGE** key .....To display the units sequentially.  
**RE-ZERO** key .....To specify a unit or mode.

The indicator **○** appears when the displayed unit or mode is specified.


Examples

| Unit          |    | Display   |
|---------------|----|-----------|
| Gram          | g  | ° Unit g  |
| Milligram     | mg | ° Unit mg |
| Counting mode | PC | ° Unit PC |
| Percent mode  | %  | ° Unit %  |
| Density mode  | DS | ° Unit DS |


- 5 Press the **PRINT** key to store the units or modes.  
The balance displays **End** and then displays the next menu item of the function table.
- 6 Press the **CAL** key to exit the function table.  
Then the balance returns to the weighing mode with the selected unit.

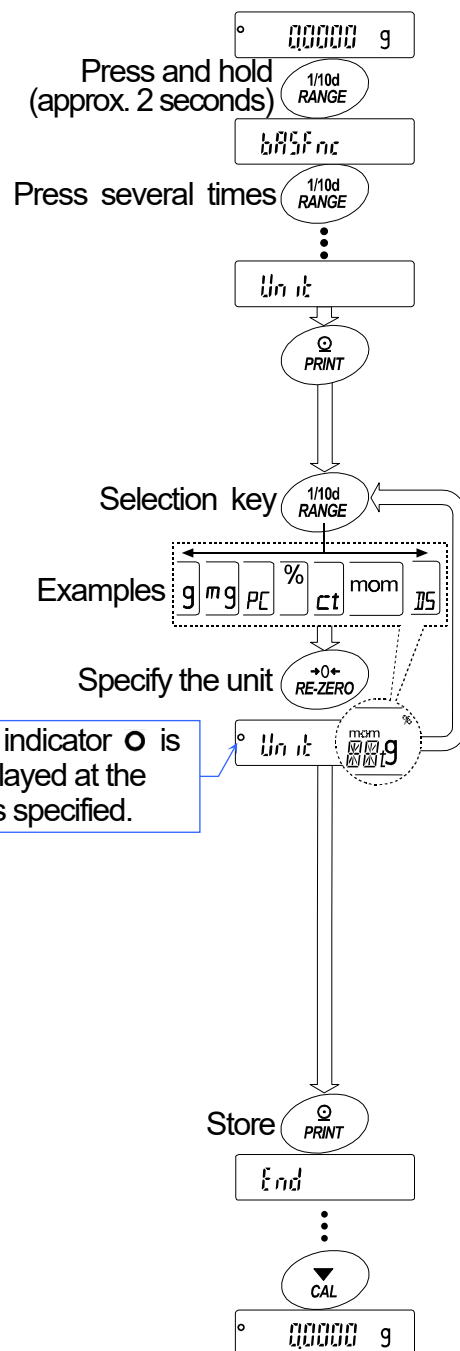

### 5.3. Selecting Units and Weighing Modes

- Preset the sequence of displaying units and weighing modes to fit the frequency of use. Refer to "5.2 Storing Active Units".
- Press the **MODE** key in the weighing mode to display the units and weighing modes in order.
- For the **BM-5**, **BM-5D**, **BM-20** and **BM-22**, the units preset at the factory, **mg** (Milligram) and **g** (Gram), are displayed in this order.
- For the **BM-252**, **BM-200**, **BM-300** and **BM-500**, the units preset at the factory, **g** (Gram) and **mg** (Milligram), are displayed in this order.
- Density mode
  - To use the density mode, it must be stored in the function table as described on page 23. Refer to "14. Density Measurement (Specific gravity)" for details about the mode.
  - To select the density mode, press the **MODE** key until the processing indicator  blinks with the unit "g" displayed.

The unit sequence of the factory settings.  
Turning the display on.

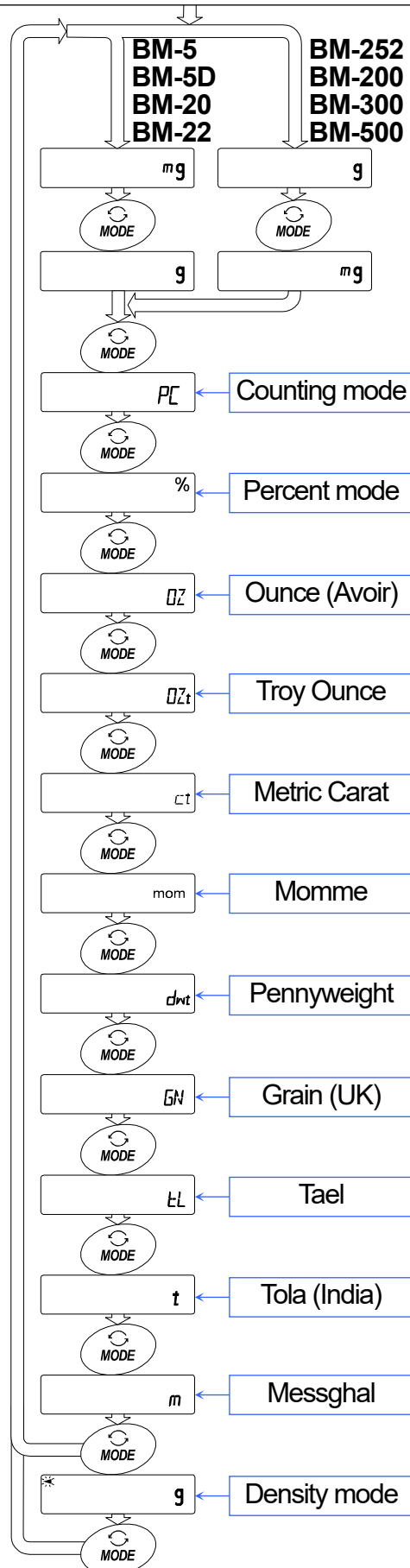




### 5.3.1. Unit Setting Example

The example below sets the units in the order with g (gram) as the first unit followed by PC (counting mode).

- 1 In the weighing mode, press and hold the **[RANGE]** key (for approx. 2 seconds) until **bASFc** is displayed, and then release the key.
- 2 Press the **[RANGE]** key several times to display **Unit**.
- 3 Press the **[PRINT]** key to enter the unit selection mode.
- 4 Press the **[RE-ZERO]** key to specify the unit of g.  
The stabilization indicator **○** appears when the unit is specified.
- 5 Press the **[RANGE]** key several times to display **Unit PC**.
- 6 Press the **[RE-ZERO]** key to specify the unit of PC.  
The stabilization indicator **○** appears when the unit is specified.
- 7 Press the **[PRINT]** key to store the units.  
The balance displays **End** and then displays the next menu item of the function table.
- 8 Press the **[CAL]** key to exit the function table. Then the balance returns to the weighing mode with g, the unit selected first.
- 9 Press the **[MODE]** key to switch between g and PC.

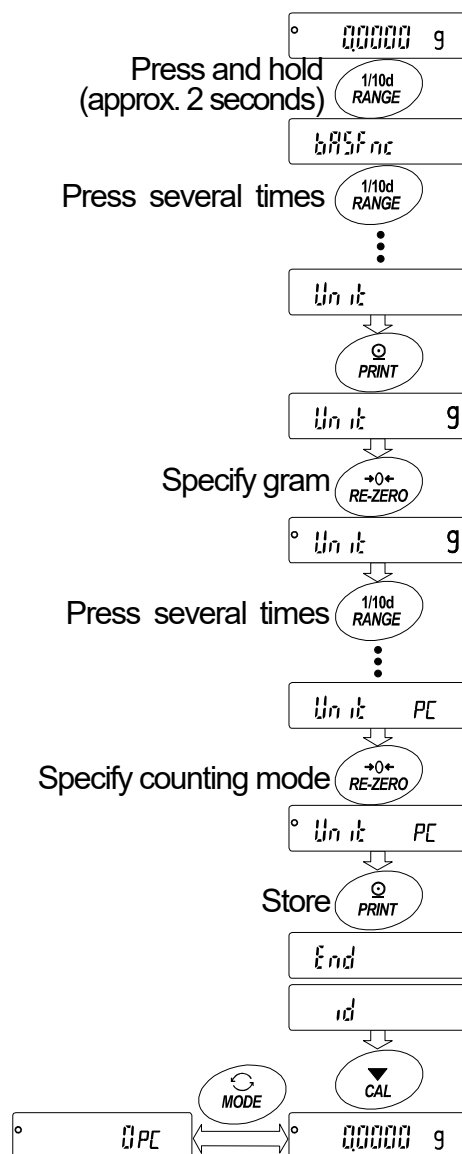

## 6. Weighing

### Precautions for the weighing operation


- Press the **[RE-ZERO]** key each time, before placing a sample on the weighing pan, to prevent possible errors.
- Place a sample in the center of the weighing pan gently.
- Temperature changes during measurement may cause weighing errors.
- Shorten the operation time as much as possible. (Opening and closing door, placing and removing sample)
- Material with an electrostatic charge or that is magnetic may cause weighing errors.
- Use a pair of tweezers to avoid a temperature change due to having your hand in the weighing chamber.
- Do not drop things on the pan, or place a weight on the pan that is beyond the weighing range.
- Keep the area clean and dry.
- Do not press keys with a sharp instrument (such as a pencil or ball point pen).
- For precision weighing, keep the AC adapter connected to the balance.
- Adjust the sensitivity periodically to maintain weighing accuracy. Refer to "[8. Sensitivity Adjustment / Calibration Test](#)".
- For the **BM-5**, **BM-5D**, **BM-20** and **BM-22**, perform pre-weighing of the container (tare) so that precision weighing can be performed.
- Consider section "[3. Installing the Balance and Precautions](#)" for the weighing operation.

### 6.1. Basic Operation (Gram Mode)

Refer to "[4. Display Symbols and Key Operation](#)" before operation.

**Note** When turning the balance on with a tare weight placed on the pan, the balance automatically displays the zero display.

- 1 Press the **[ON:OFF]** key to turn the display on.
- 2 Press the **[MODE]** key to select a unit, if necessary.
- 3 Place a container (tare) on the weighing pan, if necessary.  
Press the **[RE-ZERO]** key to cancel the tare weight.  
Then zero is displayed.  
Tare : A vessel placed on the pan, but not to be included in the weighing data.  
Example: Container.
- 4 Place a sample on the pan or in the container.  
Close the door.
- 5 Wait for the stabilization indicator **[O]** to be displayed, and then read the value.
- 6 Remove the sample and container from the pan.

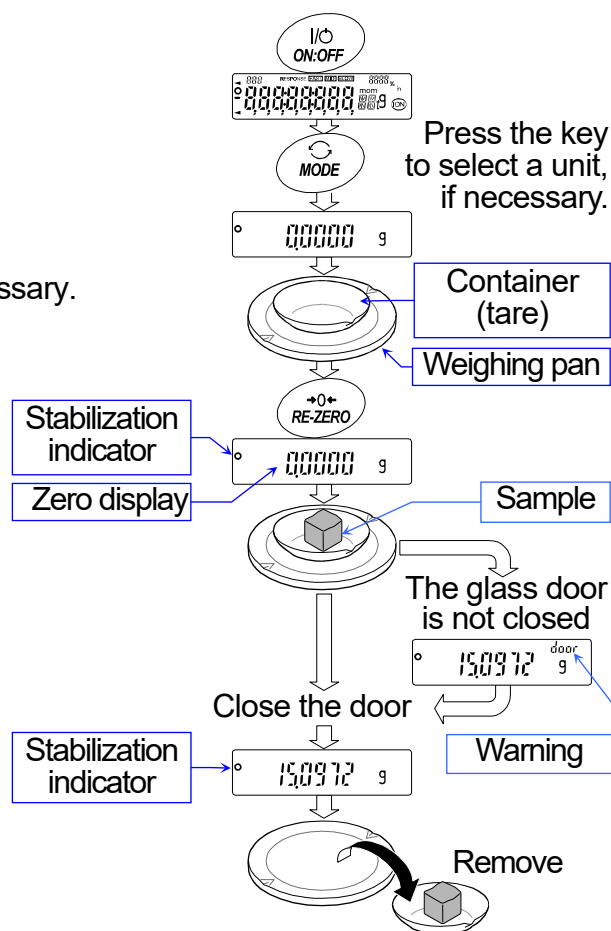

## Zero operation, tare operation and weighing range

### □ When weighing is started

The balance will determine the reference zero point when the **[ON:OFF]** key is pressed to enter the weighing mode.

Depending on the load condition at that time, the balance will automatically judge whether to perform zero or tare operation. The condition for determining which is used is "power on zero range", and when power on zero range is exceeded, the tare operation is performed.

### □ Re-zero operation

By pressing the **[RE-ZERO]** key, the display can be set to zero.

The re-zero operation with the **[RE-ZERO]** key will automatically judge whether to perform zero or tare operation.

The condition for determining which is used is "zero range", and when zero range is exceeded, the tare operation is performed.

### □ Weighing range

The range that the balance can weigh varies depending on the model.

When the total amount (net weight + tare weight) displayed for each model exceeds the maximum display, **[E]** is displayed to indicate that the weighing range is exceeded.

When exceeded in the negative direction, **[-E]** is displayed.


| Model         | Power on zero range | Zero range               | -E display range        |
|---------------|---------------------|--------------------------|-------------------------|
| <b>BM-5</b>   | Approx. $\pm 0.1$ g | Approx. -3.0 g to +0.1 g | Less than approx. -3 g  |
| <b>BM-5D</b>  |                     |                          |                         |
| <b>BM-20</b>  | Approx. $\pm 2.2$ g | Approx. -3.0 g to +0.4 g | Less than approx. -3 g  |
| <b>BM-22</b>  |                     |                          |                         |
| <b>BM-252</b> | Approx. $\pm 25$ g  | Approx. -37 g to +5.0 g  | Less than approx. -37 g |
| <b>BM-200</b> | Approx. $\pm 22$ g  | Approx. -37 g to +4.4 g  |                         |
| <b>BM-300</b> | Approx. $\pm 32$ g  | Approx. -37 g to +6.4 g  |                         |
| <b>BM-500</b> | Approx. $\pm 52$ g  | Approx. -37 g to +10.4 g |                         |

## 6.2. Smart Range

- The smart range for the **BM-5D** and **BM-22** consists of the standard range and precision range (high resolution).
- Smart range function
  - The range switches automatically, depending on the value displayed.
  - Pressing the **RE-ZERO** key allows weighing in the precision range, regardless of the tare value.
  - The range can be fixed to the standard range by pressing the **RANGE** key.

### Example

- 1 Press the **RE-ZERO** key to use the precision range.  
The balance displays zero.
- 2 Place a container (tare) on the weighing pan.  
When the value of the container (tare) exceeds the precision range, the balance automatically switches to the standard range.
- 3 Press the **RE-ZERO** key to use the precision range.  
The balance displays zero.
- 4 Place the weighing sample in the container (tare).  
When the sample is within the precision range, the weighing value can be read in high resolution.

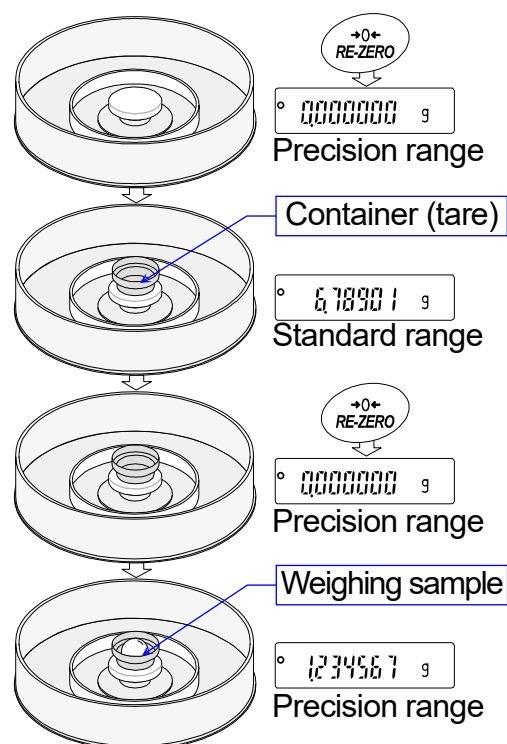

### Precision range and standard range

| Model        | Unit      | Precision range                                 | Standard range                                    |
|--------------|-----------|---|---|
| <b>BM-5D</b> | Milligram | 0.000 <sup>m</sup> g to 2100.009 <sup>m</sup> g | 2100.01 <sup>m</sup> g to 5200.08 <sup>m</sup> g  |
|              | Gram      | 0.000000 g to 2.100009 g                        | 2.10001 g to 5.20008 g                            |
| <b>BM-22</b> | Milligram | 0.000 <sup>m</sup> g to 5100.009 <sup>m</sup> g | 5100.01 <sup>m</sup> g to 22000.08 <sup>m</sup> g |
|              | Gram      | 0.000000 g to 5.100009 g                        | 5.10001 g to 22.00008 g                           |

Refer to page [20](#) and [21](#) for other units.

## 6.3. Counting Mode (PC)

- This is the mode to determine the number of objects in a sample based on the standard sample unit mass. The unit mass means an average mass of the samples. The smaller the variation in the samples is, the more accurate the count will be. The balance is equipped with the Automatic Counting Accuracy Improvement (ACAI) function to improve the counting accuracy.

### Notes

- Use samples with a unit mass of 1 mg or more for counting.
- If the sample unit mass variable is too large, it may cause a counting error.
- To improve the counting performance, use the ACAI function frequently or divide the samples into several groups and count each group.

### Selecting the counting mode

- 1 Press the **[MODE]** key to select the unit **[PC]** (counting mode).

### Storing a sample unit mass (Weighing input mode)

- 2 Press the **[RANGE]** key to enter the sample unit mass storing mode.

- 3 Select the number of samples using the **[RANGE]** key.  
It may be set to 10, 25, 50 or 100.

**Note** A greater number of samples will yield more accurate counting result.

- 4 Place a container (tare) on the weighing pan, if necessary.  
Press the **[RE-ZERO]** key to cancel the weight (tare).

The number specified in Step 3 appears.

Example: **[25] [PC]** is displayed if 25 is selected in Step 3.

- 5 Place the number of samples specified on the pan.  
In this example, 25 pieces.


- 6 Wait for the stabilization indicator to be displayed.  
Press the **[PRINT]** key to calculate and store the unit mass. Then the balance displays **[25] [PC]** and is set to count samples with this unit mass. To improve the accuracy of the unit mass, proceed to Step 8.

### Notes

- If the balance judges that the mass of the samples is too light to acquire accurate weighing, it displays an error requiring the addition of more samples to the specified number.  
Example: **[50] - PC** appears, requiring 25 more samples. Add 25 samples and press the **[PRINT]** key. When the unit mass is stored correctly, the balance proceeds to the counting mode.
- If the balance judges that the mass of the samples is too light (under 0.0001g) and cannot be stored as the unit mass, it displays **[Lo]**.
- The sample unit mass is stored in non-volatile memory, and is maintained even if the AC adapter is removed.

### Counting operation

- 7 Place the samples to be counted on the pan.  
Read the result and remove the samples from the pan.

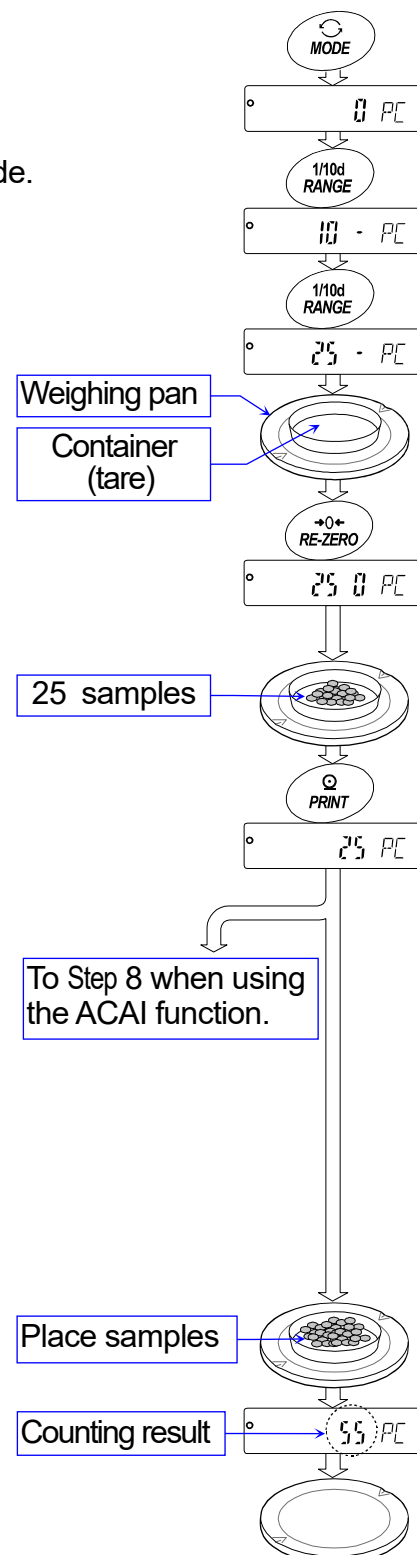


## Counting mode using the ACAI function

The ACAI is a function that improves the accuracy of the unit mass automatically by increasing the number of samples as the counting process.

**ACAI:** Automatic Counting Accuracy Improvement

Proceed to Step 8 after storing a unit mass on the previous page.

- 8 If a few more samples are added, the processing indicator turns on. To prevent an error, add three or more. The processing indicator does not turn on if overloaded. Try to add the same number of samples as displayed.
- 9 The balance re-calculates the unit mass while the processing indicator is blinking. Do not touch the balance or samples on the pan until the processing indicator turns off.
- 10 Counting accuracy is improved when the processing indicator turns off.  
Each time the above operation is performed, a more accurate unit mass will be obtained. There is no definite upper limit to the ACAI range for the number of samples exceeding 100. Try to add the same number of samples as displayed.
- 11 Remove all the samples used in ACAI and proceed with the counting operation using the improved unit mass.

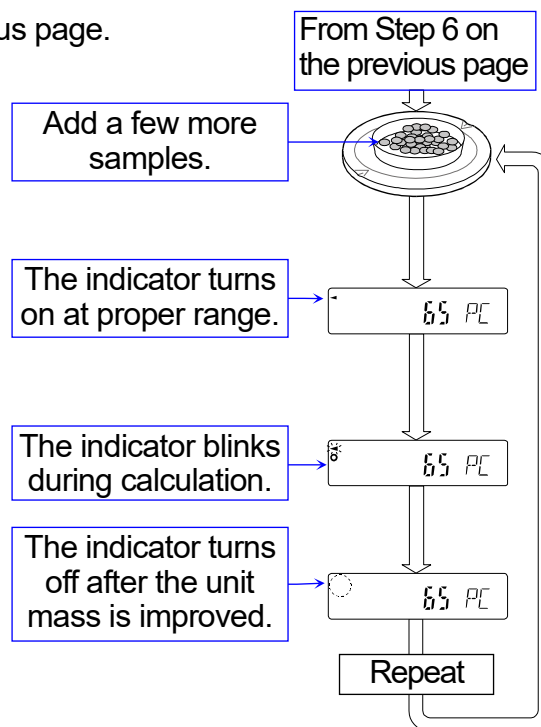

## 6.4. Percent Mode (%)

The percent mode displays the weight value in percentage compared to a 100% reference mass and is used for target weighing or checking the sample variance.

### Selecting the percent mode

- 1 Press the **[MODE]** key to select the unit **[%]** (Percent mode).  
If the percent mode cannot be selected, refer to "5. Weighing Units".

### Storing the 100% reference mass

- 2 Press the **[RANGE]** key to enter the 100% reference mass storing mode.  
The balance displays **100.0 %**.
- 3 Place a container (tare) on the weighing pan, if necessary.  
Press the **[RE-ZERO]** key to cancel the weight (tare).  
The balance displays **100.0 %**.
- 4 Place the sample to be set as the 100% reference mass on the pan or in the container.
- 5 Press the **[PRINT]** key to store the reference mass.  
The balance displays **100.00 %**.

### Notes


- The readability depends on the 100% mass.

| BM-252, BM-200<br>BM-300, BM-500 |             | BM-5, BM-5D<br>BM-20, BM-22 |             |
|----------------------------------|-------------|-----------------------------|-------------|
| 100% mass                        | Readability | 100% mass                   | Readability |
| 0.0100 g to 0.0999 g             | 1 %         | 0.00100 g to 0.0099 g       | 1 %         |
| 0.1000 g to 0.9999 g             | 0.1 %       | 0.01000 g to 0.0999 g       | 0.1 %       |
| 1.0000 g to                      | 0.01 %      | 0.10000 g to                | 0.01 %      |

- If the balance judges that the mass of the sample is too light (under 0.01g) to be used as a reference, it displays **[Lo]**.
  - The 100% reference mass can be stored in the non-volatile memory and is maintained even if the AC adapter is removed.
- 6 Remove the sample.

### Reading the percentage

- 7 Place a sample to be compared to the reference mass on the pan. The displayed percentage is based on the 100% reference mass.

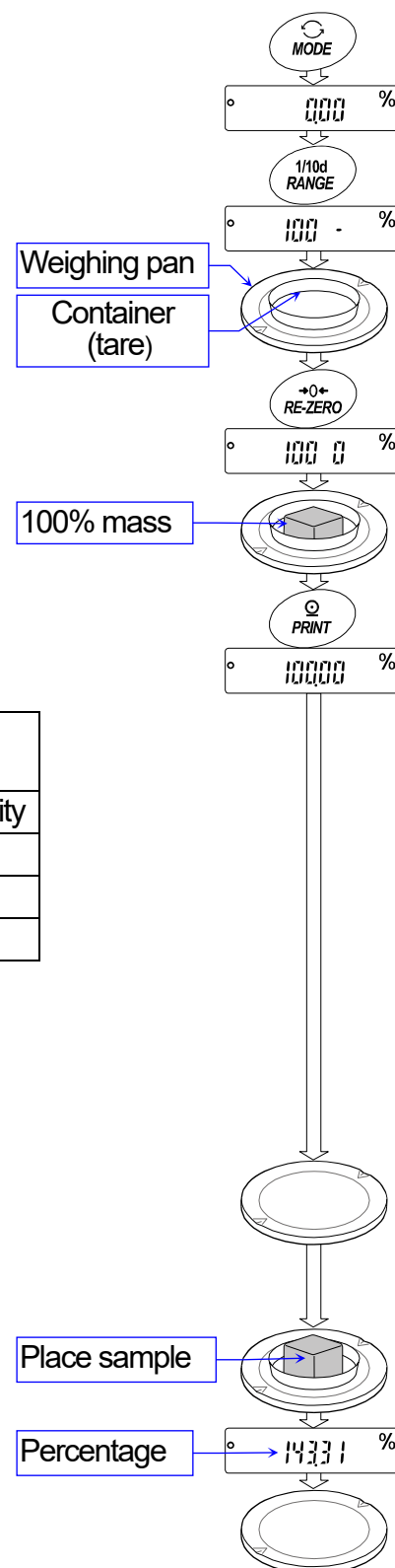

## 6.5. Ionizer (Neutralization Device)

The **BM** series is equipped with four ionizer electrode units to neutralize static electricity from a charged weighing sample.


Neutralizing static electricity before weighing will improve stability, reducing errors.

Neutralizing ion : The discharge electrodes of the ionizer continuously generate bipolar ions by corona discharge. Applying these ions to the weighing sample neutralizes the static charge.

Static electricity : In general, when the ambient humidity is less than 45% RH, nonconductors such as powders, paper and plastics easily become charged with static electricity. The influence of the static electricity may cause a weighing error of several milligrams. The ionizer effectively neutralizes the electrical charge.

### Operation

- 1 Place the sample to be weighed on the center (of the metallic circle) of the separation plate to neutralize the static charge.
  - 2 Press the **ION** key to start the neutralization.  
The **ION** mark and the LED blink.  
The neutralization will stop after the preset time of "Neutralization time ( **ION** )" in the function table (the factory setting is 3 seconds).
- The readability of the **BM-5**, **BM-5D**, **BM-20**, **BM-22** and **BM-252** is switched to 0.1 mg while performing the neutralization.
  - Pressing the **ION** key during neutralization stops the neutralization.

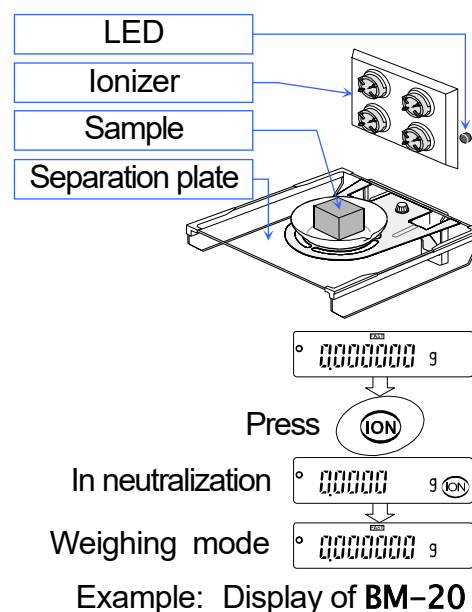

### Notes

- **Keep the weighing sample away from the electrodes while neutralizing. Placing the sample too close to the electrodes may cause the sample to become charged.**
- **Remove any obstacle between electrodes and the weighing sample.**

### 6.5.1. Maintenance of the Electrode Units


- When using the ionizer for a long time, dust and stains may stick to the electrodes. Clean them periodically to maintain performance.
- When needles of the electrodes are worn down and neutralization ability is not restored after cleaning them, replace all four electrodes with new ones. The standard life of an electrode unit is approximately 10000 hours.

### Replacing the electrode units

- 1 Turn the electrode units 45 degrees counterclockwise and remove them from the ionizer.
- 2 Replace all four units with new ones.

**Note** Do not replace, remove and clean the electrodes while the ionizer is in operation as there is a danger of electrical shock.

- 3 Turn the electrode units 45 degrees clockwise and insert them into the ionizer.

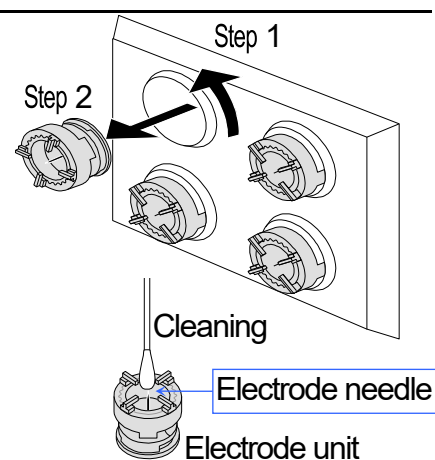




## 6.5.2. Controlling the Ionizer Externally

It is possible to control the BM ionizer function by connecting the remote controller (AD-8922A) or the foot switch (AX-SW137-PRINT, AX-SW137-REZERO).

- When specifying "on" to "AD-8922A control (8922 on)", the balance assigns the function of the **ION** key to the **MODE** key on the AD-8922A.

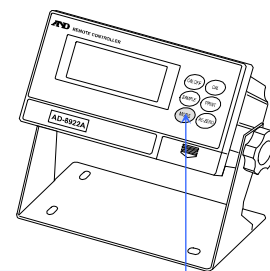

Assign the **ION** key to the **MODE** key on the AD-8922A.

- When specifying "External control input (E-5)" in "Ionizer function (ionFnc)", the following functions can be assigned to AX-SW137-PRINT (option) or AX-SW137-REZERO (option) terminals.


| Ionizer function | External control input | AX-SW137-PRINT   | AX-SW137-REZERO    |
|------------------|------------------------|------------------|--------------------|
| ▪ ionFnc         | E-5 0                  | <b>PRINT</b> key | <b>RE-ZERO</b> key |
| ionFnc           | E-5 1                  | <b>PRINT</b> key | <b>ION</b> key     |
| ionFnc           | E-5 2                  | <b>ION</b> key   | <b>RE-ZERO</b> key |

- Factory settings

- By using a  $\phi 3.5$  stereo plug MP-013LC (Marushin Electric Mfg. Co., Ltd.) or an equivalent product, you can create your own switch.  
(To use a switch, you need to solder the  $\phi 3.5$  stereo plug and the switch that you obtain yourself. )


External contact input terminal

Compatible plug:  $\phi 3.5$  mm 3-pole stereo plug  
MP-013LC or the equivalent

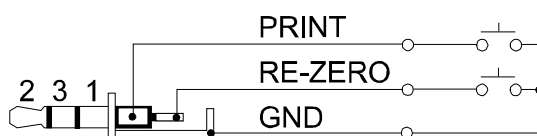

Example of external contact input terminal

### External contact input circuit

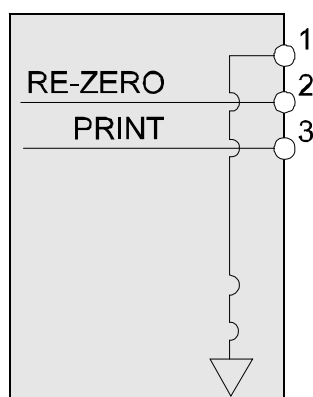

### Pin assignments

| Pin No. | Description                    |
|---------|--------------------------------|
| 1       | GND                            |
| 2       | RE-ZERO External contact input |
| 3       | PRINT External contact input   |

When pin 1 and pin 3 or pin 1 and pin 2 are shorted 100 ms or longer, the switch can operate the balance in the same way as the balance **PRINT** key or the **RE-ZERO** key, respectively.

- AX-SW137-PRINT or AX-SW137-REZERO (optional foot switches sold separately) can be used.

## 6.6. Monitoring the Clock and Environmental Sensors

In the weighing mode, the built-in clock and environmental sensors (temperature, humidity, barometric pressure) can be monitored and temperature data can be output in the function settings.

### Specifications of the environmental sensors


| Sensor              | Resolution | Measurement range   | Applicable range |
|---------------------|------------|---------------------|------------------|
| Temperature         | ±1.5 °C    | 5 °C to 40 °C       | 5 °C to 40 °C    |
| Humidity            | ±10 %      | 0 % to 100 %        |                  |
| Barometric pressure | ±10 hPa    | 300 hPa to 1100 hPa |                  |

### Notes

- The environmental sensors are built into the balance casing. Therefore, these values are not in accordance with the environmental condition.
- Refer to "20. Specifications" for the balance operating conditions.

### Operation

- Press the **SELECT** key in the weighing mode. The current time, date and sensor values (temperature, humidity, barometric pressure) are displayed for several seconds in order. The balance returns to the weighing mode automatically.
- When pressing the **SELECT** key while the current data is displayed, the display proceeds to the next data.
- When pressing the **CAL** key while the current data is displayed, the balance returns to the weighing mode.

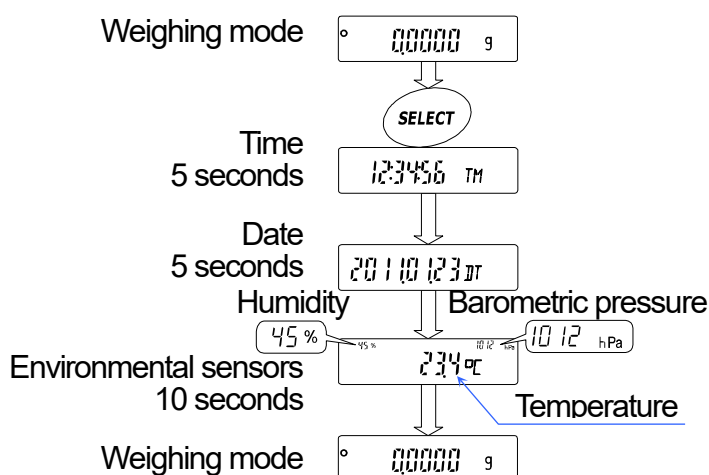

### Notes

- When adjusting the built-in clock, refer to "10.7. Clock and Calendar Function".
- When adding the time and date to the output data, refer to "Time/Date output (5-td)" of "10.2. Details of the Function Table".
- When adding temperature data to the output data, refer to "Temperature output (5-tp)" of "10.2. Details of the Function Table".

## 6.7. Repeatability Test

Repeatability is an index of variation in weighing values when the same mass is repeatedly loaded and unloaded. Standard deviation is generally used for the indicator of repeatability.

The repeatability test measures the internal mass 10 times and displays the standard deviation.

Use the repeatability test function to inspect the performance of the balance when changing the environment or when moving the balance to a new location.

Example: "Standard deviation = 0.2 mg" means that the results of repeated measurements of the same mass fall within the range of  $\pm 0.2$  mg at a frequency of about 68%.

### Notes

- When the password function is used, the repeatability test function is available only when logged in as an administrator (ADMIN).
- The repeatability test uses the internal mass. Therefore, the obtained standard deviation may disagree with that of the repeatability of "20. Specifications". Regard the standard deviation as a reference.


| Model                          | Internal mass |
|--------------------------------|---------------|
| BM-5, BM-5D                    | Approx. 5 g   |
| BM-20, BM-22                   | Approx. 20 g  |
| BM-252, BM-200, BM-300, BM-500 | Approx. 190 g |

- Press and hold the **RANGE** key (for approx. 4 seconds) until **REP TEST** is displayed, and then release the key.
- When **REP TEST** is displayed, the measurement is automatically started, **REP** blinks and the measurement count (0 to 10) is displayed.  
To cancel the current measurement, press the **CAL** key.

**Note** Avoid vibration and breezes for precise measurement during this test.



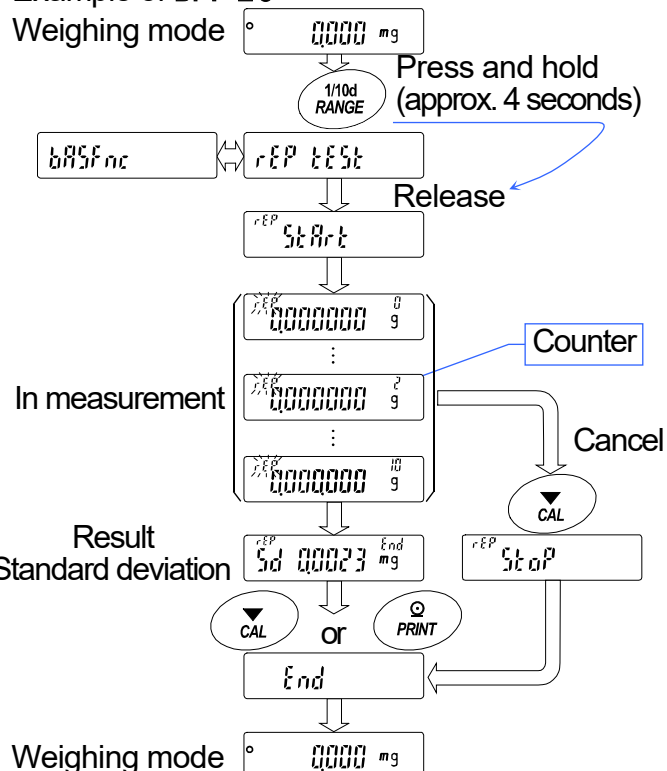

- At the end of the measurement, the standard deviation of the repeatability is displayed.
- Press the **CAL** key or **PRINT** key to return to the weighing mode.

### Example of BM-20




## 7. Response Adjustment

This function stabilizes the weight value by reducing the influence on weighing that is caused by drafts and/or vibration at the place where the balance is installed. The adjustment of the function can be performed by analyzing the environment automatically or hand-operation. The state of the function has three stages as follows :

| Indicator | Parameter | Response   | Stability   |
|-----------|-----------|--|---|
| FAST      | [End 0]   | Fast response,  | Sensitive value  |
| MID.      | [End 1]   |  |   |
| SLOW      | [End 2]   | Slow response,  | Stable value     |

Response indicator

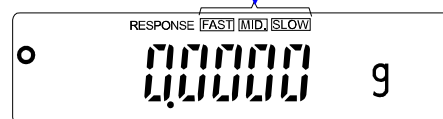


### 7.1. Automatic Response Adjustment

This function automatically updates the response adjustment by analyzing the influence of the environment using the internal mass.

- 1 Press and hold the [MODE] key (for approx. 2 seconds) until [RESPONSE] is displayed, and then release the key.
- 2 The balance automatically sets the response characteristic.

**Caution** Do not allow vibration or drafts to affect the balance during adjustment.

- 3 After automatic adjustment, the balance displays [End], returns to the weighing mode and displays the updated response indicator for about thirty seconds.

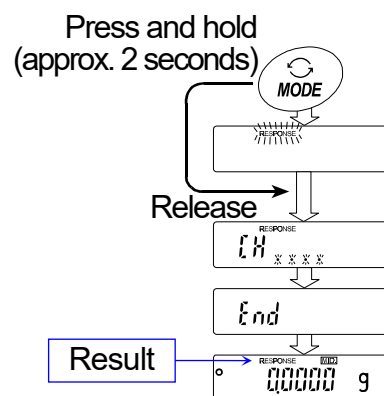

#### Notes

- If the automatic response adjustment fails, the balance displays [CH nG]. Check the ambient conditions such as breeze and vibration, also check the weighing pan. Then, perform adjustment again. Press the [CAL] key to return to the weighing mode.
- If there is anything on the weighing pan, the balance displays [CH E]. Remove the substance from the pan. Press the [CAL] key to return to the weighing mode.
- If the automatic response adjustment is not helpful, try ["7.2. Manual Response Adjustment"](#).

## 7.2. Manual Response Adjustment


- 1 Press and hold the **MODE** key (for approx. 2 seconds) until **RESPONSE** is displayed, and then release the key. Press the **MODE** key again immediately.
- 2 Select a setting for the response adjustment using the **MODE** key. Either **FAST**, **MID.** or **SLOW** can be selected.
- 3 The balance displays **End**, returns to the weighing mode and displays the updated response indicator for about thirty seconds.

### Notes

- When the response adjustment is selected, the "Condition (*Cond*)" parameter of "Environment, Display (*bRSFnc*)" in the function table will be changed accordingly as shown below.

| Display | Cond(Condition) |
|---------|-----------------|
| FAST    | 0               |
| MID.    | 1               |
| SLOW    | 2               |

- The response adjustment setting can be changed by the "Condition (*Cond*)" parameter of "Environment, Display (*bRSFnc*)" in the function table. For details, refer to "10. Function Table".

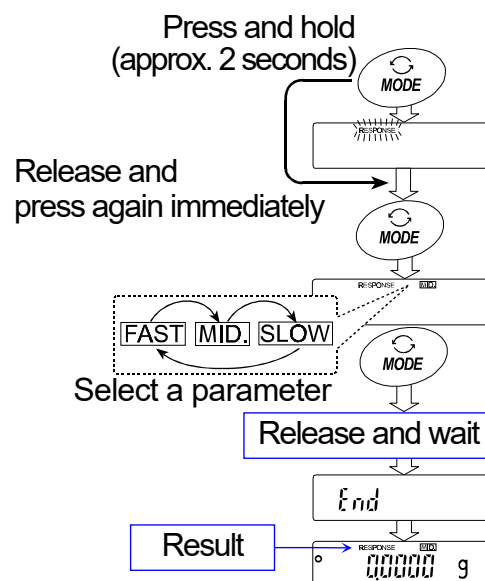

## 8. Sensitivity Adjustment / Calibration Test

Since the balance's resolution is high, weighing values may change due to gravity and daily environmental changes. It is necessary to perform sensitivity adjustment with the weight in order to keep the weighing values from changing even if gravity or the environment changes.

Adjust the balance if it is installed for the first time or relocated, or when the weighing values change significantly in daily inspection.

Sensitivity adjustment means to adjust the weighing value of the balance using the reference weight or internal mass. Calibration test means to weigh with the reference weight and compare how much the result deviates from the reference value. (Calibration test does not perform adjustment.)

### Sensitivity adjustment

Automatic sensitivity adjustment — Automatically adjusts the balance using the internal mass depending on the temperature change of the operating environment.

Sensitivity adjustment using the internal mass — Using the internal mass, adjusts the balance with a single touch.

Sensitivity adjustment using an external weight — Using an external weight, adjusts the balance.

### Calibration test

Calibration test using the internal mass — Checks the accuracy of weighing using the internal mass and outputs the result.

\*No adjustment is made.

Calibration test using an external weight — Checks the accuracy of weighing using an external weight (target weight) and outputs the result.

\*No adjustment is made.

### Notes

- Do not allow vibration or drafts to affect the balance during sensitivity adjustment.
- When using the data output for GLP / GMP using the RS-232C interface, set "GLP output (INF0)" of the function table. Refer to "10. Function Table". The time and date can be added to the GLP report. If the time and date are incorrect, refer to "10.7. Clock and Calendar Function" and adjust them.
- Calibration test is available only when "GLP output (INF0)" is set to "1", "2" or "3".
- The sensitivity adjustment and calibration test data can be stored in memory. When using memory, set "Data memory (dRtR)" of the function table. Refer to "12. Data Memory" for details.

### Note on using an external weight

- The accuracy of an external weight can influence the accuracy of weighing.
- Select a weight for sensitivity adjustment and calibration test from the following table.

| Model         | Usable weight |       |               |       |               | Adjustable range       |
|---------------|---------------|-------|---------------|-------|---------------|------------------------|
| <b>BM-5</b>   | 1 g           | 2 g   | <b>5 g*</b>   |       |               | -3.000 mg to +3.099 mg |
| <b>BM-5D</b>  | 1 g           | 2 g   | <b>5 g*</b>   |       |               |                        |
| <b>BM-20</b>  | 1 g           | 2 g   | 5 g           | 10 g  | <b>20 g*</b>  |                        |
| <b>BM-22</b>  | 1 g           | 2 g   | 5 g           | 10 g  | <b>20 g*</b>  |                        |
| <b>BM-252</b> | 10 g          | 20 g  | 50 g          | 100 g | <b>200 g*</b> | -15.00 mg to +15.99 mg |
| <b>BM-200</b> | 50 g          | 100 g | <b>200 g*</b> |       |               | -30.0 mg to +30.9 mg   |
| <b>BM-300</b> | 50 g          | 100 g | <b>200 g*</b> | 300 g |               |                        |
| <b>BM-500</b> | 50 g          | 100 g | <b>200 g*</b> | 300 g | 500 g         |                        |

**Bold type\*:** Factory settings. The weight value can be adjusted within the range above.

## About the internal mass

- The internal mass may change due to corrosion or other damage caused by the operating environment, or due to aging. Check the internal mass periodically and correct the internal mass value if necessary. Refer to "[8.6. Correcting the Internal Mass Value](#)".

## Display



This indicator means that the balance is measuring sensitivity adjustment data. Do not allow vibration or drafts to affect the balance while the indicator is displayed.

## 8.1. Automatic Sensitivity Adjustment

### Automatic sensitivity adjustment due to changes in temperature

This function automatically adjusts the balance when the balance detects an ambient temperature change. If GLP output is selected in the function table, the balance outputs the sensitivity adjustment report or stores the data in memory. Automatic sensitivity adjustment functions even if the display is turned off (standby mode). Refer to "[9.1. Permit or Inhibit](#)" for the operation.


### Caution

- **If the balance detects a load on the weighing pan, it will determine that it is in use and will not perform automatic sensitivity adjustment. The criteria for performing automatic sensitivity adjustment are as follows.**


#### The criteria for performing automatic sensitivity adjustment

|                                   |                 |
|-----------------------------------|-----------------|
| BM-5 / BM-5D / BM-20 / BM-22      | Less than 0.5 g |
| BM-252 / BM-200 / BM-300 / BM-500 |                 |

- **To maintain the correct sensitivity adjustment of the balance, do not place anything on the weighing pan while not in use.**
- **When performing a long term weighing with something placed on the pan, installing the balance in a system or considering the importance of continuous data, turn off automatic sensitivity adjustment.**

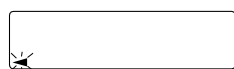

Indicates that the balance detects changes in ambient temperature and automatic sensitivity adjustment will start. If the balance is not used for several minutes with this indicator (◀) blinking, the balance performs automatic sensitivity adjustment. The blinking duration depends on the operating environment.

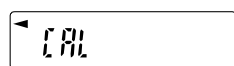


Indicates that the balance is measuring sensitivity adjustment data. Do not allow vibration or drafts to affect the balance while this indicator is displayed. After sensitivity adjustment, the balance returns to indicate the previous display.

- **The balance can be used while the indicator is blinking. However, for accurate weighing, ensure the pan is empty and allow the balance to complete automatic sensitivity adjustment.**

## 8.2. Sensitivity Adjustment Using the Internal Mass

- This function adjusts the balance using the internal mass.
  - The only operation required is to press the **[CAL]** key.
- 1 Connect the AC adapter and warm up the balance for at least one hour with nothing on the weighing pan.
  - 2 Press the **[CAL]** key to display **[CAL in]**.
  - 3 The balance performs sensitivity adjustment using the internal mass. Do not allow vibration or drafts to affect the balance.
  - 4 If GLP output is set, **[GLP]** is displayed, the sensitivity adjustment report is output to the RS-232C interface and is stored in memory. Refer to "GLP output (inFo)" and "Data memory (dAtA)" of the function table in "[11.2. GLP Report](#)" and "[12. Data Memory](#)".  
**[End]** is displayed after sensitivity adjustment.
  - 5 The balance will automatically return to the weighing mode after sensitivity adjustment.
  - 6 Confirm weighing accuracy using calibration test (**[CAL in]**).

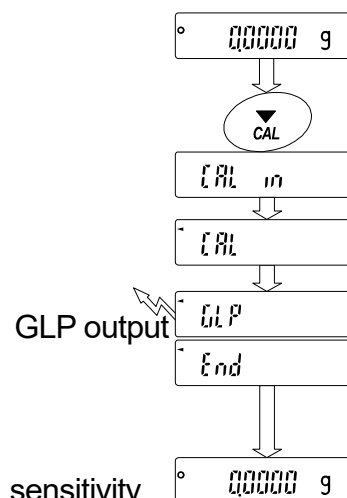



### 8.3. Calibration Test Using the Internal Mass

- This function checks the accuracy of weighing using the internal mass.

1 Connect the AC adapter and warm up the balance for at least one hour with nothing on the weighing pan.

2 Press and hold the **[CAL]** key until **[CAL in]** is displayed, and then release the key.

3 The balance measures the zero point.  
Prevent vibration and drafts to affect the balance.

4 The measured zero point data is displayed.

5 The balance measures the internal mass.  
Prevent vibration and drafts to affect the balance.


6 The value of the internal mass is displayed.  
The normal range of the value is as follows:

| Model  | Internal mass | Normal range |
|--|---------------|--------------|
| <b>BM-5, BM-5D</b>                             | 5.00000 g     | ±0.02 mg     |
| <b>BM-20, BM-22</b>                            | 20.00000 g    | ±0.02 mg     |
| <b>BM-252, BM-200</b><br><b>BM-300, BM-500</b> | 200.0000 g    | ±0.2 mg      |

7 **[GLP]** is displayed and the calibration test report is output to the RS-232C interface and is stored in memory. Refer to "GLP output (inFa)" and "Data memory (dAtA)" of the function table.

**[End]** is displayed after the calibration test.

8 The balance automatically returns to the weighing mode.

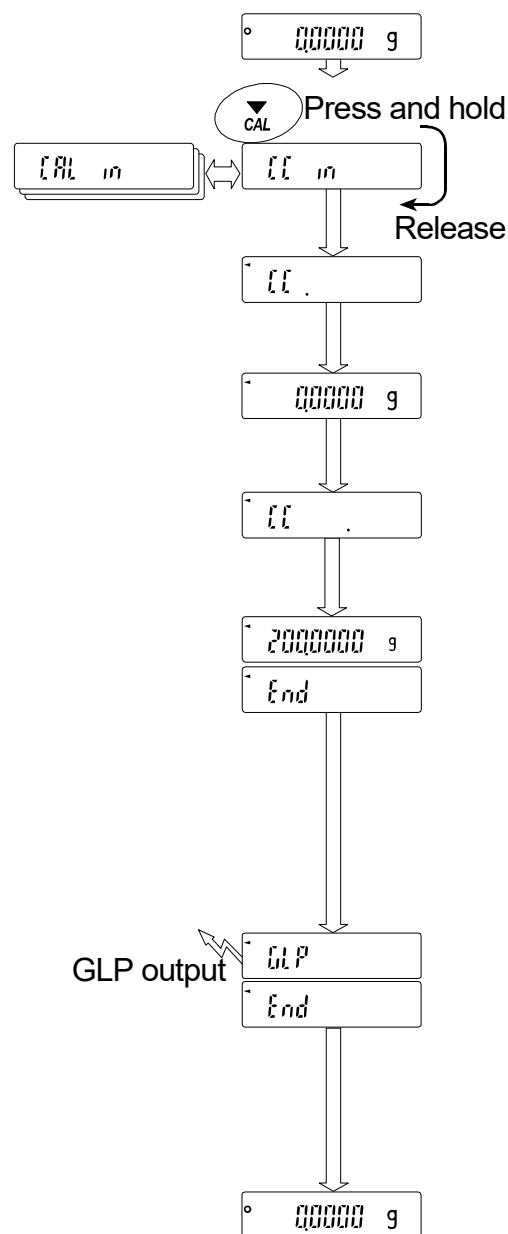


## 8.4. Sensitivity adjustment Using an External Weight

- This function adjusts the balance using an external weight.

**Note** Automatic sensitivity adjustment (due to changes in temperature) is enabled in the factory setting. Therefore, even after sensitivity adjustment using an external weight, the balance may perform automatic sensitivity adjustment whenever it detects changes in temperature.

When considering the importance of continuous data or managing the balance using the external weight only, refer to "9.1. Permit or Inhibit" to set the function switch to "inhibit".


- 1 Connect the AC adapter and warm up the balance for at least one hour with nothing on the weighing pan.
- 2 Press and hold the **[CAL]** key until **[CAL out]** is displayed, and then release the key.
- 3 When **[CAL 0]** is displayed, proceed as follows:
  - If you want to change the weight, press the **[RANGE]** key and proceed to Step 4.
  - If you use the weight value stored in the balance, proceed to Step 5.
- 4 Specify the weight value as follows:
  - [RANGE]** key ..... To switch blinking digits.
  - [RE-ZERO](+)** key ..... To select the weight or adjust the weight
  - [MODE](-)** key ..... value. Refer to page 38.
  - [PRINT]** key ..... To store the new weight value. Even if the AC adapter is removed, the data is maintained in non-volatile memory.
  - [CAL]** key ..... To cancel the operation and return to **[CAL 0]**.
- 5 Confirm that there is nothing on the pan and press the **[PRINT]** key. The balance measures the zero point. Do not allow vibration or drafts to affect the balance. The balance displays the weight value.
- 6 Place the displayed weight on the pan and press the **[PRINT]** key. The balance measures the external weight. Do not allow vibration or drafts to affect the balance.
- 7 **[End]** is displayed. Remove the weight from the pan.
- 8 If GLP output is set, **[GLP]** is displayed and the sensitivity adjustment report is output and stored. Refer to "GLP output (info)" and "Data memory (data)" of the function table in "11.2. GLP Report" and "12. Data Memory".
- 9 The balance will automatically return to the weighing mode.
- 10 Place the external weight on the pan and confirm that the value displayed is correct. If it is not within the range, check the ambient conditions such as breeze and vibration, and also check the weighing pan. Then, repeat Steps 2 to 10.

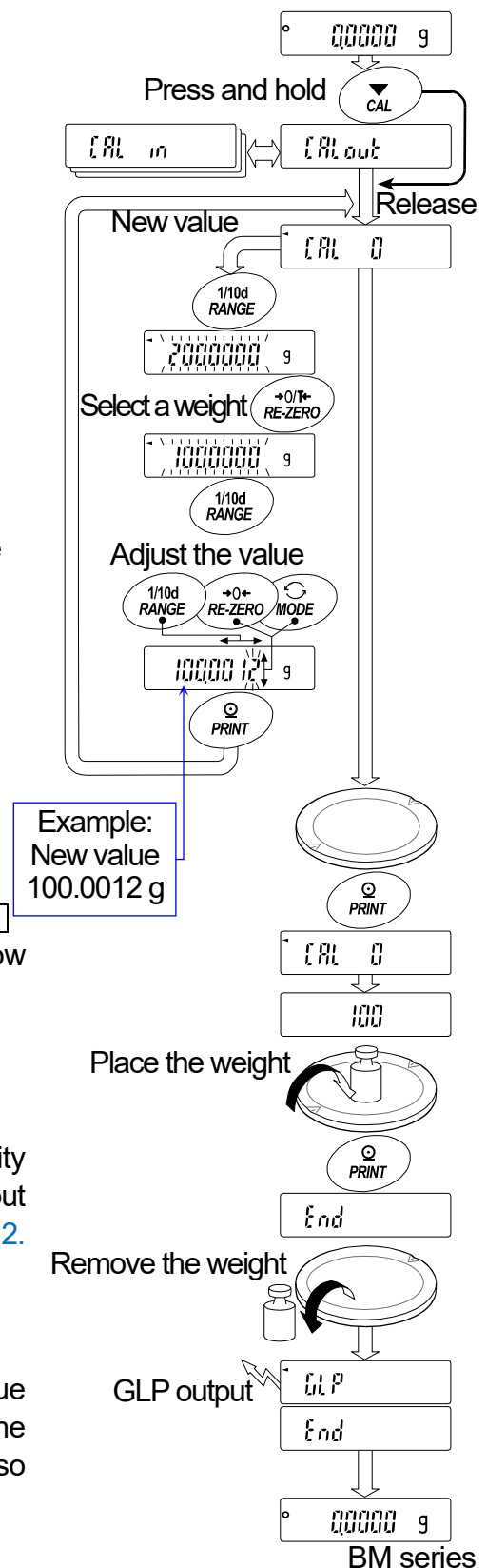

## 8.5. Calibration Test Using an External Weight


- This function checks the accuracy of weighing using an external weight.
- This is available only when the "GLP output ( *inf*o )" parameter is set to "1", "2" or "3".
- Calibration test does not perform adjustment.

- 1 Connect the AC adapter and warm up the balance for at least one hour with nothing on the weighing pan.
- 2 Press and hold the **[CAL]** key until **[EE out]** is displayed, and then release the key.
- 3 When **[EE 0]** is displayed, proceed as follows:
  - If you want to change the target weight, press the **[RANGE]** key and proceed to Step 4. A list of usable weights is on page 38.
  - If the current target weight value is used, proceed to Step 5.
- 4 Specify the target weight value as follows:
  - [RANGE]** key ..... To switch blinking digits.
  - [RE-ZERO](+)**key ... To select the target weight or adjust the weight value. Refer to page 38.
  - [MODE](-)**key ..... weight value. Refer to page 38.
  - [PRINT]** key ..... To store the new target weight value. Even if the AC adapter is removed, the data is maintained in non-volatile memory.
  - [CAL]** key ..... To cancel the operation and return to **[EE 0]**.
- 5 Confirm that there is nothing on the pan and press the **[PRINT]** key. The balance measures the zero point. Do not allow vibration or drafts to affect the balance. The balance displays the target weight value.
- 6 Place the displayed target weight on the pan and press the **[PRINT]** key. The balance measures the target weight. Do not allow vibration or drafts to affect the balance.
- 7 **[End]** is displayed. Remove the weight from the pan.
- 8 **[GLP]** is displayed and the calibration test report is output and stored. Refer to "GLP output ( *inf*o )" and "Data memory ( *data* )" of the function table in "11.2. GLP Report" and "12. Data Memory".
- 9 The balance will automatically return to the weighing mode.

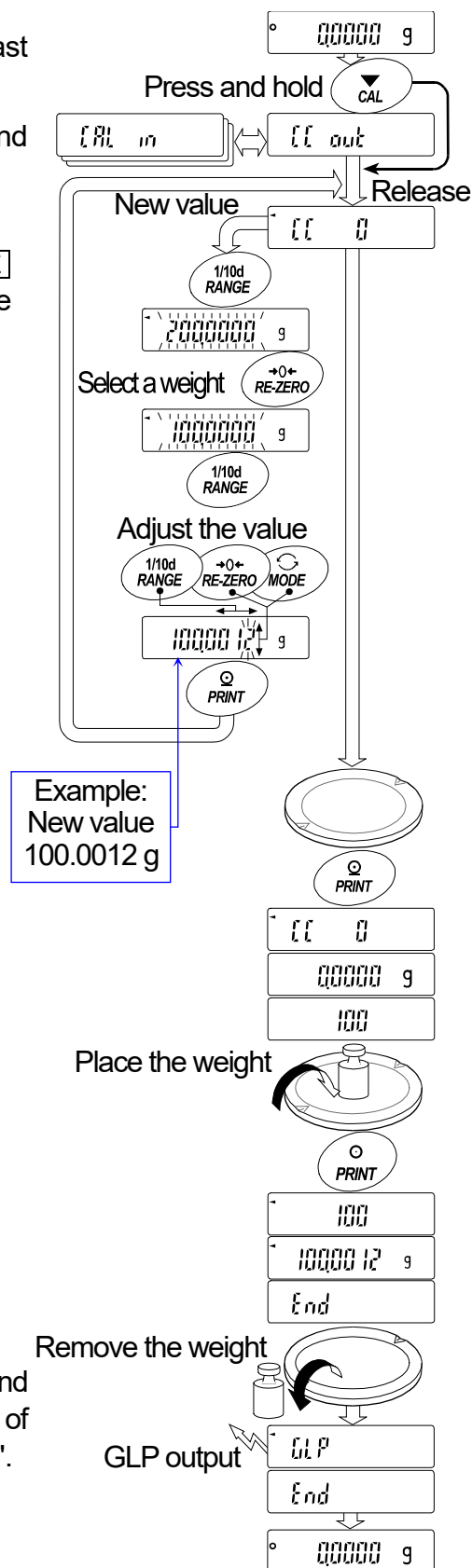

## 8.6. Correcting the Internal Mass Value

- The balance can correct the internal mass value within the range shown below. This function corrects the internal mass value to conform to an external weight. The corrected mass value is maintained in non-volatile memory even if the AC adapter is removed. The internal mass value is corrected as follows:

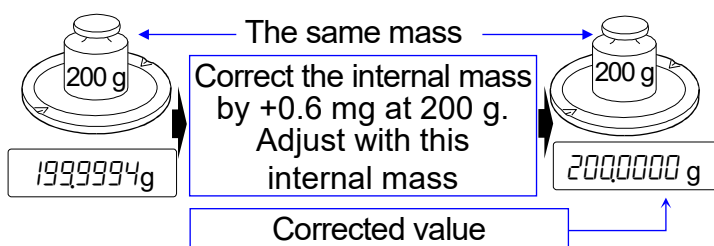

| Model                                 | Correction reference value | Correction range |
|---------------------------------------|----------------------------|------------------|
| <b>BM-5, BM-5D</b>                    | 5.00000 g                  | ±1.5 mg          |
| <b>BM-20, BM-22</b>                   | 20.00000 g                 | ±1.5 mg          |
| <b>BM-252, BM-200, BM-300, BM-500</b> | 200.0000 g                 | ±1.5 mg          |

**Example:** With the **BM-300**, 200.0000 g is corrected by +0.6 mg (200.0000 g). When performing a +0.6 mg correction with a 100 g external weight, the correction value will be +1.2 mg in accordance with the reference value of 200 g.

- Perform sensitivity adjustment using the internal mass. Weigh an external weight and get the correction value.
- Press the **ON:OFF** key to turn the display off.
- While pressing and holding the **PRINT** key and the **RANGE** key, press the **ON:OFF** key. **P5** is displayed.
- Press the **PRINT** key to display the function switches. Set the function table switch and internal mass correction switch to "1" as shown to the right using the following keys.


**RANGE** key ..... To select the blinking digit.  
**RE-ZERO** key ..... To change the value of the blinking digit.  
**PRINT** key ..... To store the new value and return to the weighing mode.  
**CAL** key ..... To cancel the current operation.

- Press and hold the **RANGE** key (for approx. 2 seconds) to enter the function table and release the key when **bR5Fnc** is displayed.
- Press the **RANGE** key several times to display **[5] in**.
- Press the **PRINT** key.

Correct the internal mass value using the following keys.

**RE-ZERO**(+)**key** .... To increase the value.  
**MODE**(-)**key** ..... To decrease the value.  
**PRINT** key ..... To store the new value and display the next menu item of the function table.  
**CAL** key ..... To cancel the current operation and display the next menu item of the function table.

- Press the **CAL** key to return to the weighing mode.
- Press the **CAL** key to adjust the balance using the internal mass.
- Check that the correction has been performed properly with the external weight. If the value is incorrect, repeat the correction.

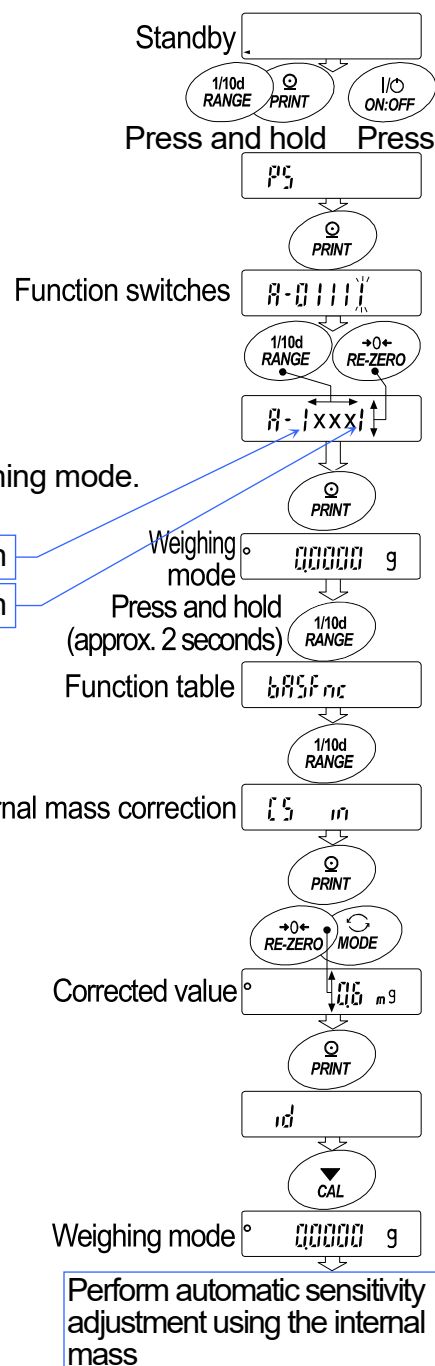

## 9. Function Switch and Initialization

### 9.1. Permit or Inhibit

- The balance stores parameters that must not be changed unintentionally.  
Example: Sensitivity adjustment history for accurate weighing, Data for adapting to the operating environment, Control data for the RS-232C interface.
- The balance is equipped with five switches for the purpose of protecting parameters as shown above. Each switch can select either "permit" or "inhibit". "Inhibit" protects parameters against unintentional operations.  
In addition to the above, when using balances with software version 1.660 or later, the password function can be used to restrict sensitivity adjustment functions.

- 1 Press the **ON:OFF** key to turn the display off.
- 2 While pressing and holding the **PRINT** key and the **RANGE** key, press the **ON:OFF** key to display **P5**.
- 3 Press the **PRINT** key. Then the balance displays the function switches.
- 4 Specify the switches using the following keys.

**RANGE** key..... To select the blinking digit.

**RE-ZERO** key..... To change the parameter of the selected switch.

0 To inhibit changes. (Cannot be used.)

1 To permit changes. (Can be used.)

**PRINT** key ..... To store the new parameter and return to the weighing mode.

**CAL** key ..... To cancel the operation ( **Err** is displayed ). To return to the weighing mode, press the **CAL** key once again.

(The display shown left indicates the factory settings)

**A-01111**

#### Function table

□ To inhibit changes to the function table.

/ To permit changes to the function table.

#### Sensitivity adjustment using the internal mass

□ Software ver. 1.00 to 1.64 To inhibit sensitivity adjustment using the internal mass.

Software ver. 1.660 or later **PASS 0** To inhibit sensitivity adjustment using the internal mass.

**PASS 1,2** To inhibit sensitivity adjustment using the internal mass

See Note 1. when logged in as a user (USER) or a guest (GUEST).

/ To permit sensitivity adjustment using the internal mass.

#### Sensitivity adjustment using an external weight

□ Software ver. 1.00 to 1.64 To inhibit sensitivity adjustment using an external weight.

Software ver. 1.660 or later **PASS 0** To inhibit sensitivity adjustment using an external weight.

**PASS 1,2** To inhibit sensitivity adjustment using an external weight

See Note 1. when logged in as a user (USER) or a guest (GUEST).

/ To permit sensitivity adjustment using an external weight.

#### Automatic sensitivity adjustment (for variation of ambient temperature)

□ To inhibit automatic sensitivity adjustment.

/ To permit automatic sensitivity adjustment.

#### Internal mass correction

□ To inhibit correction.

/ To permit correction.


**Note 1:** The operation is permitted when logged in as an administrator (Admin).

## 9.2. Initializing the Balance

- This function restores the following parameters to the factory settings.
  - Sensitivity adjustment history
  - Function table (excluding the time and date)
  - Sample unit mass value (counting mode),  
100% reference mass value (percent mode)
  - The data that is stored in the balance using the data memory function
  - External weight value
  - Function switch settings ("9.1. Permit or Inhibit")
  - Liquid density and temperature in the density mode

**Note** Be sure to adjust the balance after initialization.

- 1 Press the **ON:OFF** key to turn the display off.
- 2 While pressing and holding the **PRINT** key and the **RANGE** key, press the **ON:OFF** key to display **P5**.
- 3 Press the **RANGE** key to display **CLr**.
- 4 Press the **PRINT** key to display **CLr 00**.  
To cancel the operation, press the **CAL** key.
- 5 Press the **RE-ZERO** key to display **CLr 00**.
- 6 Press the **PRINT** key to initialize the balance.  
The balance will automatically return to the weighing mode.

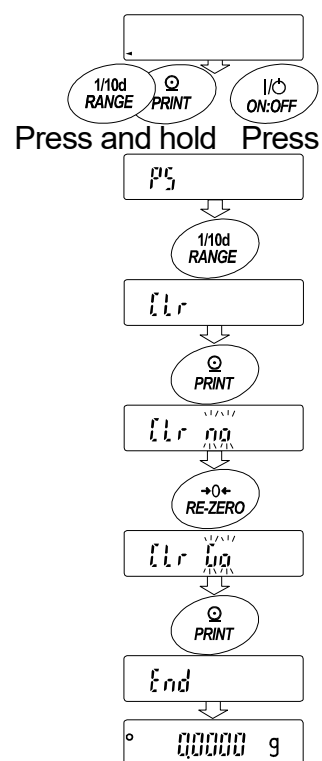






## 10. Function Table

The function table reads or rewrites the parameters that are stored in the balance.  
 These parameters are maintained in non-volatile memory, even if the AC adapter is removed.  
 The function table menu consists of two layers. The first layer is the "Class" and the second layer is the "Item".

### 10.1. Setting the Function Table

#### Display symbol and keys


|   |   |
|---|---|
|    | The symbol "o" indicates that the parameter displayed is in effect.   |
|    | When pressed and held (for approx. 2 seconds) in the weighing mode, enters the function table mode.<br>Selects the class or item in the function table mode.                    |
|    | Changes the parameter.  |
|  | When a class is displayed, enters an item in the class.<br>When an item is displayed, stores the new parameter and displays the next class.                                     |
|  | When an item is displayed, cancels the new parameter and displays the next class.<br>When a class is displayed, exits the function table mode and returns to the weighing mode. |

#### Setting procedure

- 1 In the weighing mode, press and hold the **RANGE** key (for approx. 2 seconds) until **bASFnC** is displayed, and then release the key.
- 2 Press the **RANGE** key to select a class.
- 3 Press the **PRINT** key to enter the class.
- 4 Press the **RANGE** key to select an item.
- 5 Press the **RE-ZERO** key to select a parameter for the selected item.
- 6 Repeat Steps 4 and 5 to select other items of the class.
- 7 To store the parameters of the selected class, press the **PRINT** key.  
 Then the next class is displayed.  
 To cancel the current operation, press the **CAL** key. Then the next class is displayed.
- 8 To specify parameters for another class, proceed to Step 2.  
 To finish the setting, press the **CAL** key to return to the weighing mode.

## Setting example

This example sets the "Data memory (*data*)" parameter to "1" (To store the weighing data) and the "Interval time (*int*)" parameter to "5" (1 minute).

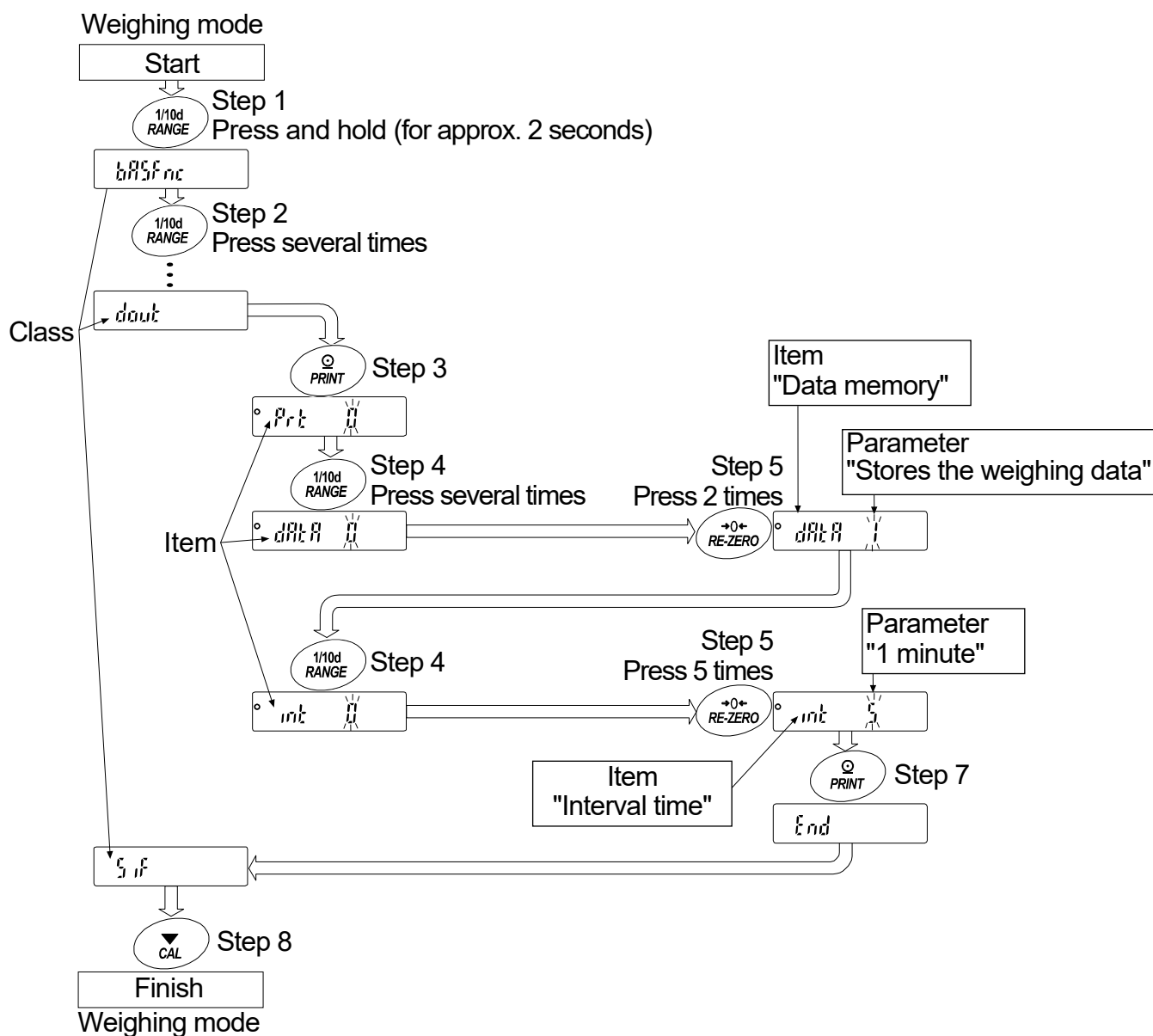



## 10.2. Details of the Function Table

| Class                            | Item and Parameter                            |   | Description  |  |
|----------------------------------|---|---|--|--|
| bRSFnc<br>Environment<br>Display | Cond<br>Condition                             | 0   | Fast response, sensitive value                                       | Can be changed by response adjustment.   |
|                                  |   | 1   |  |  |
|                                  |   | 2   | Slow response, stable value  |  |
|                                  | St-b<br>Stability band width                  | 0   | Stable range is ±1 digit   | #1<br>The stabilization indicator illuminates when the display fluctuation is within the range.            |
|                                  |   | 1   | ±2 digit   |  |
|                                  |   | 2   | Stable range is ±3 digits  |  |
|                                  | PLoc<br>Display lock function                 | OFF                                       | OFF  | Display lock function to measure the volume dispensed from the pipette.                                    |
|                                  |   | ON  | ON   |  |
|                                  | trc<br>Zero tracking                          | 0   | OFF  | Keeps zero display by tracking zero drift.   |
|                                  |   | 1   | Normal   |  |
|                                  |   | 2   | Strong   |  |
|                                  | SPd<br>Display refresh rate                   | 0   | 5 times/second   | Period to refresh the display  |
| 1                                |   | 10 times/second                           |  |  |
| Pnt<br>Decimal point             | 0   | Point (.)                                 | Decimal point format   |  |
|                                  | 1   | Comma (,)                                 |  |  |
| P-on<br>Auto display-ON          | OFF   | OFF                                       | Turns on the weighing mode display when the AC adapter is connected. |  |
|                                  | ON  | ON  |  |  |
| bEEP<br>Beep                     | OFF   | OFF                                       | Sounds a beep when a key is touched.                                 |  |
|                                  | ON  | ON  |  |  |
| CL Add<br>Clock                  | Refer to "10.7. Clock and Calendar Function " |   | The time and date are added to the output data.                      |  |
| dout<br>Data output              | Prt<br>Data output mode                       | 0   | Key mode   | Accepts the <span>PRINT</span> key only when the display is stable.  |
|                                  |   | 1   | Auto print mode A<br>Reference = zero                                | Outputs data when the display is stable and conditions of $RP-P$ , $RP-b$ and the reference value are met. |
|                                  |   | 2   | Auto print mode B<br>Reference = last stable value                   |  |
|                                  |   | 3   | Stream mode / Interval mode  | With $dAtR$ 0, outputs data continuously; with $dAtR$ 1, uses interval memory.                             |
|                                  | RP-P<br>Auto print polarity                   | 0   | Plus only  | Displayed value > Reference  |
|                                  |   | 1   | Minus only   | Displayed value < Reference  |
|                                  |   | 2   | Both   | Regardless of displayed value  |
|                                  | RP-b<br>Auto print difference                 | 0   | 10 digits  | #2<br>Difference between reference value and displayed value   |
|                                  |   | 1   | 100 digits   |  |
|                                  |   | 2   | 1000 digits  |  |
|                                  | dAtR<br>Data memory                           | 0   | Not used   | Related items:<br>$Prt$ , $int$ , $d-no$ , $S-td$ , $info$   |
|                                  |   | 1   | Stores the weighing data   |  |
| 2                                |   | Stores the sensitivity adjustment history |  |  |

▪ : Factory settings.

#1 The unit of readability is digit.

Example: If 1 mg display is selected using the RANGE key for the **BM-300**, 1 mg is one digit.

#2 Usable readability of the balance is one digit.

Example: In gram display, one digit is 0.01 mg for the **BM-252** and 0.1 mg for the **BM-300**.

| Class                   | Item and Parameter          |       | Description                             |   |
|-------------------------|-----------------------------|-------|---|---|
| dout<br>Data output     | int<br>Interval time        | ▪ 0   | Every measurement                       | Interval time in the interval memory mode when using <i>Prt 3, dAtA 1</i>   |
|                         |                             | 1     | 2 seconds                               |   |
|                         |                             | 2     | 5 seconds                               |   |
|                         |                             | 3     | 10 seconds                              |   |
|                         |                             | 4     | 30 seconds                              |   |
|                         |                             | 5     | 1 minute                                |   |
|                         |                             | 6     | 2 minutes                               |   |
|                         |                             | 7     | 5 minutes                               |   |
|                         |                             | 8     | 10 minutes                              |   |
|                         | d-no<br>Data number output  | ▪ OFF | No output                               | Refer to "12. Data Memory"  |
|                         |                             | on    | Output                                  |   |
|                         | S-t-d<br>Time/Date output   | ▪ 0   | No output                               | Selects whether or not the time or date is added to the weighting data. Refer to "10.7. Clock and Calendar Function" for details. |
|                         |                             | 1     | Time only                               |   |
|                         |                             | 2     | Date only                               |   |
|                         |                             | 3     | Time and date                           |   |
|                         | S-id<br>ID number output    | ▪ OFF | No output                               | Selects whether or not the ID number is output.   |
|                         |                             | on    | Output                                  |   |
|                         | PUSE<br>Data output pause   | ▪ OFF | No pause                                | Selects the data output interval.   |
|                         |                             | on    | Pause (1.6 seconds)                     |   |
|                         | At-F<br>Auto feed           | ▪ OFF | Not used                                | Selects whether or not automatic feed is performed.   |
|                         |                             | on    | Used                                    |   |
|                         | info<br>GLP output          | ▪ 0   | No output                               | Refer to "11. ID Number and GLP Report" for details.<br>#3 Available for balances with software version 1.660 or later.           |
|                         |                             | 1     | AD-8121 format                          |   |
|                         |                             | 2     | General (balance clock data)            |   |
|                         |                             | 3     | General (external device clock data) #3 |   |
|                         | Pr-d<br>Zero after output   | ▪ OFF | Not used                                | Adjusts zero automatically after data is output   |
|                         |                             | on    | Used                                    |   |
|                         | S-t-P<br>Temperature output | ▪ OFF | Not output                              | Adds temperature to data output   |
|                         |                             | on    | Output                                  |   |
| SIF<br>Serial interface | bps<br>Baud rate            | 0     | 600 bps                                 |   |
|                         |                             | 1     | 1200 bps                                |   |
|                         |                             | ▪ 2   | 2400 bps                                |   |
|                         |                             | 3     | 4800 bps                                |   |
|                         |                             | 4     | 9600 bps                                |   |
|                         |                             | 5     | 19200 bps                               |   |
|                         | bPr<br>Data bit, parity bit | ▪ 0   | 7 bits, even                            |   |
|                         |                             | 1     | 7 bits, odd                             |   |
|                         |                             | 2     | 8 bits, none                            |   |
|                         | CrLF<br>Terminator          | ▪ 0   | CR LF                                   | CR: ASCII 0Dh<br>LF: ASCII 0Ah  |
|                         |                             | 1     | CR                                      |   |
|                         | tYPE<br>Data format         | ▪ 0   | A&D standard format                     | Refer to "10.5. Description of the Item "Data Format".  |
|                         |                             | 1     | DP format                               |   |
|                         |                             | 2     | KF format                               |   |
|                         |                             | 3     | MT format                               |   |
|                         |                             | 4     | NU format                               |   |
|                         |                             | 5     | CSV format                              |   |

▪ : Factory settings.

**Caution** The balance may not transmit the data completely at the specified refresh rate, depending on the baud rate or data added to the weighing data such as time, date and ID number.

| Class                    | Item and Parameter     |     | Description   |  |  |
|--------------------------|------------------------|-----|---|--|--|
| Serial interface         | Timeout                | OFF | No limit  | Selects the wait time to receive a command.      |  |
|                          |                        | ON  | 1 second  |  |  |
|                          | AK, Error code         | OFF | No output   | AK: ASCII 06h                                    |  |
|                          |                        | ON  | Output  |  |  |
| Density function         | Liquid density input   | 0   | Water temperature   | Available only when density mode is selected     |  |
|                          |                        | 1   | Liquid density  |  |  |
| Unit                     |                        |     | Refer to "5. Weighing Units".   |  |  |
| Internal mass correction |                        |     | Displayed only when the internal mass value correction switch is set to 1. Refer to "8. Sensitivity Adjustment / Calibration Test". |  |  |
| ID number setting        |                        |     | Refer to "11. ID Number and GLP Report".  |  |  |
| Ionizer function         | Neutralization time    | 0   | 3 seconds   | Refer to "6.5. Ionizer (Neutralization Device)". |  |
|                          |                        | 1   | 10 seconds  |  |  |
|                          |                        | 2   | Manual operation (max. 10 min.)   |  |  |
|                          | AD8922A control        | OFF | The [ION] key is not used.  |  |  |
|                          |                        | ON  | Assigns the [ION] key to the [MODE] key on AD-8922A.  |  |  |
|                          | External control input | 0   | Assigns the [PRINT] key and [RE-ZERO] key.  |  |  |
|                          |                        | 1   | Assigns the [PRINT] key and [ION] key.  |  |  |
|                          |                        | 2   | Assigns the [ION] key and [RE-ZERO] key.  |  |  |
| Door sensor function     | Door sensor            | OFF | Not used  | Refer to "6.1. Basic Operation"                  |  |
|                          |                        | ON  | Used  |  |  |
| Password function        | #4 Password function   | 0   | Not used  | Refer to "15. Password Function".                |  |
|                          |                        | 1   | Used (Limits weighing operations)   |  |  |
|                          |                        | 2   | Used (Allows basic weighing)  |  |  |
| Password                 | Admin                  |     | Password input by the administrator   |  |  |
|                          | USER 01                |     | Password input by user 1  |  |  |
|                          | USER 10                |     | Password input by user 10 #5  |  |  |

0 : Factory settings.

#4 Only "Not used (0)" and "Used (1)" are available for balances with software version 1.50 to 1.64. All the parameters are available for balances with software version 1.660 or later.

#5 Available for balances with software version 1.660 or later.

For balances with software version 1.50 to 1.64, only the password input by the administrator is available.

## 10.3. Description of the Class "Environment, Display"

### Condition (Cond)

Cond 0 This parameter is for sensitive response to the fluctuation of a mass value. Used for powder target weighing, weighing a very light sample or when quick response weighing is required. After setting, the balance displays [FAST].

Cond 2 This parameter is for stable weighing with slow response. Used to prevent a weight value from drifting due to vibration or drafts. After setting, the balance displays [SLOW].

**Note** In automatic response adjustment, this parameter is selected automatically.

## Stability band width ( $Sb$ )

This item controls the width to regard a mass value as a stable value. When the fluctuation per second is less than the parameter, the balance displays the stabilization indicator and outputs or stores the data. The parameter influences the "Auto print mode".

$Sb$  0 This parameter is used for sensitive response of the stabilization indicator. Used for exact weighing.

$Sb$  2 This parameter ignores slight fluctuations of a mass value. Used to prevent a mass value from drifting due to vibration or drafts.


## Display lock function ( $PLoc$ )

□ This function is used to lock the display when measuring the volume dispensed from the pipette. With "Display lock function is used (  $PLoc$  on )" selected in the function table, when the liquid is dispensed into the weighing bottle from the pipette and a stable weighing value is obtained, the balance calculates an average value and displays the result (locked).

□ When using this function for the measurement of a small volume, the influence of liquid evaporation can be reduced.

**Caution** Even if the weighing bottle (evaporator trap) to reduce the water evaporation is used, at least 0.05 mg (0.05  $\mu\text{l}$ ) evaporation will occur.

□ When canceling this function, press the RE-ZERO key.

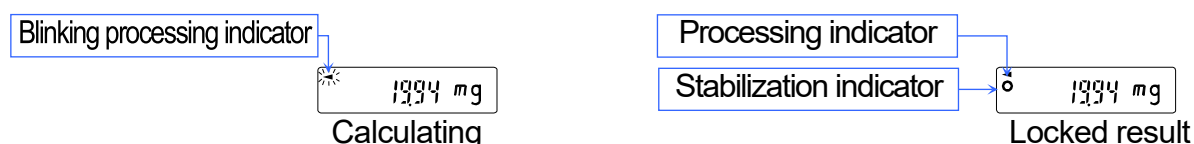

## Zero tracking ( $trc$ )

This function tracks zero point drift caused by changes in the environment and stabilizes the zero point. When the weighing data is only a few digits, turn the function off for accurate weighing.

$trc$  0 The tracking function is not used. Used for weighing a very light sample.

$trc$  1 The normal tracking function is used.

$trc$  2 The strong tracking function is used.

$trc$  3 The very strong tracking function is used. Used for stable zero display.

## Display refresh rate ( $SPd$ )

The periodic time to refresh the display. This parameter influences "Baud rate", "Data output pause" and the data output rate of "Stream mode".

## Decimal point ( $Pnt$ )

The decimal point format can be selected.

## Auto display-ON ( $P-on$ )

When the AC adapter is plugged in, the display is automatically turned on without pressing the ON:OFF key, to display the weighing mode. Used when the balance is built into an automated system. One hour warm up is necessary for accurate weighing.

## 10.4. Description of the Item "Data Output Mode"

The parameter setting of "Data output mode (*Prt*)" applies to the performance when the "Data memory (*dRtR*)" parameter is set to "1" (to store the weighing data) and when the data is transmitted using the RS-232C interface.

### Key mode

When the PRINT key is pressed with the stabilization indicator turned on, the balance outputs or stores the weighing data and the display blinks one time.

Required setting    *dout*    *Prt* 0    Key mode

### Auto print modes A and B

When the displayed value is stable and the conditions of "Auto print polarity", "Auto print difference" and the reference value are met, the balance outputs or stores the weighing data.

When the PRINT key is pressed with the stabilization indicator turned on, the balance outputs or stores the data and the display blinks one time.

#### Auto print mode A

Example                      For weighing each time a sample is placed and removed, with "*Rr-d*" set to "1" (to adjust zero after the data is output).

Required setting    *dout*    *Prt* 1    Auto print mode A (reference = zero)  
                          *dout*    *RP-P*    Auto print polarity  
                          *dout*    *RP-b*    Auto print difference  
                          *dout*    *Rr-d* on    Zero after output

#### Auto print mode B

Example                      For weighing while a sample is added.

Required setting    *dout*    *Prt* 2    Auto print mode B (reference = last stable value)  
                          *dout*    *RP-P*    Auto print polarity  
                          *dout*    *RP-b*    Auto print difference

### Stream mode

The balance outputs the weighing data continuously regardless of the display condition. When the display refresh rate is set to 5 times / second (*SPd* 0), the data output rate is also set to the same 5 times / second. The display does not blink in this mode. The interval memory mode is used when the "Data memory (*dRtR*)" parameter is set to "1" (to store the weighing data).

Example                      For monitoring data on a computer.

Required setting    *dout*    *Prt* 3    Stream mode  
                          *dout*    *dRtR* 0    Data memory function is not used  
                          *bRSFnc* *SPd*    Display refresh rate  
                          *SIF*    *bPS*    Baud rate

**Caution**    The balance may not transmit the data completely at the specified refresh rate, depending on the baud rate or data added to the weighing data such as time, date and ID number.

### Interval mode

The balance outputs the weighing data periodically regardless of the display condition and the display blinks one time. When the "Data memory (*dRtR*)" parameter of "Data output (*dout*)" is set to "1" (to store the weighing data), the weighing data is periodically stored in memory.

Pressing the PRINT key starts output and pressing the PRINT key again during output stops the output.

Example                      For periodical weighing without a personal computer command and to output all of the data, to a computer, at one time.

Required setting    *dout*    *Prt* 3    Interval mode  
                          *dout*    *dRtR* 1    Data memory function is used  
                          *dout*    *int*    Interval time


## 10.5. Description of the Item "Data Format"

### A&D standard format

5 IF TYPE 0

This format is used when the peripheral equipment can receive the A&D format. With an AD-8121B printer, use MODE 1 or 2. With an AD-8127 printer or a thermal printer, AD-8129TH, use the external key printing mode, manual printing mode, auto printing mode or interval printing mode.

- This format consists of fifteen or sixteen characters excluding the terminator.
- A header of two characters indicates the balance condition.
- The polarity sign is placed before the data with the leading zeros. If the data is zero, the plus sign is applied.
- The unit, consisting of three characters, follows the data.

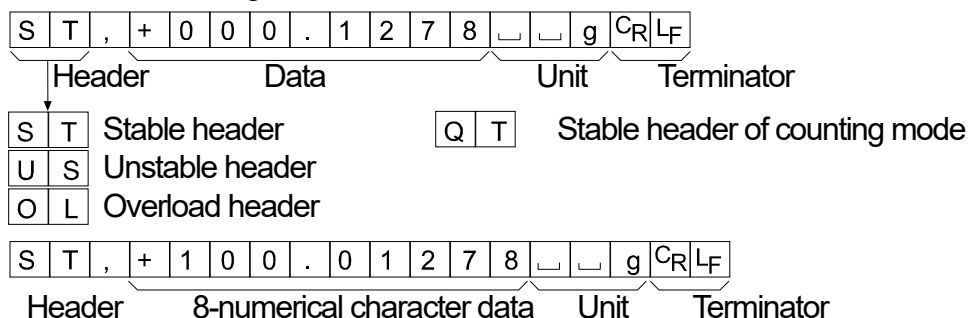


### DP (Dump print) format

5 IF TYPE 1

This format is used when the peripheral equipment cannot receive the A&D format.

With an AD-8121B printer, use MODE 3. With an AD-8127 printer or a thermal printer, AD-8129TH, use the dump printing mode.

- This format consists of sixteen characters excluding the terminator.
- A header of two characters indicates the balance condition. No overload header is used.
- The polarity sign is placed before the data, with spaces in place of leading zeros, if the data is not zero or overloaded.
- The unit, consisting of three characters, follows the data.

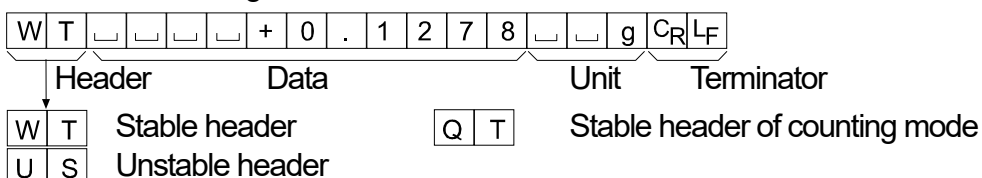


### KF format

5 IF TYPE 2

This is the Karl-Fischer moisture meter format and is used when the peripheral equipment can only communicate using this format.

- This format consists of fourteen characters excluding the terminator.
- This format has no header characters.
- The polarity sign is placed before the data, with spaces in place of leading zeros, if the data is not zero or overloaded.
- This format outputs the unit only for a stable value.

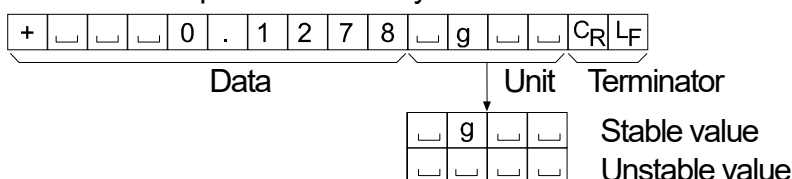


## MT format

5 IF TYPE 3

This format is used when the peripheral equipment of other manufacturer is connected. Please note that the connection is not guaranteed.

- A header of two characters indicates the balance condition.
- The polarity sign is used only for negative data.
- The weighing data uses spaces in place of the leading zeros.
- The character length of this format changes dependent upon the unit

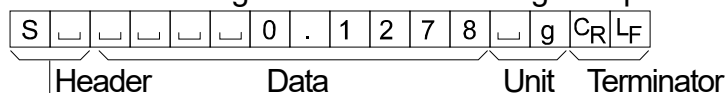


|   |   |   |
|---|---|---|
| S |   | Stable header   |
| S | D | Unstable header   |
| S | I | Overload header   |
|   |   | Stable header (When data is output using the <span style="border: 1px solid black; padding: 0 2px;">PRINT</span> key) |

## NU (numerical) format

5 IF TYPE 4

This format outputs only numerical data.

- This format consists of ten characters excluding the terminator.
- The polarity sign is placed before the data with the leading zeros. If the data is zero, the plus sign is used.

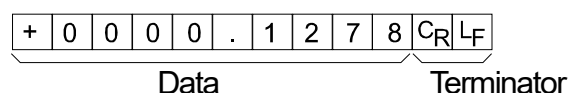

## CSV format

5 IF TYPE 5

- This format separates the data of A&D standard format and the unit by a comma ( , ).
- This format outputs the unit even when the data is overloaded.
- When a comma ( , ) is selected for the decimal point, the separators are set to semicolon ( ; ).
- When the ID number, data number, time and date are added in "Data output (*dout*)" of the function table, outputs ID number, data number, date, time and weighing data in this order and separates each item by a comma and treats all the items as one group of data.

LAB-0123, No,012, 2010/11/01, 12:34:56, ST,+0000.1278, g<CR><LF>

ID number
Data number
Date
Time
Weighing data

|   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |                |                |   |                |                |
|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|----------------|----------------|---|----------------|----------------|
| S | T | , | + | 0 | 0 | 0 | . | 1 | 2 | 7 | 8 | , |   |   | g | C <sub>R</sub> | L <sub>F</sub> |   |                |                |
| O | L | , | + | 9 | 9 | 9 | 9 | 9 | 9 | 9 | E | + | 1 | 9 | , |                |                | g | C <sub>R</sub> | L <sub>F</sub> |

**Note** When the balance is connected to a personal computer via the USB interface, the ID number, data number, time and date cannot be added to the weighing data.

## ID number

*dout 5- id on*

The number to identify a specific balance.

- This format consists of eight characters excluding the terminator.

|   |   |   |   |   |   |   |   |                |                |
|---|---|---|---|---|---|---|---|----------------|----------------|
| L | A | B | - | 0 | 1 | 2 | 3 | C <sub>R</sub> | L <sub>F</sub> |
|---|---|---|---|---|---|---|---|----------------|----------------|

## Data number

*dout d-no on*

This format outputs the data number just before the data is transmitted using the RS-232C interface.

- This format consists of six characters excluding the terminator.
- When CSV format (*5 IF TYPE 5*) is selected, the period ( . ) is replaced with a comma ( , ).

|   |   |   |   |   |   |                |                |
|---|---|---|---|---|---|----------------|----------------|
| N | O | . | 0 | 0 | 1 | C <sub>R</sub> | L <sub>F</sub> |
|---|---|---|---|---|---|----------------|----------------|

Data number

Terminator

## Date

*dout S-td 2 or 3*

- The date output order can be changed in "Time/Date output (*S-td*)" and "Clock (*CL Adj*)".

This format outputs the year in a four-digit format.

|   |   |   |   |   |   |   |   |   |   |                |                |
|---|---|---|---|---|---|---|---|---|---|----------------|----------------|
| 2 | 0 | 1 | 0 | / | 1 | 1 | / | 0 | 1 | C <sub>R</sub> | L <sub>F</sub> |
|---|---|---|---|---|---|---|---|---|---|----------------|----------------|

## Time

*dout S-td 1 or 3*

- This format outputs the time in a 24-hour format.

|   |   |   |   |   |   |   |   |                |                |
|---|---|---|---|---|---|---|---|----------------|----------------|
| 1 | 2 | : | 3 | 4 | : | 5 | 6 | C <sub>R</sub> | L <sub>F</sub> |
|---|---|---|---|---|---|---|---|----------------|----------------|

## Temperature

*dout S-tP on*

- This format outputs the temperature. The unit is "C".

**Note** Humidity and barometric pressure are not output.

|   |   |   |   |   |   |   |   |   |                |                |
|---|---|---|---|---|---|---|---|---|----------------|----------------|
| + | 0 | 2 | 3 | . | 4 | □ | □ | C | C <sub>R</sub> | L <sub>F</sub> |
|---|---|---|---|---|---|---|---|---|----------------|----------------|



## 10.6. Data Format Examples

### Stable

° 0.1278 g

|   |   |   |   |   |   |   |   |   |   |   |                |                |   |                |                |                |                |
|---|---|---|---|---|---|---|---|---|---|---|----------------|----------------|---|----------------|----------------|----------------|----------------|
| S | T | , | + | 0 | 0 | 0 | . | 1 | 2 | 7 | 8              | □              | □ | g              | C <sub>R</sub> | L <sub>F</sub> |                |
| S | T | , | + | 1 | 0 | 0 | . | 0 | 1 | 2 | 7              | 8              | □ | □              | g              | C <sub>R</sub> | L <sub>F</sub> |
| W | T | □ | □ | □ | □ | + | 0 | . | 1 | 2 | 7              | 8              | □ | □              | g              | C <sub>R</sub> | L <sub>F</sub> |
| + | □ | □ | □ | 0 | . | 1 | 2 | 7 | 8 | □ | g              | □              | □ | C <sub>R</sub> | L <sub>F</sub> |                |                |
| S | □ | □ | □ | □ | □ | 0 | . | 1 | 2 | 7 | 8              | □              | g | C <sub>R</sub> | L <sub>F</sub> |                |                |
| + | 0 | 0 | 0 | 0 | 0 | . | 1 | 2 | 7 | 8 | C <sub>R</sub> | L <sub>F</sub> |   |                |                |                |                |

### Unstable

- 18.3690 g

|   |   |   |   |   |   |   |   |   |   |                |                |   |   |                |                |                |                |
|---|---|---|---|---|---|---|---|---|---|----------------|----------------|---|---|----------------|----------------|----------------|----------------|
| U | S | , | - | 0 | 1 | 8 | . | 3 | 6 | 9              | 0              | □ | □ | g              | C <sub>R</sub> | L <sub>F</sub> |                |
| U | S | , | - | 1 | 0 | 1 | . | 8 | 3 | 6              | 9              | 0 | □ | □              | g              | C <sub>R</sub> | L <sub>F</sub> |
| U | S | □ | □ | □ | - | 1 | 8 | . | 3 | 6              | 9              | 0 | □ | □              | g              | C <sub>R</sub> | L <sub>F</sub> |
| - | □ | □ | 1 | 8 | . | 3 | 6 | 9 | 0 | □              | □              | □ | □ | C <sub>R</sub> | L <sub>F</sub> |                |                |
| S | D | □ | □ | - | 1 | 8 | . | 3 | 6 | 9              | 0              | □ | g | C <sub>R</sub> | L <sub>F</sub> |                |                |
| - | 0 | 0 | 1 | 8 | . | 3 | 6 | 9 | 0 | C <sub>R</sub> | L <sub>F</sub> |   |   |                |                |                |                |

### Overload

Positive error

g

|   |   |   |                |                |   |   |   |   |   |   |                |                |   |   |                |                               |
|---|---|---|----------------|----------------|---|---|---|---|---|---|----------------|----------------|---|---|----------------|-------------------------------|
| O | L | , | +              | 9              | 9 | 9 | 9 | 9 | 9 | 9 | E              | +              | 1 | 9 | C <sub>R</sub> | L <sub>F</sub>                |
| ␣ | ␣ | ␣ | ␣              | ␣              | ␣ | ␣ | ␣ | ␣ | E | ␣ | ␣              | ␣              | ␣ | ␣ | ␣              | C <sub>R</sub> L <sub>F</sub> |
| ␣ | ␣ | ␣ | ␣              | ␣              | ␣ | H | ␣ | ␣ | ␣ | ␣ | ␣              | ␣              | ␣ | ␣ | C <sub>R</sub> | L <sub>F</sub>                |
| S | I | + | C <sub>R</sub> | L <sub>F</sub> |   |   |   |   |   |   |                |                |   |   |                |                               |
| + | 9 | 9 | 9              | 9              | 9 | 9 | 9 | 9 | 9 | 9 | C <sub>R</sub> | L <sub>F</sub> |   |   |                |                               |

### Overload

Negative error

-g

|   |   |   |                |                |   |   |   |   |   |   |                |                |   |   |                |                               |
|---|---|---|----------------|----------------|---|---|---|---|---|---|----------------|----------------|---|---|----------------|-------------------------------|
| O | L | , | -              | 9              | 9 | 9 | 9 | 9 | 9 | 9 | E              | +              | 1 | 9 | C <sub>R</sub> | L <sub>F</sub>                |
|   |   |   |                |                |   |   |   | - | E |   |                |                |   |   |                | C <sub>R</sub> L <sub>F</sub> |
|   |   |   |                |                |   |   | L |   |   |   |                |                |   |   | C <sub>R</sub> | L <sub>F</sub>                |
| S | I | - | C <sub>R</sub> | L <sub>F</sub> |   |   |   |   |   |   |                |                |   |   |                |                               |
| - | 9 | 9 | 9              | 9              | 9 | 9 | 9 | 9 | 9 | 9 | C <sub>R</sub> | L <sub>F</sub> |   |   |                |                               |

Space, ASCII 20h  
 C<sub>R</sub> Carriage Return, ASCII 0Dh  
 L<sub>F</sub> Line Feed, ASCII 0Ah

## Units

|                              |     | A&D   | D.P.  | KF   | MT   |
|------------------------------|-----|---|---|--|--|
| g                            | g   | <input type="text"/> <input type="text"/> g     | <input type="text"/> <input type="text"/> g     | <input type="text"/> g <input type="text"/> <input type="text"/>     | <input type="text"/> <input type="text"/> g                          |
| mg                           | mg  | <input type="text"/> <input type="text"/> m g   | <input type="text"/> <input type="text"/> m g   | <input type="text"/> <input type="text"/> m g <input type="text"/>   | <input type="text"/> <input type="text"/> m g                        |
| Counting mode                | PC  | <input type="text"/> <input type="text"/> P C   | <input type="text"/> <input type="text"/> P C   | <input type="text"/> <input type="text"/> p c s                      | <input type="text"/> <input type="text"/> P C S                      |
| Percent mode                 | %   | <input type="text"/> <input type="text"/> %     | <input type="text"/> <input type="text"/> %     | <input type="text"/> % <input type="text"/> <input type="text"/>     | <input type="text"/> %   |
| Ounce (Avoir)                | oz  | <input type="text"/> <input type="text"/> o z   | <input type="text"/> <input type="text"/> o z   | <input type="text"/> <input type="text"/> o z <input type="text"/>   | <input type="text"/> <input type="text"/> o z                        |
| Troy Ounce                   | ozt | <input type="text"/> o z <input type="text"/> t | <input type="text"/> o z <input type="text"/> t | <input type="text"/> <input type="text"/> o z <input type="text"/> t | <input type="text"/> <input type="text"/> o z <input type="text"/> t |
| Metric Carat                 | ct  | <input type="text"/> <input type="text"/> c t   | <input type="text"/> <input type="text"/> c t   | <input type="text"/> <input type="text"/> c t <input type="text"/>   | <input type="text"/> <input type="text"/> c t                        |
| Momme                        | mom | <input type="text"/> m o m                      | <input type="text"/> m o m                      | <input type="text"/> <input type="text"/> m o m                      | <input type="text"/> <input type="text"/> m o                        |
| Pennyweight                  | dwt | <input type="text"/> d w t                      | <input type="text"/> d w t                      | <input type="text"/> <input type="text"/> d w t                      | <input type="text"/> <input type="text"/> d w t                      |
| Grain                        | GN  | <input type="text"/> <input type="text"/> G N   | <input type="text"/> <input type="text"/> G N   | <input type="text"/> <input type="text"/> g r <input type="text"/>   | <input type="text"/> <input type="text"/> G N                        |
| Tael (HK general, Singapore) | tL  | <input type="text"/> <input type="text"/> t l   | <input type="text"/> <input type="text"/> t l   | <input type="text"/> <input type="text"/> t l s                      | <input type="text"/> <input type="text"/> t l                        |
| Tael (HK, jewelry)           | tL  | <input type="text"/> <input type="text"/> t l   | <input type="text"/> <input type="text"/> t l   | <input type="text"/> <input type="text"/> t l h                      | <input type="text"/> <input type="text"/> t l                        |
| Tael (Taiwan)                | tL  | <input type="text"/> <input type="text"/> t l   | <input type="text"/> <input type="text"/> t l   | <input type="text"/> <input type="text"/> t l t                      | <input type="text"/> <input type="text"/> t l                        |
| Tael (China)                 | tL  | <input type="text"/> <input type="text"/> t l   | <input type="text"/> <input type="text"/> t l   | <input type="text"/> <input type="text"/> t l c                      | <input type="text"/> <input type="text"/> t l                        |
| Tola (India)                 | t   | <input type="text"/> <input type="text"/> t     | <input type="text"/> <input type="text"/> t     | <input type="text"/> <input type="text"/> t o l                      | <input type="text"/> <input type="text"/> t                          |
| Messghal                     | m   | <input type="text"/> m e s                      | <input type="text"/> m e s                      | <input type="text"/> <input type="text"/> M S <input type="text"/>   | <input type="text"/> <input type="text"/> m                          |
| Density                      | D5  | <input type="text"/> <input type="text"/> D S   | <input type="text"/> <input type="text"/> D S   | <input type="text"/> <input type="text"/> D S <input type="text"/>   | <input type="text"/> <input type="text"/> D S                        |

Space, ASCII 20h

## 10.7. Clock and Calendar Function

The balance is equipped with a clock and calendar function. When "GLP output (Info)" and "Time/Date output (Setd)" are used, the time and date can be added to the output data. The operation is as follows:

### Operation

- 1 In the weighing mode, press and hold the **RANGE** key (for approx. 2 seconds) until **bRSFnc** is displayed, and then release the key.
- 2 Press the **RANGE** key several times to display **CL Adj.**
- 3 Press the **PRINT** key to enter the time and date setting mode.

### Checking the current time

- 4 The current time is displayed with all the digits blinking.
  - To change the time, press the **RE-ZERO** key and proceed to Step 5.
  - When the time is correct, press the **RANGE** key and proceed to Step 6.
  - To exit the mode, press the **CAL** key and proceed to Step 8.

### Storing a new time

- 5 Set the time in a 24-hour format using the following keys.
  - RE-ZERO**(+)key... To increase the value by one.
  - MODE**(-)key... To decrease the value by one.
  - RANGE** key... To select the digit to change the value.
  - PRINT** key... To store the new time and proceed to Step 6.
  - CAL** key... To cancel the operation and proceed to Step 6.

### Checking the current date

- 6 The current date is displayed with all the digits blinking.
  - To change the display order of year (y), month (m) and day (d), press the **MODE** key. The date is output in the order as specified.
  - To change the date, press the **RE-ZERO** key and proceed to Step 7.
  - When the date is correct, press the **CAL** key and proceed to Step 8.
  - To change the time again, press the **RANGE** key and proceed back to Step 4.


### Storing a new date

- 7 Set the date using the following keys.
  - RE-ZERO**(+)key... To increase the value by one.
  - MODE**(-)key... To decrease the value by one.
  - RANGE** key... To select the digits to change the value.
  - PRINT** key... To store the new date and proceed to Step 8.
  - CAL** key... To cancel the operation and proceed to Step 8.

### Quitting the operation

- 8 The balance displays the next menu item of the function table.  
Press the **CAL** key to exit the mode and return to the weighing mode.

**Note** Do not enter invalid values such as non-existing time and date in this function. When the clock backup battery has been depleted, the balance displays **btc PF**. Under this condition, press any key and set the time and date. The dead battery only affects the clock and calendar function. Even so, the function works normally as long as the AC adapter is connected to the balance.

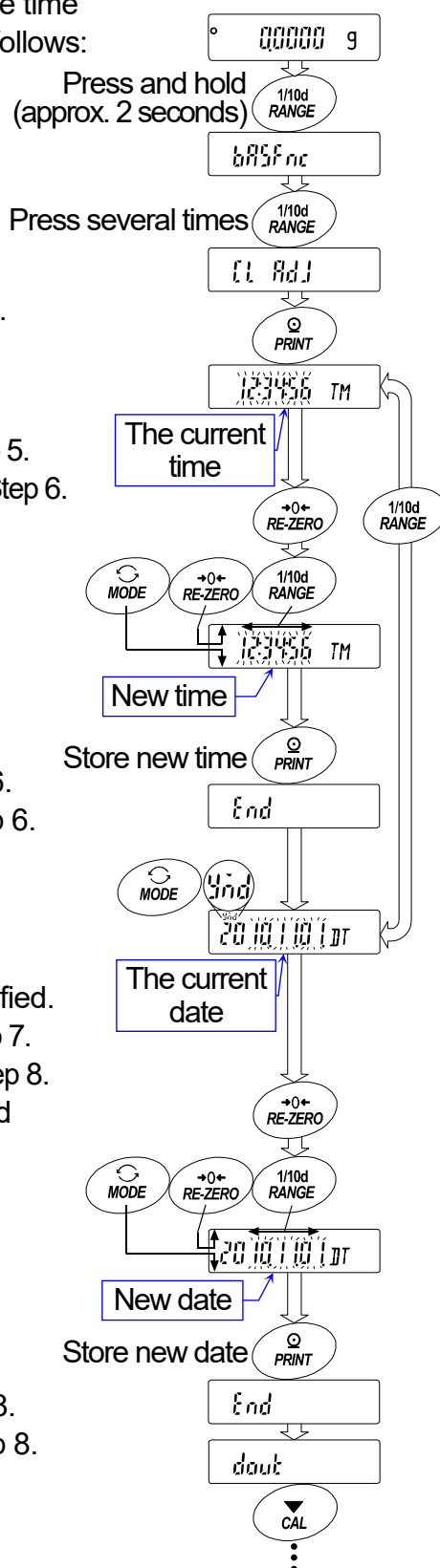

# 11. ID Number and GLP Report

- The ID number is used to identify the balance when Good Laboratory Practice (GLP) or Good Manufacturing Practice (GMP) is used.
- The output format for GLP/GMP compliant report is selected in "GLP output ( *inFo* )" of the function table and can be output to a personal computer or printer using the RS-232C serial interface.
- The GLP/GMP compliant report includes the balance manufacturer, model, serial number, ID number, date, time and space for signature. The sensitivity adjustment report and the calibration test report include the weight used and the result.
- The balance can output the following reports for GLP / GMP using the RS-232C serial interface.
  - "Sensitivity adjustment report" of the automatic sensitivity adjustment and the sensitivity adjustment using the internal mass.
  - "Sensitivity adjustment report" of the sensitivity adjustment using an external weight.
  - "Calibration test report" of the calibration test using the internal mass.
  - "Calibration test report" of the calibration test using an external weight.
  - "Title block" and "End block" for the weighing data.
- Sensitivity adjustment report and calibration test report can be stored in memory and several reports can be output at the same time. Refer to "12. Data Memory" for details.
- The ID number is maintained in non-volatile memory even if the AC adapter is removed.
- Refer to "10.7. Clock and Calendar Function" for checking and storing the time and date.
- When an AD-8127 printer or AD-8129TH printer is connected to the balance and outputs the GLP report, the time and date can be added to the report using the clock function of the printer. (Available for balances with software version 1.660 or later) This is very useful for managing the time and date using the AD-8127 or AD-8129TH password lock function to prevent them from being falsified. To output the GLP/GMP compliant report, set the AD-8127 printer or AD-8129TH printer to the dump printing mode.

## 11.1. Setting the ID Number

- 1 Press and hold the RANGE key (for approx. 2 seconds) until bASFnC is displayed, and then release the key.
- 2 Press the RANGE key several times to display id.
- 3 Press the PRINT key. Set the ID number using the following keys.
  - RE-ZERO key.....To set the character of the digit selected.  
Refer to the display character table shown below.
  - RANGE key.....To select the digit to change the character.
  - PRINT key.....To store the new ID number and display bASFnC.
  - CAL key.....To cancel the new ID number and display bASFnC.
- 4 When bASFnC is displayed, press the CAL key to return to the weighing mode.

### Display character table

|   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |   |
|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|---|
| 0 | 1 | 2 | 3 | 4 | 5 | 6 | 7 | 8 | 9 | - | ␣ | A | B | C | D | E | F | G | H | I | J | K | L | M | N | O | P | Q | R | S | T | U | V | W | X | Y | Z |
| 0 | 1 | 2 | 3 | 4 | 5 | 6 | 7 | 8 | 9 | - | ␣ | A | B | C | D | E | F | G | H | I | J | K | L | M | N | O | P | Q | R | S | T | U | V | W | X | Y | Z |

␣ Space

## 11.2. GLP Report

To output the GLP/GMP compliant report to an AD-8121B printer, AD-8127 printer, thermal printer AD-8129TH or to a personal computer, set the "GLP output ( *inFo* )" parameter of the function table to "1" (AD-8121 format), "2" (General format using the balance built-in clock data) or "3" (General format using the external device clock data).

### Notes on outputting the GLP/GMP compliant report to a printer

- Refer to "16.2. Connecting to Peripheral Equipment" for connection to a printer.
- With an AD-8121B printer, use MODE 3. If MODE1 is used, select temporarily the dump print mode by pressing the **[STAT.]** key of the AD-8121B printer.  
With an AD-8127 printer or thermal printer AD-8129TH, use the dump printing mode.  
If the external key printing mode is used, press and hold the **[ENT]** key of the AD-8127 printer or thermal printer AD-8129TH (for approx. 2 seconds) to switch between the external key printing mode and the dump printing mode.  
If the data memory function is used (except *dAtA* 0), the GLP output cannot be output automatically.
- If the time and date are not correct when the balance built-in clock data is output ( *inFo* 1, 2), set the correct time and date in "Clock (CL Adj)" of the function table.
- " *inFo* 3" is available for balances with software version 1.660 or later.

### Sensitivity adjustment report using the internal mass

**Note** If the data memory function is used (except *dAtA* 0), the "Title block" and "End block" cannot be output.

#### Key operation

- 1 Press the **[CAL]** key to display **[CAL in]**. The balance performs sensitivity adjustment using the internal mass.
- 2 If GLP output is used, **[GLP]** is displayed and the sensitivity adjustment report is output.
- 3 The balance returns to the weighing mode.

Function setting " *inFo* 1"  
AD-8121 format

```

      A & D
MODEL  BM-300
S/N    01234567
ID     LAB-0123
DATE   2011/02/20
TIME   12:34:56
CALIBRATED (INT.)
SIGNATURE
-----

```

← Manufacturer →  
 ← Model →  
 ← Serial number →  
 ← ID number →  
 ← Date →  
 ← Time →  
 ← Sensitivity adjustment with internal weight →  
 ← Signature →

Function setting " *inFo* 2"  
General format using the balance built-in clock data

```

.....A.&D<TERM>
MODEL.....BM-300<TERM>
S/N.....01234567<TERM>
ID.....LAB-0123<TERM>
DATE<TERM>
.....2011/02/20<TERM>
TIME<TERM>
.....12:34:56<TERM>
CALIBRATED (INT.)<TERM>
SIGNATURE<TERM>
<TERM>
<TERM>
-----<TERM>
<TERM>
<TERM>

```


─ Space, ASCII 20h  
 <TERM> Terminator, CR, LF or CR  
 CR Carriage return, ASCII 0Dh  
 LF Line feed, ASCII 0Ah

## Outputting the clock data of the external device “ inf 3 ”

When the "GLP output ( inf )" parameter of the function table is set to "3", the clock data of external devices such as a printer or a personal computer can be used instead of the balance built-in clock data for the GLP/GMP compliant report. This is useful for standardizing the clock data using the clock function of an external device.

### Notes

- To output the clock data of an external device, the external device such as an AD-8127 printer or thermal printer AD-8129TH or Windows communication tools software WinCT RsCom must have a function to output the time and date when <ESC>D or <ESC>T is received.
- Even with inf 3 set, the data memory function saves the balance built-in clock data for the sensitivity adjustment report.

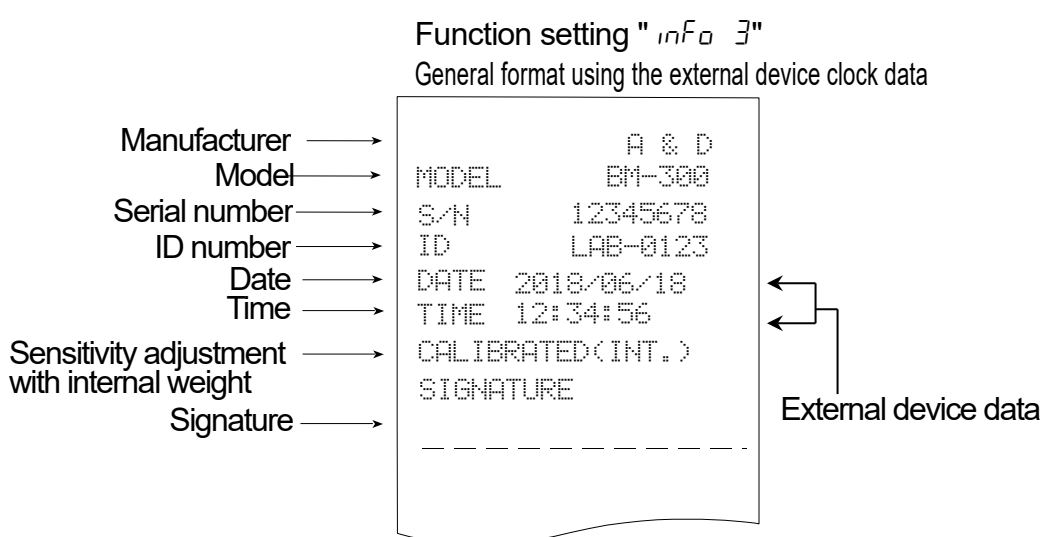

Calibration test report using the internal mass

- Note**
- Calibration test does not perform adjustment.
  - If the data memory function is used (except *data 0*), the "Title block" and "End block" cannot be output.

Key operation

- Press and hold the **CAL** key to display **[[ in ]]** and release the key.
- [[ ]]** is displayed and the balance is checked automatically.
- The zero point is measured and the weight value is displayed for a few seconds.
- The internal mass is weighed and the weight value is displayed for a few seconds.
- If GLP output is used, **[GLP]** is displayed and the calibration test report is output.
- The balance returns to the weighing mode.

Command

This calibration test can be performed with the TST command.

Function setting "inFa 1"  
AD-8121 format

```

      A & D
MODEL    BM-300
S/N      01234567
ID        LAB-0123
DATE     2011/02/20
TIME     12:34:56
CAL. TEST<INT.>
ACTUAL
      0.0000 g
    +200.0002 g
TARGET
    +200.0000 g
SIGNATURE
-----
```

Function setting "inFa 2"  
General format using the balance built-in clock data

```

.....A.&.D<TERM>
MODEL.....BM-300<TERM>
S/N.....01234567<TERM>
ID.....LAB-0123<TERM>
DATE<TERM>
.....2011/02/20<TERM>
TIME<TERM>
.....12:34:56<TERM>
CAL. TEST<INT.><TERM>
ACTUAL<TERM>
.....0.0000...g<TERM>
....+200.0002...g<TERM>
TARGET<TERM>
....+200.0000...g<TERM>
SIGNATURE<TERM>
<TERM>
<TERM>
-----<TERM>
<TERM>
<TERM>
```

- Space, ASCII 20h
- <TERM> Terminator, CR , LF or CR
- CR Carriage return, ASCII 0Dh
- LF Line feed, ASCII 0Ah

## Sensitivity adjustment report using an external weight

**Note** If the data memory function is used (except *data 0*), the "Title block" and "End block" cannot be output.

### Key operation

- Press and hold the **[CAL]** key to display **[CAL out]** and release the key.
- [CAL 0]** is displayed.
- To update the weight value, press the **[RANGE]** key and proceed to Step 4.  
To use the preset weight value, proceed to Step 5.
- Specify the weight value using the following keys.  
**[RANGE]** key.....To select the blinking digit.  
**[RE-ZERO](+)**key...To increase the value of the blinking digit.  
**[MODE](-)**key .....To decrease the value of the blinking digit.  
**[PRINT]** key .....To store the new value.
- Press the **[PRINT]** key. The zero point is measured and the weight value is displayed for a few seconds.
- Place the displayed weight on the pan. Press the **[PRINT]** key to measure the weight and the weight value is displayed for a few seconds.
- [End]** is displayed. Remove the weight from the pan.
- If GLP output is used, **[GLP]** is displayed and the sensitivity adjustment report is output.
- The balance returns to the weighing mode.

Function setting "info 1"  
AD-8121 format

```

          A & D
MODEL    BM-300
S/N      01234567
ID        LAB-0123
DATE     2011/02/20
TIME     12:34:56
CALIBRATED(EXT.)
CAL.WEIGHT
          +200.0000 g
SIGNATURE
          - - - - -
    
```

\_ Space, ASCII 20h  
 <TERM> Terminator, CR , LF or CR  
 CR Carriage return, ASCII 0Dh  
 LF Line feed, ASCII 0Ah

Function setting "info 2"  
General format using the balance built-in clock data

```

          A & D<TERM>
MODEL_____BM-300<TERM>
S/N_____01234567<TERM>
ID_____LAB-0123<TERM>
DATE<TERM>
          2011/02/20<TERM>
TIME<TERM>
          12:34:56<TERM>
CALIBRATED(EXT.)<TERM>
CAL.WEIGHT<TERM>
          ____+200.0000__g<TERM>
SIGNATURE<TERM>
<TERM>
<TERM>
          -----<TERM>
<TERM>
<TERM>
    
```



## Calibration test report using an external weight

**Note** □ Calibration test does not perform adjustment.

- If the data memory function is used (except *data 0*), the "Title block" and "End block" cannot be output.

### Key operation

- 1 Press and hold the **CAL** key to display **EE out** and release the key.
- 2 **CAL 0** is displayed.
- 3 To update the target weight value, press the **RANGE** key and proceed to Step 4.  
To use the preset target weight value, proceed to Step 5.
- 4 Specify the target weight value using the following keys.  
**RANGE** key ..... To select the blinking digit.  
**RE-ZERO**(+)key .... To increase the value of the blinking digit.  
**MODE**(-)key ..... To decrease the value of the blinking digit.  
**PRINT** key..... To store the new value.
- 5 Press the **PRINT** key. The zero point is measured and the weight value is displayed for a few seconds.
- 6 Place the displayed target weight on the pan. Press the **PRINT** key to measure the weight. The weight value is displayed for a few seconds.
- 7 **End** is displayed. Remove the weight from the pan.
- 8 If GLP output is used, **GLP** is displayed and the calibration test report is output.
- 9 The balance returns to the weighing mode.

Function setting "inF0 1"  
AD-8121 format

```

      A & D
MODEL    BM-300
S/N      01234567
ID       LAB-0123
DATE    2011/02/20
TIME     12:34:56
CAL.TEST(EXT.)
ACTUAL
      0.0000 g
      +200.0002 g
TARGET
      +200.0000 g
SIGNATURE
.....

```

Function setting "inF0 2"  
General format using the balance built-in clock data

```

.....A_&_D<TERM>
MODEL.....BM-300<TERM>
S/N.....01234567<TERM>
ID.....LAB-0123<TERM>
DATE<TERM>
.....2011/02/20<TERM>
TIME<TERM>
.....12:34:56<TERM>
CAL. TEST (EXT.)<TERM>
ACTUAL<TERM>
.....0.0000...g<TERM>
.....+200.0002...g<TERM>
TARGET<TERM>
.....+200.0000...g<TERM>
SIGNATURE<TERM>
<TERM>
<TERM>
-----<TERM>
<TERM>
<TERM>

```

- Space, ASCII 20h
- <TERM> Terminator, CR , LF or CR
- CR Carriage return, ASCII 0Dh
- LF Line feed, ASCII 0Ah


## Title block and end block

When the weighing data are recorded as GLP data, a "Title block" is inserted at the beginning and an "End block" is inserted at the end of a group of weight values in the GLP report.

**Note** If the data memory function is used (except *data 0*), the "Title block" and "End block" cannot be output.

### Key operation

- 1 With the weighing data displayed, press and hold the **PRINT** key (for approx. 2 seconds) until **Start** is displayed, and then release the key. The "Title block" is output.
- 2 The weighing data is output according to the parameter setting of the data output mode (*Pr*) of the function table.
- 3 Press and hold the **PRINT** key (for approx. 2 seconds) until **RecEnd** is displayed, and then release the key. The "End block" is output.

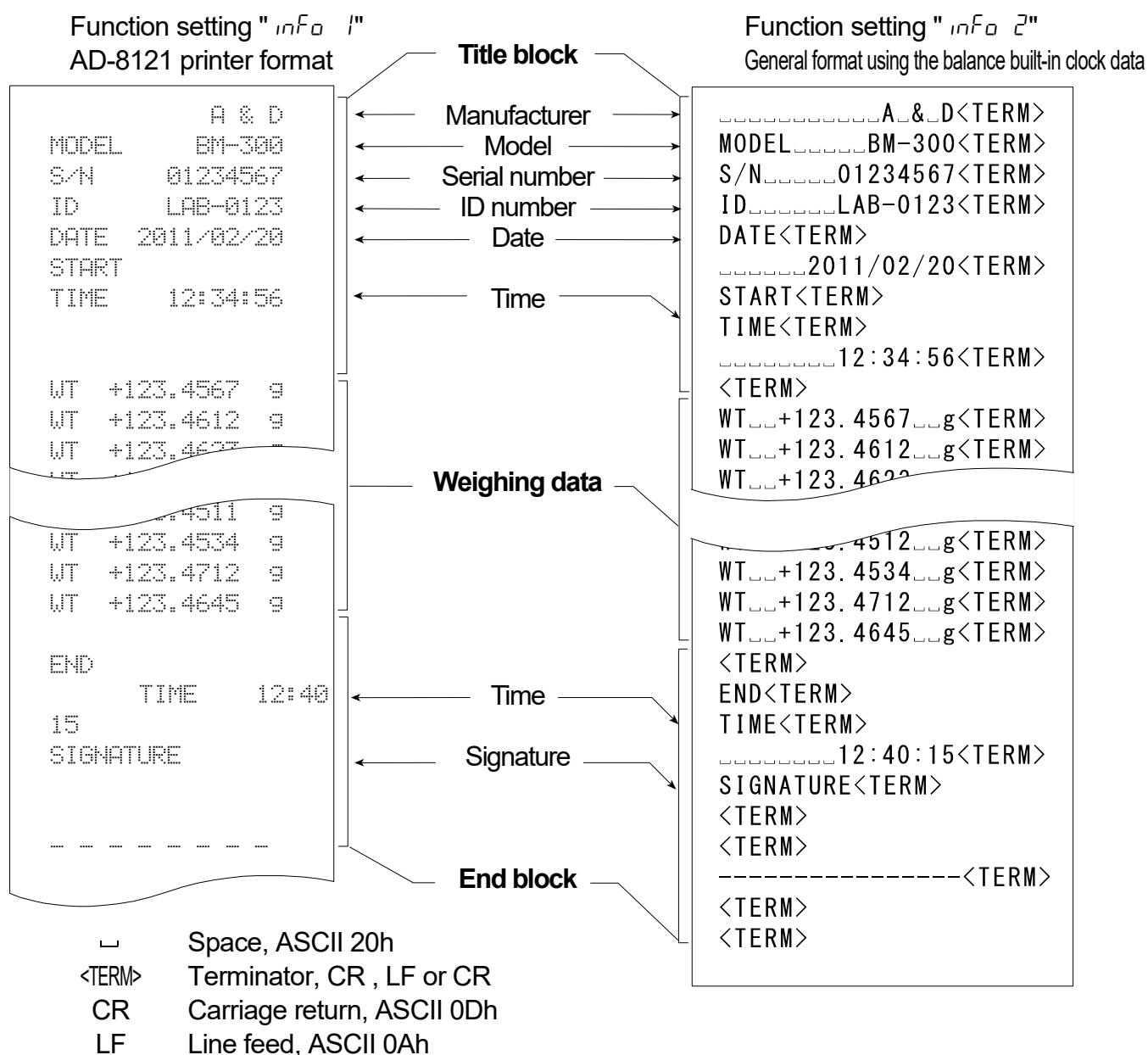

## 12. Data Memory

Data memory is a function to store weighing data, sensitivity adjustment and calibration test in memory. The data stored in memory are available for outputting at one time to a printer or personal computer.

|   |                             |                |
|---|-----------------------------|----------------|
| Weighing data   | Excluding the time and date | Up to 200 sets |
|   | Including the time and date | Up to 100 sets |
| Sensitivity adjustment report:<br>Using the internal mass and an external weight Automatic sensitivity adjustment |                             | Last 50 sets   |
| Calibration test report:<br>Using the internal mass and an external weight  |                             |                |

### 12.1. Notes on Using Data Memory

- To use the memory function, set the "Data memory (dAtA)" parameter and the "Time/Date output (5-t-d)" parameter of the function table. Refer to "10. Function Table" for details on setting the data memory.
- For weighing data, the data contents to be stored and the storage capacity depend on the "Time/Date output (5-t-d)" parameter setting.
- If a different type of data exists in memory, "Lr" blinks in the upper left of the display to prompt you to delete the data.  
For example: You want to store weighing data but sensitivity adjustment report or unit mass data remains in memory. "Lr" blinks.
- If the stored data is corrupted, "Err" blinks in the upper left of the display.  
Under such a condition, before storing data, delete the data in memory as follows:

Upper left of  
the display



#### Releasing "Lr" or "Err"

- 1 Press and hold the **PRINT** key (for approx. 2 seconds) until **Lr na** with "na" blinking is displayed, and then release the key.
- 2 Press the **RE-ZERO** key to display **Lr 0.0** with "0.0" blinking.
- 3 Press the **PRINT** key to delete all the data in memory.
- 4 The balance displays **End** and returns to the weighing mode.

### 12.2. Data Memory for Weighing Data

#### Features

- The balance can store 200 sets of weighing data in memory (if the time and date are added, the balance can store 100 sets). Even if the AC adapter is removed, the data is maintained in the memory.
- It is not necessary to connect a printer or personal computer to the balance continually, because the balance stores the weighing data in memory.
- The data memory function does not occupy a printer or personal computer while weighing.
- The data in memory can be displayed on the balance for confirmation.
- The data in memory can be output at one time to a printer or personal computer.  
Comment: Select items for addition to the output data, such as the ID number, data number, time and date, in the function settings.

## Storing the weighing data

**Note** If "Lr" blinks in the upper left of the display, delete the data in memory.

- 1 Set the "Data memory (dAtA)" parameter to "1" (To store the weighing data).
- 2 Specify whether or not to add the time and date using the "Time/Date output (5-t d)" parameter.
- 3 The storing mode depends on the "Data output mode (Prt)" parameter setting.  
Four modes are available to store data.

Key mode ..... When the **PRINT** key is pressed and the displayed value is stable, the balance stores the weighing data.


Auto print mode A ..... When the displayed value is stable and the conditions of "Auto print polarity", "Auto print difference" and zero point (reference value) are met, the balance stores the weighing data.

Auto print mode B ..... When the displayed value is stable and the conditions of "Auto print polarity", "Auto print difference" and last stable data (reference value) are met, the balance stores the weighing data.


Interval mode ..... Weighing data is stored at an interval specified in "Interval time (int)".  
Press the **PRINT** key to start and stop this mode.

## Enabling the data memory function

- 1 Press and hold the **RANGE** key (for approx. 2 seconds) until **bASFnC** is displayed, and then release the key.
- 2 Press the **RANGE** key several times to display **dout**.
- 3 Press the **PRINT** key.
- 4 Press the **RANGE** key several times to display **dAtA**.
- 5 Press the **RE-ZERO** key to display **dAtA 1**.
- 6 Press the **PRINT** key to store the setting.
- 7 Press the **CAL** key to return to the weighing mode.

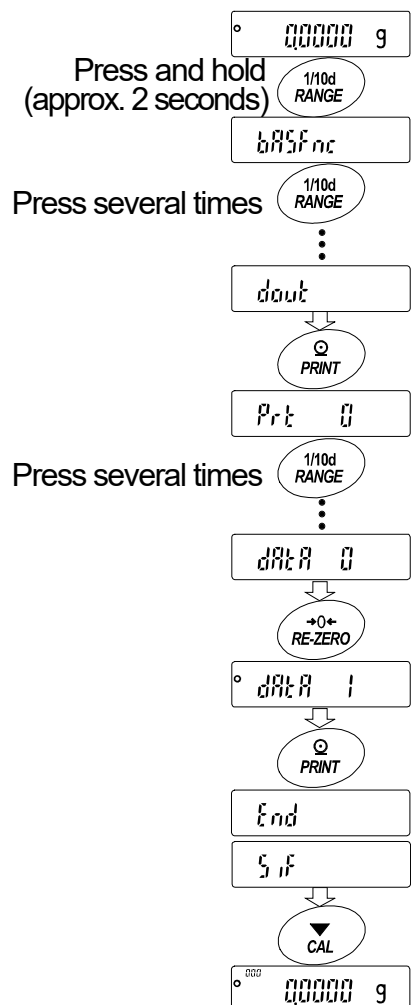

## Display symbols for data memory

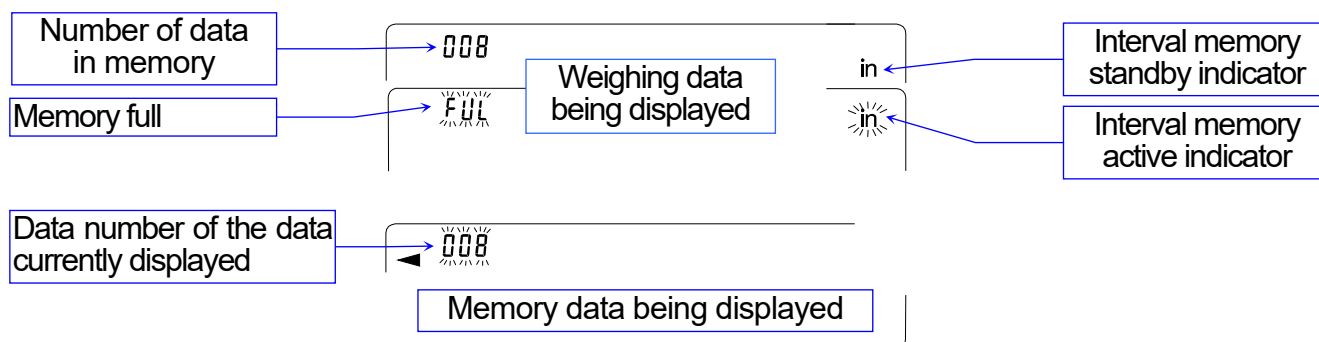

### Notes

- When weighing data is being stored in memory, the data can be output to a personal computer using the RS-232C interface at the same time.
- "FUL" indicates that memory is full or the memory capacity has been reached. More data cannot be stored unless the memory data is deleted.
- Automatic sensitivity adjustment cannot be used while the interval memory mode is active.
- The following commands cannot be used during data storage.

## Setting the function table

Parameter settings for each output mode are as follows:

| Mode \ Item       | Data output mode | Auto print polarity, difference | Data memory function | Interval time |
|-------------------|------------------|---------------------------------|----------------------|---------------|
| Key mode          | Print 0          | Not used                        | dAtA 1               | Not used      |
| Auto print mode A | Print 1          | RP-A 0 to 2                     | dAtA 1               |               |
| Auto print mode B | Print 2          | RP-b 0 to 2                     | dAtA 1               |               |
| Interval mode     | Print 3          | Not used                        | dAtA 1               | int 0 to 8    |


Additional parameter settings, as follows:

|             |     |          |               |           |        |          |
|-------------|-----|----------|---------------|-----------|--------|----------|
| Data number | No  | d-na oFF | Time and date | No        | S-td 0 | 200 sets |
|             | Yes | d-na on  |               | Time only | S-td 1 | 100 sets |
| ID number   | No  | S-id oFF |               | Date only | S-td 2 |          |
|             | Yes | S-id on  |               | Both      | S-td 3 |          |

## Recalling the memory data

Confirm that the "Data memory (dAtA)" parameter is set to "I".

- 1 Press and hold the **PRINT** key (for approx. 2 seconds) until **rEALL** is displayed, and then release the key. The type of data appears in the upper left of the display as shown to the right.
- 2 Press the **PRINT** key to enter the memory recall mode.  
Recall the data in memory using the following keys.  
**RE-ZERO** key.....To proceed to the next data set.  
**MODE** key .....To go back to the previous data set.  
**PRINT** key.....To transmit the current data using the RS-232C interface.  
 With **RANGE** held down, press the **CAL** key  
 To delete the current data.  
**CAL** key.....To exit the memory recall mode.
- 3 Press the **CAL** key to return to the weighing mode.

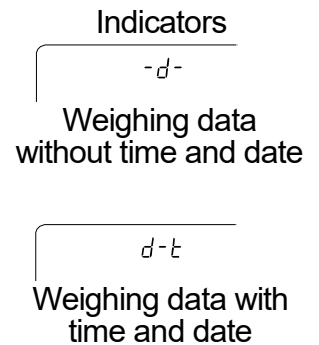


## Transmitting all memory data at one time

Confirm that the "Serial interface (SIF)" parameters are set properly.


Refer to "10. Function Table" and "16.2. Connecting to Peripheral Equipment.".

- 1 Press and hold the **PRINT** key (for approx. 2 seconds) until **rEALL** is displayed, and then release the key.
- 2 Press the **RANGE** key to display **out**.
- 3 Press the **PRINT** key to display **out na** with "na" blinking.
- 4 Press the **RE-ZERO** key to display **out 00** with "00" blinking.
- 5 Press the **PRINT** key to transmit all data using the RS-232C interface.
- 6 The balance displays **CLER** when all data is transmitted.  
Press the **CAL** key to return to the weighing mode.

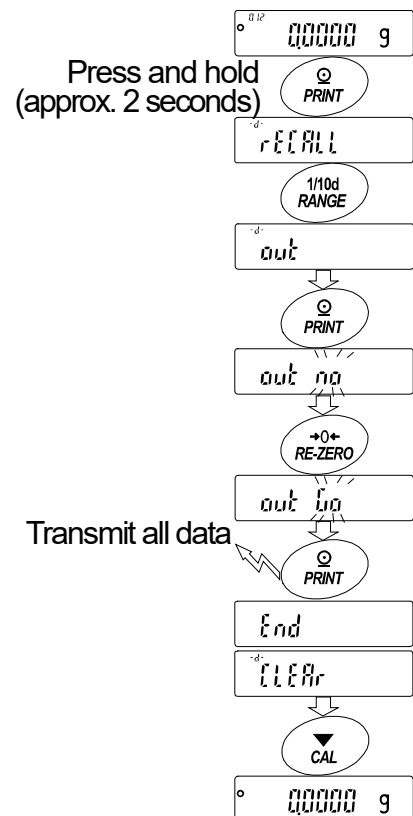

## Deleting all memory data at one time

- 1 Press and hold the **PRINT** key (for approx. 2 seconds) until **rEEALL** is displayed, and then release the key.
- 2 Press the **RANGE** key several times to display **CLER**.
- 3 Press the **PRINT** key to display **CLr na** with "na" blinking.
- 4 Press the **RE-ZERO** key to display **CLr 00** with "00" blinking.
- 5 Press the **PRINT** key to delete all data.
- 6 The balance displays **End** and returns to the weighing mode.

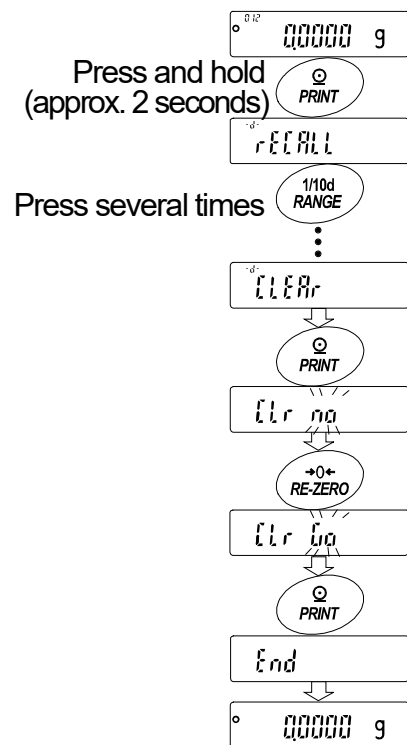

## 12.3. Data Memory for Sensitivity adjustment and Calibration Test Data

### Features

- Sensitivity adjustment history (when and how it is performed) and calibration test data can be stored in memory.
- All the data in memory is available to be output at one time to a printer or personal computer.
- Up to 50 data sets of the latest sensitivity adjustment or calibration test can be stored. When the memory capacity has been reached, "FUL" is displayed in the upper left of the display.

Indicator

FUL

### Storing the sensitivity adjustment and calibration test data

**Note** If "CLr" appears blinking in the upper left of the display, delete the data in memory.

Store the sensitivity adjustment and calibration test data as follows:


- 1 Set the "Data memory (dAtA)" parameter to "2" (to store the sensitivity adjustment and calibration test data).
- 2 Set the "GLP output (GLPo)" parameter to "1", "2" or "3".
- 3 With the settings above, each time sensitivity adjustment or calibration test is performed, the data is stored automatically.

## Transmitting the memory data

### Notes


- Confirm that the "Serial interface (S/F)" parameters are set properly.  
Refer to "10. Function Table" and "16.2. Connecting to Peripheral Equipment".
- Confirm that the "Data memory (dAtA)" parameter is set to "2".

- 1 Press and hold the **PRINT** key (for approx. 2 seconds) until **out** is displayed, and then release the key.
- 2 Press the **PRINT** key to display **out no** with "no" blinking.
- 3 Press the **RE-ZERO** key to display **out 0.0** with "0.0" blinking.
- 4 Press the **PRINT** key to transmit all memory data using the RS-232C interface.
- 5 The balance displays **Clear** when all memory data is output.  
Press the **CAL** key to return to the weighing mode.

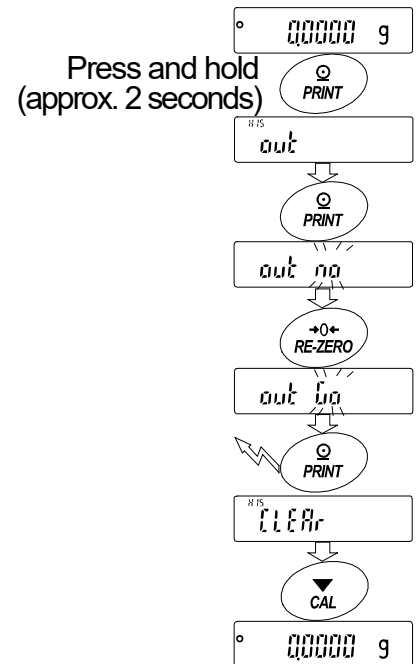

## Deleting data stored in memory

- 1 Press and hold the **PRINT** key (for approx. 2 seconds) until **out** is displayed, and then release the key.
- 2 Press the **RANGE** key to display **Clear**.
- 3 Press the **PRINT** key to display **Clr no** with "no" blinking.
- 4 Press the **RE-ZERO** key to display **Clr 0.0** with "0.0" blinking.
- 5 Press the **PRINT** key to delete all data.
- 6 The balance displays **End** and returns to the weighing mode when all the data has been deleted.

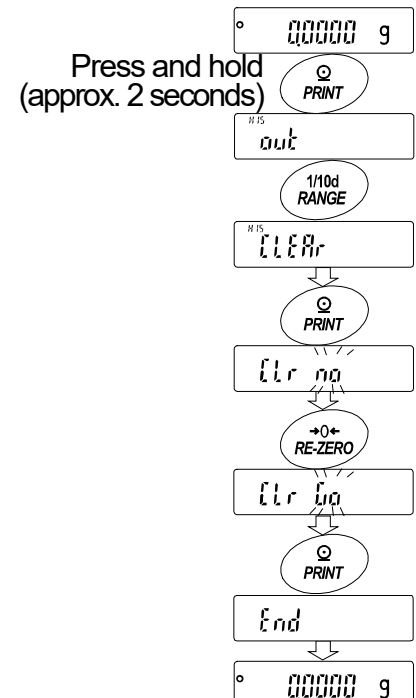





## 13. Underhook

The underhook can be used for magnetic materials or density measurement. The built-in underhook is revealed by removing the cap on the bottom of the balance. Use the underhook as shown below.

### Caution

- Do not apply excessive force to the underhook.
- When not in use, attach the cap to prevent dust from getting into the balance.
- Do not push the underhook upward
- When turning the balance over, parts and weighing sample in the weighing chamber will fall off. (Parts: the weighing pan, pan support, breeze break ring, fine range ring and dust plate ) Remove them first.
- When the cap is removed, a weighing error may occur, because of drafts entering into the internal portion of the balance. Arrange the room condition.

- 1 Remove the cap on the bottom of the balance.
- 2 Hang your weighing pan on the underhook.  
Place the balance on the solid table.

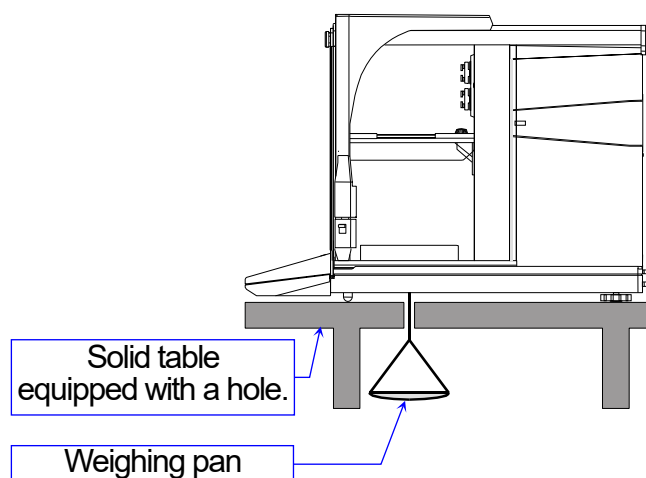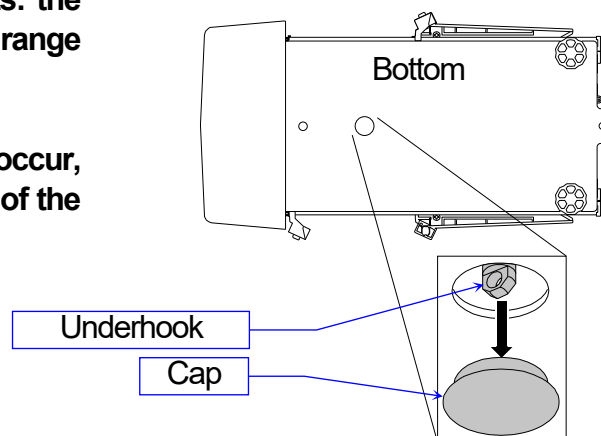

## 14. Density Measurement (Specific gravity)

The balance is equipped with a density mode. It calculates the density of a solid using the weight of a sample in air and the weight in liquid. We recommend the use of the AD-1653 Density Determination Kit.

### Notes

- The density mode was not selected for use when the balance was shipped from the factory. To use this mode, change the function table settings to activate the density mode. Refer to "[5.2. Storing Active Units](#)".
- When the density mode is selected, the response adjustment function cannot be used.
- The readability is 0.0001 g while using the density mode.

### Formula to obtain the density

The density can be obtained by the following formula.

$$\rho = \frac{A}{A - B} \times \rho_0$$

Where

- $\rho$  : Density of a sample
- A : Weight value of a sample in air
- B : Weight value of a sample in liquid
- $\rho_0$  : Density of a liquid

### Prior to measurement: Changing the function table

Prior to measurement, change the function table as follows:

- 1 Selecting the density mode.  
The density mode is not available in the factory setting. To use the mode, select it (unit of Unit 15.) in the function table. Refer to "[5.2. Storing Active Units](#)". Then, the density mode can be selected as one of the weighing units by pressing the MODE key.
- 2 Selecting a way to set the density of a liquid.  
Select the liquid density input method from the function table below.

**Note** The following function table is displayed only when the unit of Unit 15 is active. The table of the density function (d5 Fnc) is displayed after 5 rF.

| Class  | Item and Parameter  |   | Description       |
|--|---|---|-------------------|
| <span style="border: 1px solid black; padding: 0 2px;">d5 Fnc</span><br>Density function | <span style="border: 1px solid black; padding: 0 2px;">Ld in</span><br>Liquid density input | <span style="border: 1px solid black; padding: 0 2px;">■ □</span> | Water temperature |
|  |   | <span style="border: 1px solid black; padding: 0 2px;">/</span>   | Liquid density    |

▪ Factory setting.

## Entering the density of a liquid


Two ways to set the density of a liquid are available in the function table, "Liquid density input ( $Ld_{in}$ ): by entering the water temperature or by entering the density directly.

- Press the **MODE** key as necessary to select the density mode.  
Press and hold the **MODE** key (for approx. 2 seconds) to select to "water temperature" or "liquid density". In the density mode : The unit is "g (gram)". The processing indicator (◀) blinks. Response adjustment is off.  
If you cancel the density mode, press the **MODE** key.

00000 g

### Entering the water temperature ( $Ld_{in}$ °)

- Press and hold the **MODE** key (for approx. 2 seconds). The water temperature currently set (unit: °C, factory setting : 25°C) is displayed. Use the following keys to change the value.

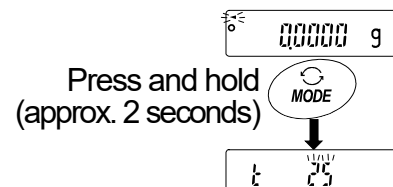

**RE-ZERO**(+)key .... To increase the temperature by one degree.  
(0°C is displayed after 99°C)

**MODE**(-)key ..... To decrease the temperature by one degree.  
(99°C is displayed after 0°C)

**PRINT** key..... To store the new water temperature, display **End** and return to the density mode. Proceed to Step 1.

**CAL** key ..... To cancel the change and return to the density mode.  
Proceed to Step 1.

### The relation between the water temperature and density


| °C | +0      | +1      | +2      | +3      | +4      | +5      | +6      | +7      | +8      | +9      |
|----|---------|---------|---------|---------|---------|---------|---------|---------|---------|---------|
| 0  | 0.99984 | 0.99990 | 0.99994 | 0.99996 | 0.99997 | 0.99996 | 0.99994 | 0.99990 | 0.99985 | 0.99978 |
| 10 | 0.99970 | 0.99961 | 0.99949 | 0.99938 | 0.99924 | 0.99910 | 0.99894 | 0.99877 | 0.99860 | 0.99841 |
| 20 | 0.99820 | 0.99799 | 0.99777 | 0.99754 | 0.99730 | 0.99704 | 0.99678 | 0.99651 | 0.99623 | 0.99594 |
| 30 | 0.99565 | 0.99534 | 0.99503 | 0.99470 | 0.99437 | 0.99403 | 0.99368 | 0.99333 | 0.99297 | 0.99259 |
| 40 | 0.99222 | 0.99183 | 0.99144 | 0.99104 | 0.99063 | 0.99021 | 0.98979 | 0.98936 | 0.98893 | 0.98849 |

g/cm<sup>3</sup>

### Entering the density directly ( $Ld_{in}$ l)

Press and hold the **MODE** key (for approx. 2 seconds). The density currently set (unit : g / cm<sup>3</sup>, factory setting : 1.0000g / cm<sup>3</sup>) is displayed. Use the following keys to change the value.

The range to set the density is 0.0000g / cm<sup>3</sup> to 1.9999g / cm<sup>3</sup>.

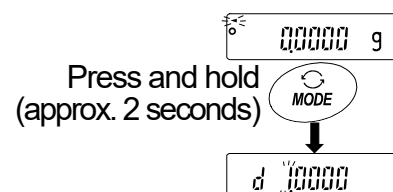

**RE-ZERO**(+)key..... To increase the value of the digit selected.

**MODE**(-)key ..... To decrease the value of the digit selected.

**RANGE** key..... To select the digit to change the value.

**PRINT** key..... To store the change, display **End** and return to the density mode. (Step 1).

**CAL** key ..... To cancel the change and return to the density mode. (Step 1)

## Measuring the density ( specific gravity )

**Note** If the liquid temperature or the type of liquid is changed during measurement, input the value of the liquid density again as necessary. The density display is four decimal places. The readability cannot be changed with the **[RANGE]** key.

The density (or specific gravity) is displayed after "Weighing in air" and "Weighing in liquid". The procedure of each measurement is as follows:

g : gram. ◀ : the processing indicator

1 Enter the density mode. (g lights. ◀ blinks.)  
Place nothing on both pans and press the **[RE-ZERO]** key to display zero.

2 Place the sample on the pan in air (g lights. ◀ blinks.)  
and wait for the weight value to be stable.  
To store or output the weight value, press the **[PRINT]** key.  
Press the **[RANGE]** key to confirm the weight value in air and proceed to the next step.

**Note** If negative value or **[E]** (out of range) is displayed, the **[RANGE]** key is inactive.


3 Move the sample to the pan in liquid (g lights. ◀ lights.)  
and wait for the weight value to be stable.  
To store or output the weight value, press the **[PRINT]** key.  
Press the **[RANGE]** key to confirm the weight value in liquid and proceed to the next step.

**Note** If **[E]** (out of range) is displayed, the **[RANGE]** key is inactive.

4 To store or output the density value, press the **[PRINT]** key. (Unit : 1/15 g turns off. ◀ lights.)  
Press the **[RANGE]** key to measure another sample and proceed to Step 2.

5 If the liquid temperature or the type of liquid is changed during measurement, input the value of the liquid density again as necessary.

6 Press the **[MODE]** key to proceed to other modes.

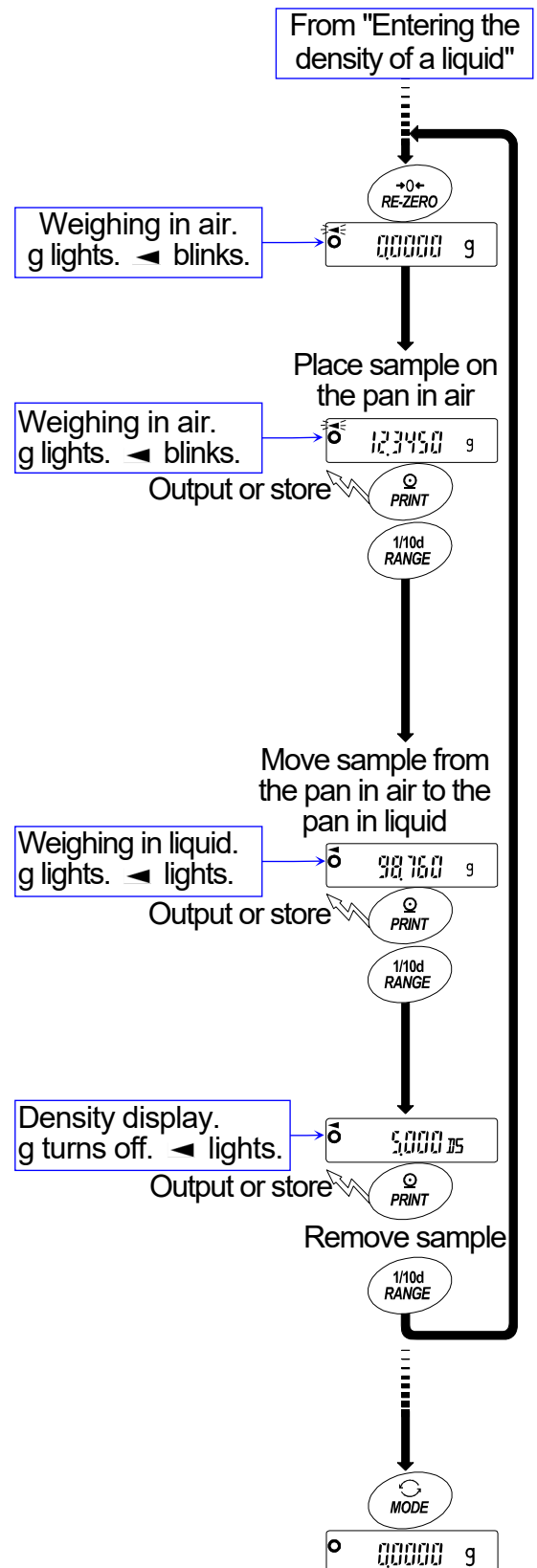

## 15. Password Function

The password function can be used to restrict balance use or functions. It is useful for preventing the time and date setting from being falsified and preventing users from changing other function settings.

Password: Four digits ( $4 \times 4 \times 4 \times 4 = 256$  outcomes) using the **MODE**, **RANGE**, **PRINT** and **RE-ZERO** keys.

At the factory setting, the password function has been disabled.

Settings related to the password function are performed in the function table.

**Note** The password function varies depending on the balance software version.

### 15.1. Balance Software Version 1.50 to 1.64

Two settings are available in “Password function (**LocFunc**)” of the function table.

|               |   |
|---------------|---|
| <b>PASS 0</b> | No password function.                                     |
| <b>PASS 1</b> | Entering a password is required before starting weighing. |

**PASS 0** (No password function)

No password is required. Anybody can perform weighing, use all the balance functions and make changes to them.

**PASS 1** (Entering a password is required before starting weighing.)

The administrator (**Admin**) sets a unique password to restrict balance users. Only one password by the administrator can be set.

The correct password must be entered for the balance to enter the weighing mode.

## 15.2. Balance Software Version 1.660 or later

Three settings are available in “Password function (*LoCFunc*)” of the function table.

|               |   |
|---------------|---|
| <i>PASS 0</i> | No password function.   |
| <i>PASS 1</i> | Entering a password is required before starting weighing.                   |
| <i>PASS 2</i> | Logging in by the administrator password is required before making changes. |

### *PASS 0* (No password function)

No password is required. Anybody can perform weighing, use all the balance functions and make changes to them.

### *PASS 1* (Entering a password is required before starting weighing.)

The administrator (*Admin*) sets unique passwords to restrict balance users. Pressing the ☐ ON:OFF key before weighing will prompt you to enter a password.

The correct password must be entered for the balance to enter the weighing mode.

Two login levels are available: administrator (*Admin*) and user (*USER 01 to 10*).

|                                |   |
|--------------------------------|---|
| Administrator ( <i>Admin</i> ) | All the functions and settings can be used.   |
|                                | A password can be set for each of 10 users.   |
| User ( <i>USER 01 to 10</i> )  | Initialization and changes to the settings are restricted, including the clock setting. |
| No password                    | The balance cannot be used.   |

### *PASS 2* (Logging in by the administrator password is required before making changes.)

Anybody can perform weighing, but initialization and changes to the settings are restricted, including the clock setting. Even if the ☐ ON:OFF key is pressed to start weighing, a password will not be required.

Two login levels are available: administrator (*Admin*) and guest (*GUEST*).

|  |   |
|--|---|
| Administrator ( <i>Admin</i> )         | All the functions and settings can be used.   |
| Guest ( <i>GUEST</i> )<br>*No password | Initialization and changes to the settings are restricted, including the clock setting. |

When weighing is started by pressing the ☐ ON:OFF key while holding down the ☐ CAL key in the standby mode, the balance will prompt you to enter the administrator password.

## Items restricted depending on the login level

| Login level                    | Weighing                          |                            |                            |
|--------------------------------|-----------------------------------|----------------------------|----------------------------|
|                                | Entering password before weighing | Sensitivity adjustment     | Changes to the settings *1 |
| Administrator ( <i>Admin</i> ) | Required                          | Available                  | Available                  |
| User ( <i>USER 01 to 10</i> )  |                                   | Available or not available | Not available              |
| Guest ( <i>GUEST</i> )         | Not required                      | *2                         |                            |

\*1: Response adjustment, repeatability test, function switch and initialization, function settings (clock setting)


\*2: Available under ordinary circumstances, but not available to users and guests when the administrator sets the function switch to “inhibit”.

## 15.3. Using the Password Function

The password function can be switched among “Not used(0)”, “Used(1)” and “Used(2)” in “Password function (LocFnc)” of the function table.

- 1 In the weighing mode, press and hold the **RANGE** key (for approx. 2 seconds) until **bASFn** is displayed, and then release the key.
- 2 Press the **RANGE** key several times to display **LocFnc**.
- 3 Press the **PRINT** key to display **PASS 0**.  
To cancel the operation, press the **CAL** key.
- 4 Press the **RE-ZERO** key to switch the numbers. Display “1” or “2”.
- 5 Press the **PRINT** key to display **SurE YES No** with “No” blinking (selected).
- 6 Press the **RE-ZERO** key to switch between “YES” and “No”.
- 7 Display **SurE YES No** with “YES” blinking (selected).
- 8 Press the **PRINT** key to store the new setting. The password function can now be used.
- 9 **Loc no** is displayed. To store or change the password, proceed to Step 3 in “15.6. Storing or Changing the Password”.  
Not to store or change the password, press the **CAL** key to return to the weighing mode.

**Note** When the balance software version is 1.50 to 1.64, only “Not used(0)” or “Used(1)” is available.

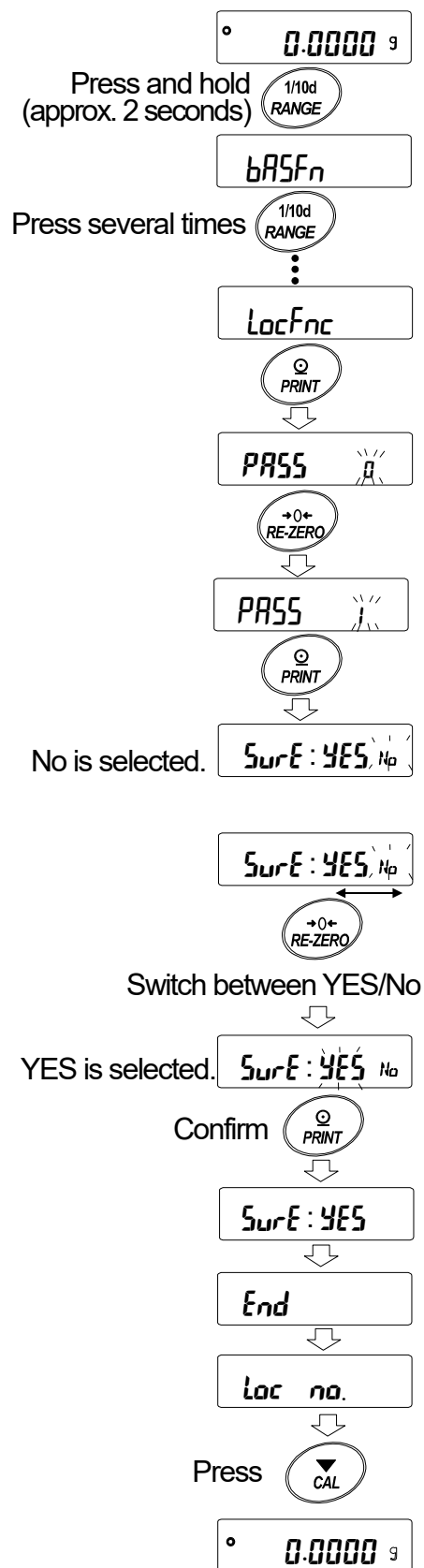


## 15.4. Entering the Password Before Weighing

### 15.4.1. *PASS* 1 (Entering a password is required before starting weighing)

- 1 In the standby mode, press the **ON/OFF** key.
- 2 After **USER** **PASS** **WT**, **---** **PASS** is displayed to prompt you to enter a password.
- 3 Enter a four-digit password using the following keys.  
The display will turn off automatically after ten minutes of non-operation.  

|                    |       |                          |
|--------------------|-------|--------------------------|
| <b>MODE</b> key    | ..... | To enter $\bar{n}$       |
| <b>RANGE</b> key   | ..... | To enter $r$             |
| <b>PRINT</b> key   | ..... | To enter $P$             |
| <b>RE-ZERO</b> key | ..... | To enter $Z$             |
| <b>CAL</b> key     | ..... | To go back one character |

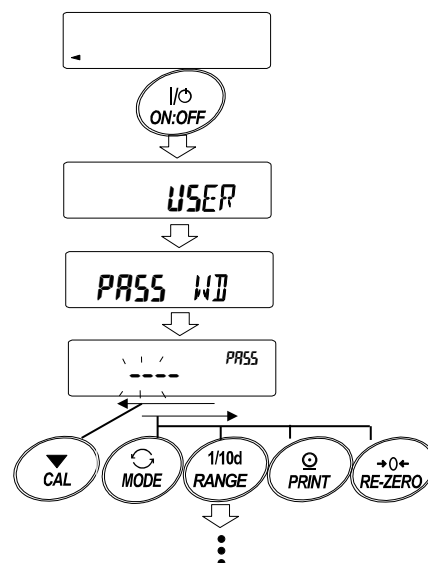


- 4 When the password is correct, the balance displays the login level, turns all the display segments on, and then enters the weighing mode.

When the administrator password is entered, you are logged in as an administrator.

The default password set at the factory is **2222** (press the **RE-ZERO** key four times), for administrator level.

When the password is incorrect, the balance displays **FAIL**, sounds the buzzer three times and turns the display off.

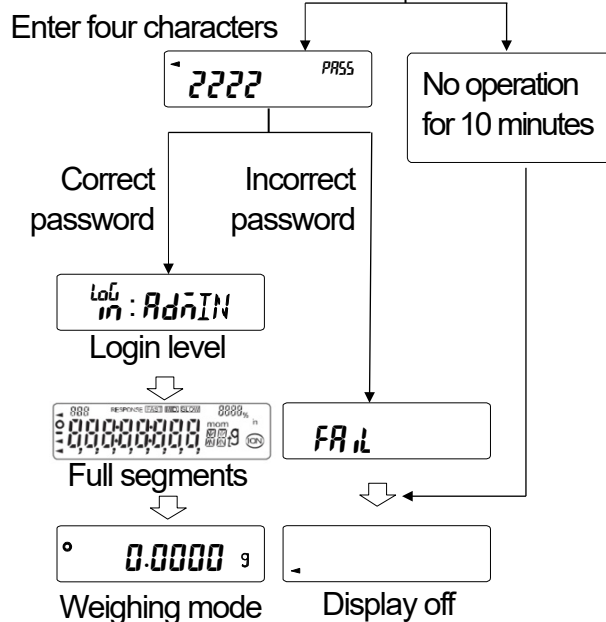

**Note** When the balance software version is 1.50 to 1.64, the login level is not displayed when logged in.  
When the balance software version is 1.660 or later, enter the administrator password to log in as an administrator.



## 15.4.2. PASS 2 (Logging in by the administrator password is required before making changes)

### Logging in as an administrator (AdmIN)

- 1 In the standby mode, while holding down the **CAL** key, press the **ON:OFF** key.

While holding down

- 2 Enter a four-digit password using the following keys.  
The display will turn off automatically after ten minutes of non-operation.

**MODE** key ..... To enter  $\bar{n}$

**RANGE** key ..... To enter  $r$

**PRINT** key ..... To enter  $P$


**RE-ZERO** key ..... To enter  $Z$

**CAL** key ..... To go back one character

- 3 When the password is correct, the balance displays the login level, turns all the display segments on, and then enters the weighing mode.


The default password set at the factory is **2222** (press the **RE-ZERO** key four times), for administrator level.

When the password is incorrect, the balance displays **FAIL**, sounds the buzzer three times and turns the display off.

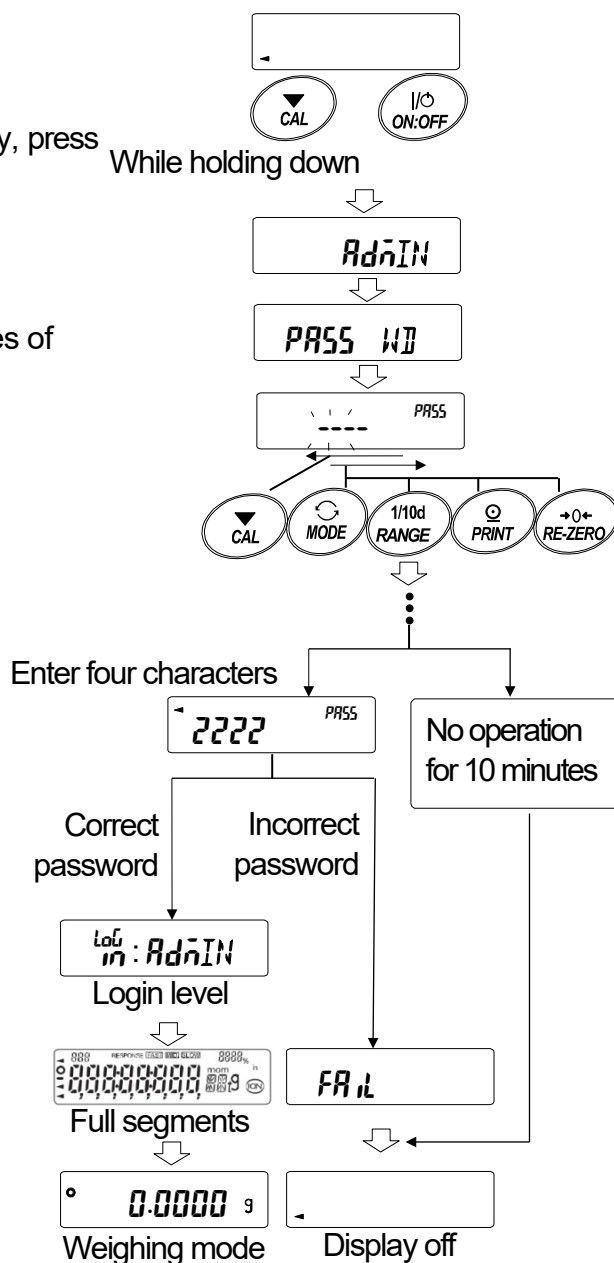


### Logging in as a guest (GUEST)

- 1 In the standby mode, press the **ON:OFF** key.
- 2 After **Log in: GUEST**, the balance enters the weighing mode.

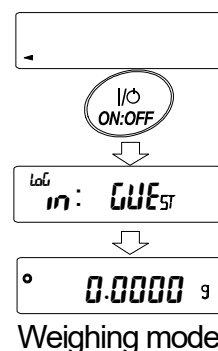

## 15.5. Logging Out


- 1 To log out, press the **ON:OFF** key to turn the display off.  
With the function **PASS 1** set, the password will need to be entered again to enter the weighing mode from the standby mode.

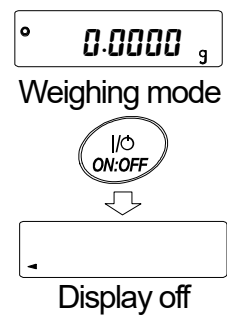

## 15.6. Storing or Changing the Password

The password can be changed in "Password (Loc no.)" of the function table.

- 1 In the weighing mode, press and hold the **RANGE** key (for approx. 2 seconds) until **bPASSn** is displayed, and then release the key.
- 2 Press the **RANGE** key several times to display **Loc no.**
- 3 Press the **PRINT** key to display the login level (**AdnIN**).
- 4 Press the **RANGE** key to display the login level to be changed (**AdnIN/USER 01 to 10**).  
The stabilization indicator appears at a login level where a password has been stored. (The password can be changed.)
- 5 To change the password, press the **PRINT** key. For details, refer to "15.7. Changing the Password".

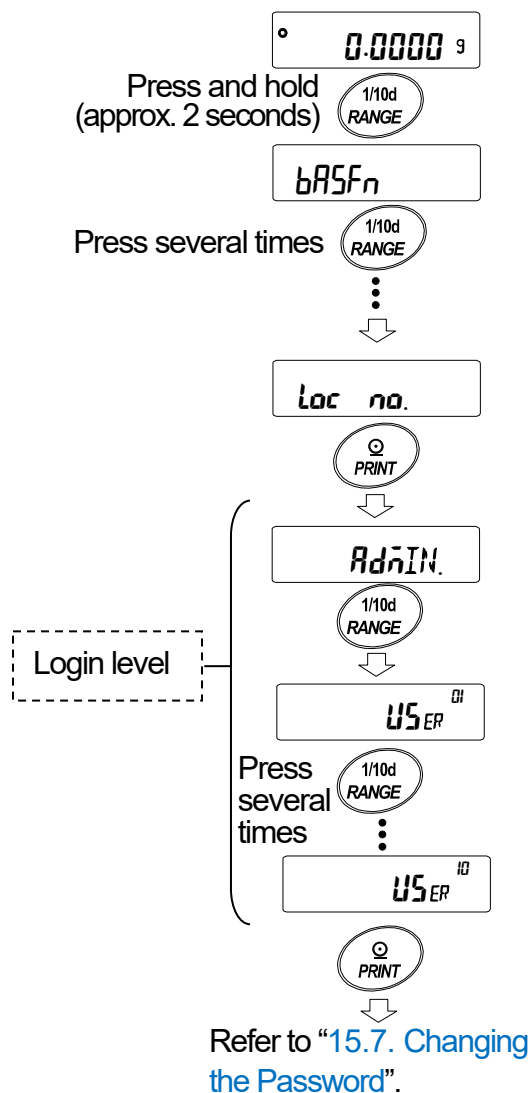

### Notes

- To log out, press the **ON:OFF** key to turn the display off.
- With the function **PASS 2** set, the administrator password is required to log in as an administrator. No password needs to be stored for **USER 01 to 10**.
- When the balance software version is 1.50 to 1.64, only one password can be stored. While the balance displays **Loc no.**, press the **PRINT** key to display the current password.

## 15.7. Changing the Password

- 1 Refer to “15.6. Storing or Changing the Password” to display the login level to be changed.
- 2 Press the **PRINT** key to display the current password.  
The default password set at the factory is **2222** (press the **RE-ZERO** key four times), for administrator level.
- 3 Enter a new four-digit password using the following keys.  
The display will turn off automatically after ten minutes of non-operation.

**MODE** key ..... To enter **7**

**RANGE** key ..... To enter **7**

**PRINT** key ..... To enter **7**

**RE-ZERO** key ..... To enter **7**


**CAL** key ..... To go back one character

**CAL** key (press and hold for approx. 2 seconds)

..... To delete the password

Refer to “15.8. Deleting the Password (USER 01~10)”.

- 4 The new password is displayed.
- 5 **SURE :YES No** is displayed with “No” blinking (selected).  
Press the **CAL** key to return to the fourth character entry position.
- 6 Press the **RE-ZERO** key to display **SURE :YES No** with “YES” blinking (selected).
- 7 While “YES” is blinking, press the **PRINT** key to store the new password.
- 8 When setting is complete, **End** and then the next login level is displayed.  
To set a password for the next level, follow the procedure starting from Step 2.  
To finish the operation, press the **CAL** key twice to return to the weighing mode.

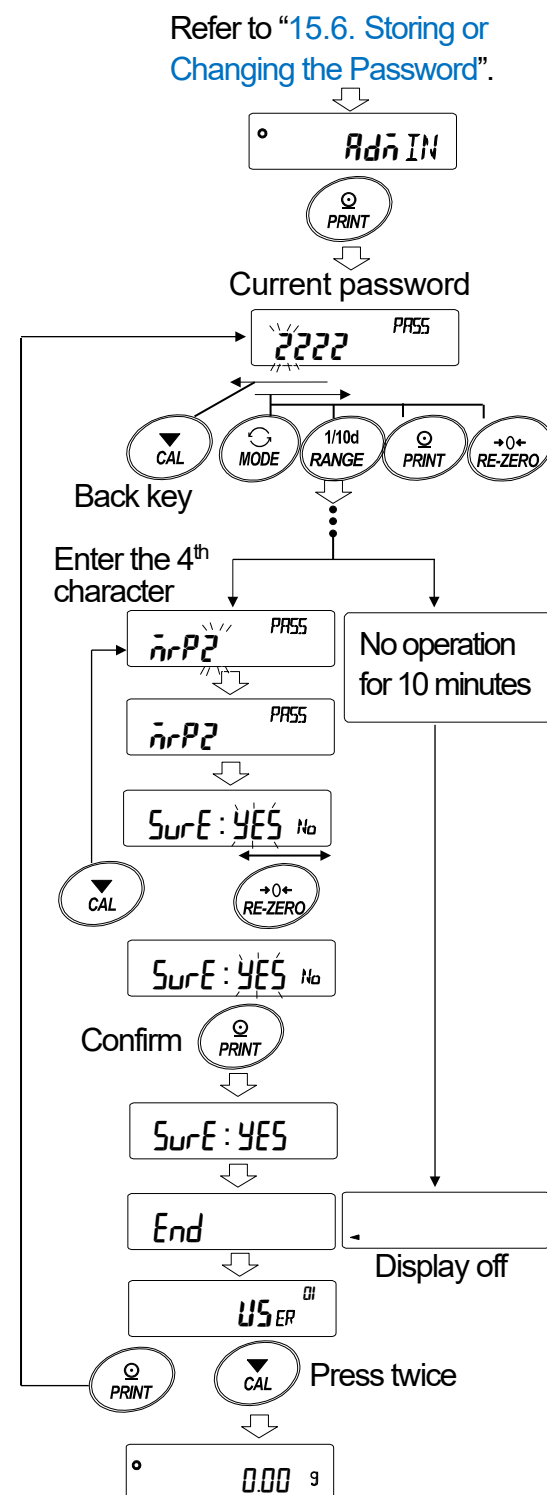


### Notes

- If you have forgotten the password, the balance cannot be used. Make a record of the password and keep it where you can refer to it if needed.
- The same password as the administrator password cannot be used for **USER 01 to 10**.

## 15.8. Deleting the Password (USER 01~10)

- 1 Refer to “15.7. Changing the Password” to select the user whose password to be deleted and display the password.
- 2 Press and hold the **CAL** key (for approx. 2 seconds) to display **CLEAR** (blinking).
- 3 Press the **PRINT** key to display **CLEAR No**.
- 4 Press the **RE-ZERO** key to switch between “No” and “Yes”.
- 5 While **CLEAR Yes** is displayed, press the **PRINT** key to delete the password. The balance displays **End** and then displays the login level.

Refer to “15.7. Changing the Password”.

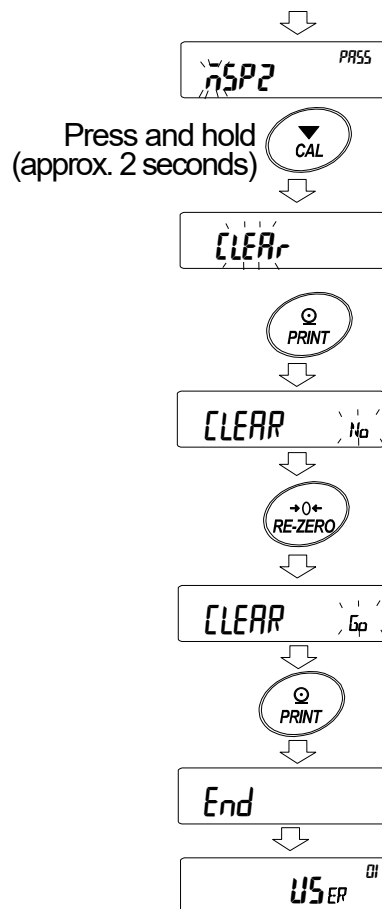

**Note** The administrator password cannot be deleted, but it can be changed.

Refer to “15.6 Storing or Changing the Password” and “15.7 Changing the Password” to change it.

## 15.9. Forgotten Password


If you have forgotten the password, the balance cannot be used.  
Contact your local A&D dealer to undo the password.

# 16. Standard Input and Output Interface

## 16.1. RS-232C Interface


The balance is a DCE device. Connect the balance to a personal computer (DTE) using a straight through cable.

Transmission system : EIA RS-232C  
 Transmission form : Asynchronous, bi-directional, half duplex  
 Transmission rate : 10 times/second or 5 times/second (same as data refresh rate)  
 Data format : Baud rate : 600, 1200, 2400, 4800, 9600, 19200 bps  
                   Data bits : 7 or 8 bits  
                   Parity : Even, Odd (Data bits 7 bits)  
                               None (Data bits 8 bits)  
 Stop bit : 1 bit  
 Code : ASCII

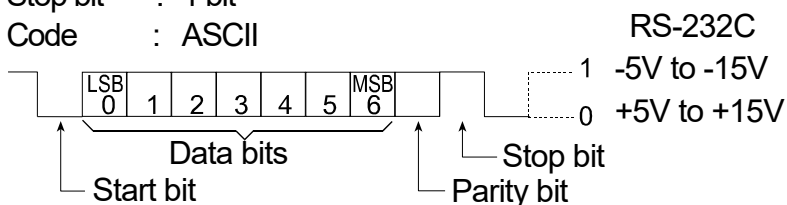


### D-Sub 9 pin assignments

| Pin No. | Signal name | Direction | Description     |
|---------|-------------|-----------|-----------------|
| 1       | —           | —         | No connection   |
| 2       | RXD         | Output    | Transmit data   |
| 3       | TXD         | Input     | Receive data    |
| 4       | —           | —         | No connection   |
| 5       | SG          | —         | Signal ground   |
| 6       | DSR         | Output    | Data set ready  |
| 7       | RTS         | Input     | Request to send |
| 8       | CTS         | Output    | Clear to send   |
| 9       | —           | —         | No connection   |

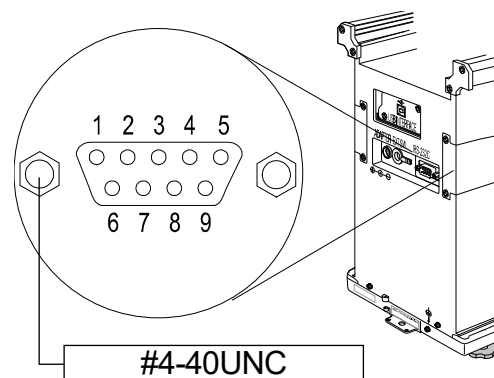

Signal names other than TXD and RXD are DTE-side names.

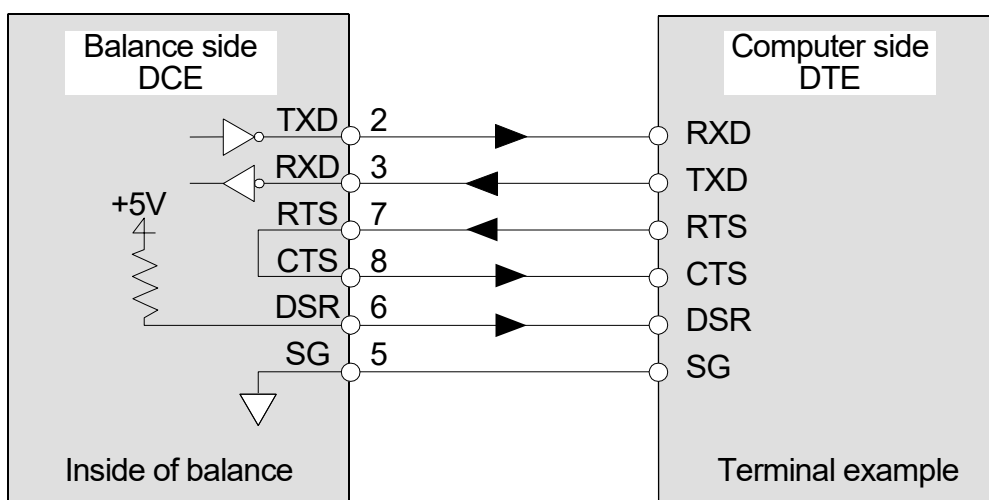

## 16.2. Connecting to Peripheral Equipment

### 16.2.1. Connecting to the AD-8121B Printer

When connecting the AD-8121B printer to the balance for printing weighing data, configure the printer and the balance as follows




| Class            | Item and Parameter                   | Factory settings | AD-8121B MODE 1 | AD-8121B MODE 2 | AD-8121B MODE 3 |
|------------------|--------------------------------------|------------------|-----------------|-----------------|-----------------|
| Data output      | <i>Prt</i><br>Data output mode       | 0                | 0, 1, 2         | 3               | 0, 1, 2         |
|                  | <i>RP-P</i><br>Auto print polarity   | 0                | #1              | No setting      | #1              |
|                  | <i>RP-b</i><br>Auto print difference | 1                |                 |                 |                 |
|                  | <i>S-id</i><br>ID number output      | off              | off             | off             | off, on         |
|                  | <i>d-no</i><br>Data number output    | off              | off             | off             | off, on         |
|                  | <i>S-t-d</i><br>Time/Date output     | 0                | 0               | 0               | 0, 1, 2, 3      |
|                  | <i>PUSE</i><br>Data output pause     | off              | off             | off             | off, on #2      |
|                  | <i>Rt-F</i><br>Auto feed             | off              | off             | off             | off, on         |
| Serial interface | <i>bPS</i><br>Baud rate              | 2                | 2               | 2               | 2               |
|                  | <i>btPr</i><br>Data bit, parity bit  | 0                | 0               | 0               | 0               |
|                  | <i>ErLF</i><br>Terminator            | 0                | 0               | 0               | 0               |
|                  | <i>tYPE</i><br>Data format           | 0                | 0               | 0               | 1               |

#1 Set the parameters when auto print mode A or B (*Prt* 1 or 2) is selected.

#2 Set on when multiple lines are printed. Example: When adding ID number, set on.

#### Notes

- Refer to "11.2. GLP Report" for print samples.
- Settings of AD-8121B

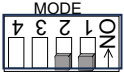
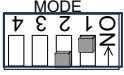
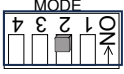
| MODE   | AD-8121B DIP switch   | Description  |
|--------|---|--|
| MODE 1 |  | Print at receiving data.<br>Standard mode, statistical calculation mode  |
| MODE 2 |  | Print by <span style="border: 1px solid black; padding: 0 2px;">DATA</span> key operation or built-in timer.<br>Standard mode, interval mode, chart mode |
| MODE 3 |  | Print at receiving data.<br>Dump print mode  |

DIP switch No.3 : Handling unstable data

- ON To print
- OFF Not to print

DIP switch No.4 : Data input specifications (Interface selection)

- ON Current loop
- OFF RS-232C

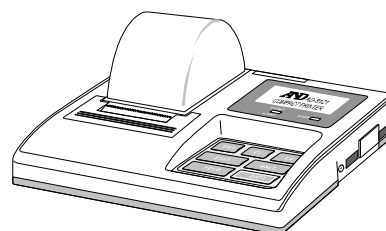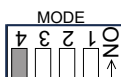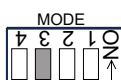

## 16.2.2. Connecting to the AD-8127 Printer or thermal printer AD-8129TH

When connecting the AD-8127 printer or thermal printer AD-8129TH to the balance for printing weighing data, configure the printer and the balance as follows according to the usage examples.

### Function settings of the printer

| Usage example  | Multi printer AD-8127<br>Thermal printer AD-8129TH<br>Printing mode settings |
|--|--|
| When printing the balance weighing data with the balance's <span>PRINT</span> key or its "Auto print" mode.  | EXT.KEY  |
| When printing the balance weighing data with the printer's "Printing" key or its "Interval printing" mode.<br>When printing charts with the printer. | MANUAL<br>AUTO<br>TIMER<br>CHART   |
| When printing the balance's statistical calculation results.<br>When printing the balance's GLP report.  | DUMP   |

- Refer to the instruction manual of the AD-8127 printer or thermal printer AD-8129TH for how to change the function settings of each printer.

### Function settings of the balance

| Usage example  | Balance<br><i>Prt</i><br>Data output mode | Balance<br><i>TYPE</i><br>Data format |
|--|---|---------------------------------------|
| When printing the balance weighing data with the balance's <span>PRINT</span> key or its "Auto print" mode.  | 0, 1, 2                                   | 0                                     |
| When printing the balance weighing data with the printer's "Printing" key or its "Interval printing" mode.<br>When printing charts with the printer. | 3   | 0                                     |
| When printing the balance's statistical calculation results.<br>When printing the balance's GLP report.  | 0, 1, 2                                   | 1                                     |

- Refer to "[10. Function Table](#)" for how to change the function settings of the balance.



## 16.2.3. Connecting to a Computer Using the RS-232C Interface

---

### Features

The balance is of the DCE type (Data Communication Equipment), which can be connected to a personal computer using the RS-232C interface. Before connection, read the personal computer manual thoroughly. Use a standard DCE cable for connection (cable type: straight-through).

### Using Windows Communication Tools Software (WinCT)

When Windows is used as an operating system in a personal computer, the WinCT software can be used to transmit the weighing data to the personal computer. The current version of the WinCT can be downloaded from the A&D website. Confirm the windows version supported by this software when downloading. Refer to the WinCT instruction manual in the A&D website. The WinCT software has three communication methods: "RsCom", "RsKey" and "RsWeight".

#### RsCom

- RsCom can transmit commands to control the balance.
- RsCom can make bi-directional communication between the balance and a personal computer using the RS-232C interface.
- RsCom can display or store the data using a text file format. RsCom can also print the data using a printer connected to the personal computer.
- When several ports of a personal computer have balances connected, the personal computer can communicate with each balance simultaneously.
- RsCom can share a personal computer with other application software.
- RsCom can receive the balance GLP report.

#### RsKey

- RsKey can transmit the weighing data output from the balance directly to other application software such as Microsoft Excel.
- RsKey can be used with most application software.
- RsKey can receive the balance GLP report.

#### RsWeight

- Can retrieve the weight data from the balance and display the data as a graph in real time.
- Can calculate and display the maximum, minimum, average, standard deviation and coefficient of variation.

### Using the WinCT software, the balance can do the following:

- **Analyzing the weighing data and the statistics input by "RsKey"**  
The weighing data can be input directly into an Excel worksheet. Then, Excel can analyze the data to obtain total, average, standard deviation, maximum and minimum values, and display them in a graph.
- **Controlling the balance using commands from a personal computer**  
By using "RsCom", the personal computer sends commands such as "re-zero" or "send weighing data" to the balance and controls the balance.
- **Printing the balance GLP report using your printer**  
The balance GLP report can be printed using a printer connected to the personal computer.
- **Receiving weighing data at a certain interval**  
The weighing data can be received at a certain interval and data characteristic with elapsed time can be obtained.
- **Using the balance memory function**  
The data can be stored in the balance's memory. Of the data stored, the weighing data, sensitivity adjustment data and calibration test data can be transmitted to a personal computer at one time.
- **Using a personal computer as an external indicator**  
With the "RsKey" test mode function, a personal computer can be used as an external weight indicator for the balance. (To do this, set the balance data output mode to stream mode.)

## 16.2.4. Connecting to a Computer Using the USB Interface

### Features

- The weighing data can be transmitted to a personal computer when connecting the balance and personal computer via the USB interface. The applicable operating system is Windows 7 or later, as of December 2018.
- The USB interface uses standard device driver software and does not require installation of special driver software. By simply connecting, weighing data can be transmitted.
- The data can be transmitted to most application software (ex.: EXCEL, WORD, Notepad).

### Notes

- **The balance can transmit the weighing data only without the header and unit.**
- **The USB interface of the balance cannot receive the command that controls the balance. When using the command, connect to the RS-232C interface of the balance. When connecting the USB interface of a personal computer to the RS-232C interface of the balance, use the USB converter (AX-USB-9P).**
- **Do not use the screensaver or suspend mode of a personal computer.**
- **Even if stream mode is used, output will be performed approximately 2 times/second due to key input operation.**
- **To add the ID number, data number, time and date to the weighing data or to output the GLP report, set the “Data output pause (PUSE)” parameter of “Data output (dout)” to “on” (Pause 1.6 seconds). The time and the date are separated by a dot when output.**
- **When the “Data format (LDF)” parameter is set to “5” (CSV format), the ID number, data number, time and date cannot be added to the weighing data.**

### Using the USB interface

- 1 Start the Windows of the personal computer.  
Connect the AC adapter to the balance and press the  key to turn the display on.
- 2 Connect the personal computer to the balance using the USB cable.  
**Note** Do not use stream mode of data output mode (dout: Prt 3).
- 3 At the first connection only, the USB driver software is automatically installed.
- 4 Open the application (ex.: EXCEL) that is to receive the weighing data.  
Move the cursor onto the position to input the weighing data.
- 5 Press the  key to input the weighing data at the position of the cursor.
- 6 To finish the operation, remove the USB cable.

### Examples

- **Case 1: Transmitting the weighing data with key operation**  
Data output mode of the balance: Key mode (factory setting), dout: Prt 0  
Press the  key to transmit the weighing data. (basic use)
- **Case 2: Transmitting the stable weighing data**  
Data output mode of the balance: Auto print mode A and B, dout: Prt 1 or 2  
When a sample is placed and the stabilization indicator  is displayed, the weighing data is transmitted. This mode is useful when repeating the weighing.
- **Case 3: Transmitting the weighing data periodically**  
Data output mode of the balance: Interval mode, dout: Prt 3, dAtA 1 and int  
The weighing data is transmitted periodically.  
The interval time: 2, 5, 10, 30 seconds, 1, 2, 5, 10 minute(s).

## 16.3. Commands

### 16.3.1. Command List

**Note** Before transmitting a command, add the terminator that is specified in the “Terminator (CRLF)” parameter of “Serial interface (SIF)” in the function table.

| Commands to query weighing data |   |
|---------------------------------|---|
| C                               | Cancels the S or SIR command.               |
| Q                               | Requests the weighing data immediately.     |
| S                               | Requests the weighing data when stabilized. |
| SI                              | Requests the weighing data immediately.     |
| SIR                             | Requests the weighing data continuously.    |

| Commands to control the balance |   |
|---------------------------------|---|
| CAL                             | Same as the <b>CAL</b> key. Sensitivity adjustment using the internal mass. |
| OFF                             | Turns the display off.  |
| ON                              | Turns the display on.   |
| P                               | Same as the <b>ON:OFF</b> key. Turning the display on or off.               |
| PRT                             | Same as the <b>PRINT</b> key.   |
| R                               | Same as the <b>RE-ZERO</b> key. Zero display. #1                            |
| RNG                             | Same as the <b>RANGE</b> key. Changing the readability.                     |
| TR                              | Tares the balance. Canceling the container's weight. #1                     |
| TST                             | Performs calibration test. Calibration test using the internal mass.        |
| U                               | Same as the <b>MODE</b> key. Changing the unit.                             |
| KL                              | Sets locking or unlocking of the key lock (function to disable all keys).   |
| LK                              | Sets the specified to be locked (disabled) #1                               |

| Commands to request data |   |
|--------------------------|---|
| ?ID                      | Requests the identification number.   |
| ?PT                      | Outputs the tare value. #1  |
| ?SN                      | Requests the serial number of the balance.                                  |
| ?TN                      | Requests the model name of the balance.                                     |
| ?KL                      | Outputs the status of the key lock (function to disable all keys) settings. |
| ?LK                      | Outputs the status of the specified key lock (disable) setting.             |

| Command to set data  |                         |
|----------------------|-------------------------|
| PT:***.**** <b>g</b> | Sets the tare value. #1 |

- #1
- Command **R** assumes the point as zero and sets the display to zero.
  - Command **TR** cancels the tare weight when the displayed value is greater than zero. The canceled tare value can be read using the command **?PT**.
  - Command **PT:\*\*\*.\*\*\*\* g** sets the tare value digitally.  
An example of this command sets a negative target value and fills with the sample until the display becomes zero. For details, refer to page 95.
- Step 1 Place a container.
- Step 2 Set the display to zero using the **R** command.
- Step 3 Set a target value using the **PT:** command.

### 16.3.2. Acknowledge Code and Error Codes

---

When the "AK, Error code (*ErrCd*)" parameter of "Serial interface function (*SIF*)" is set to "on", the balance outputs <AK> code or error code to each command as follows:

<AK> (06h)                      Acknowledge in ASCII.

- When the balance receives a command to request data and cannot process it, the balance transmits an error code (EC, *Err*).

When the balance receives a command to request data and can process it, the balance outputs the data.

- When the balance receives a command to control the balance and cannot process it, the balance transmits an error code (EC, *Err*).

When the balance receives a command to control the balance and can process it, the balance transmits the acknowledge code.

<AK> (06h)                      Acknowledge in ASCII.

Among commands to control the balance, the following transmit the acknowledge code both when the balance receives the command and when the balance has accomplished the command. If the command cannot be processed properly, the balance transmits an error code (EC, *Err*). This error can be resolved using the CAL command. *xx* is error code number.

|             |  |
|-------------|--|
| CAL command | (Sensitivity adjustment command using internal mass) |
| ON command  | (Display ON command)                                 |
| P command   | (Display ON/OFF command)                             |
| R command   | (RE-ZERO command)                                    |
| TR command  | (Tare command)                                       |
| TST command | (Calibration test command)                           |

- When a communication error has occurred due to external noise, or a parity error has occurred due to transmission error, the balance transmits an error code (EC, *Err*). In this case, send the command again. *xx* is error code number.

### 16.3.3. Settings Related to RS-232C


---

Concerning the RS-232C, the balance has two functions: "Data output (*dout*)" and "Serial interface (*SIF*)". Set each function as necessary.


### 16.3.4. Command Examples

This example uses "ErCd on" of "S iF" so that the <AK> (06h) code is output.


#### ON command (turning the display on)

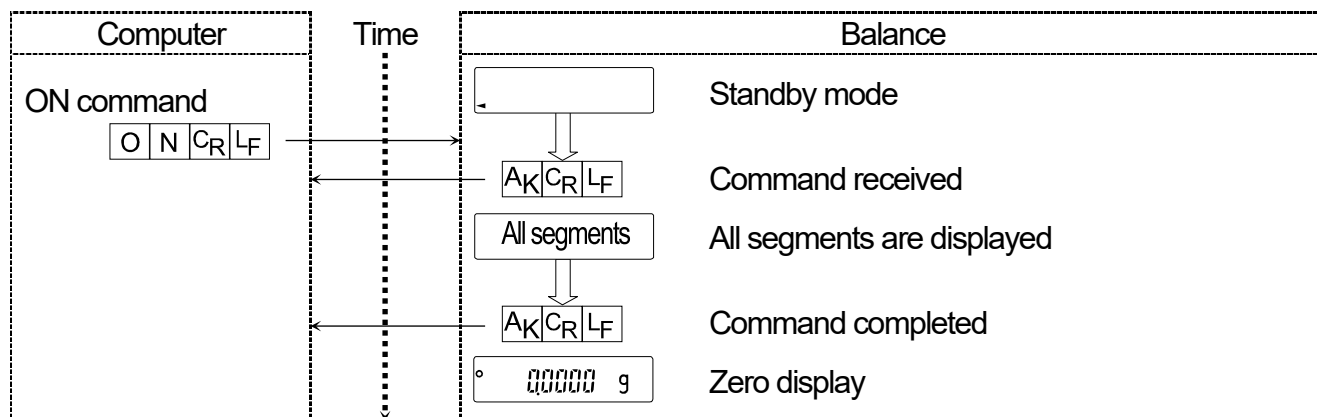

#### R command (Re-zeroing the display)

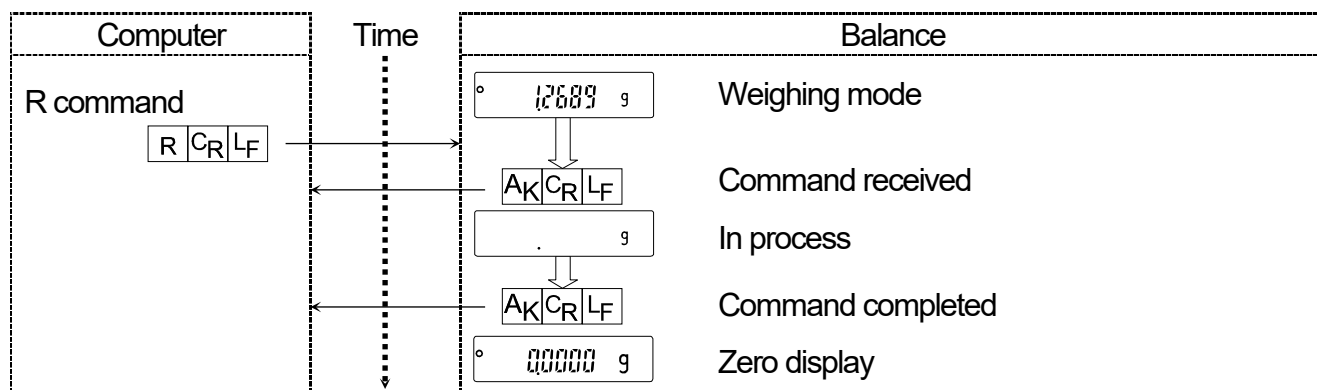


#### CAL command (Sensitivity adjustment with internal mass)

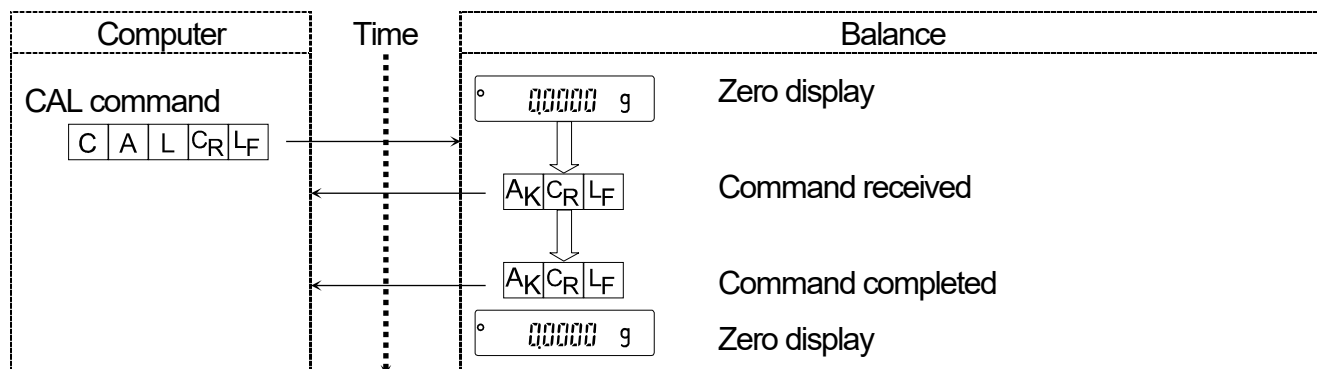


## Error code

This example is of an error using the R command. "Error" is used. The balance transmits an error code when the received command cannot be achieved.

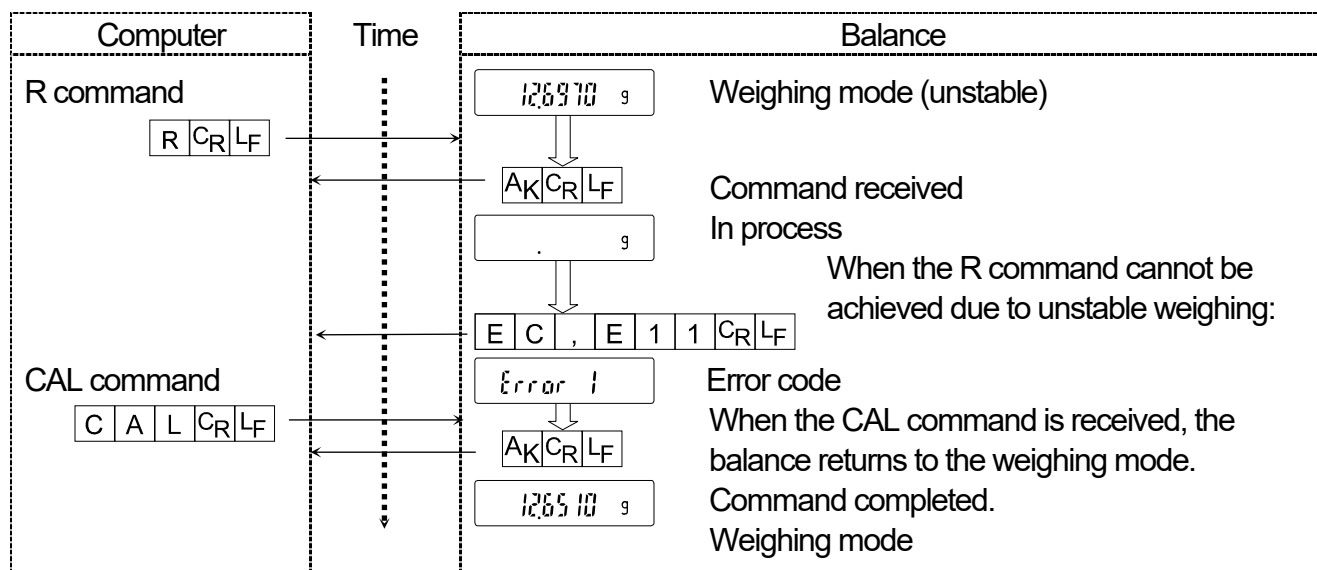


## Weighing with a tare

This example uses "Error" of "S if" so that the <AK> (06h) code is output.

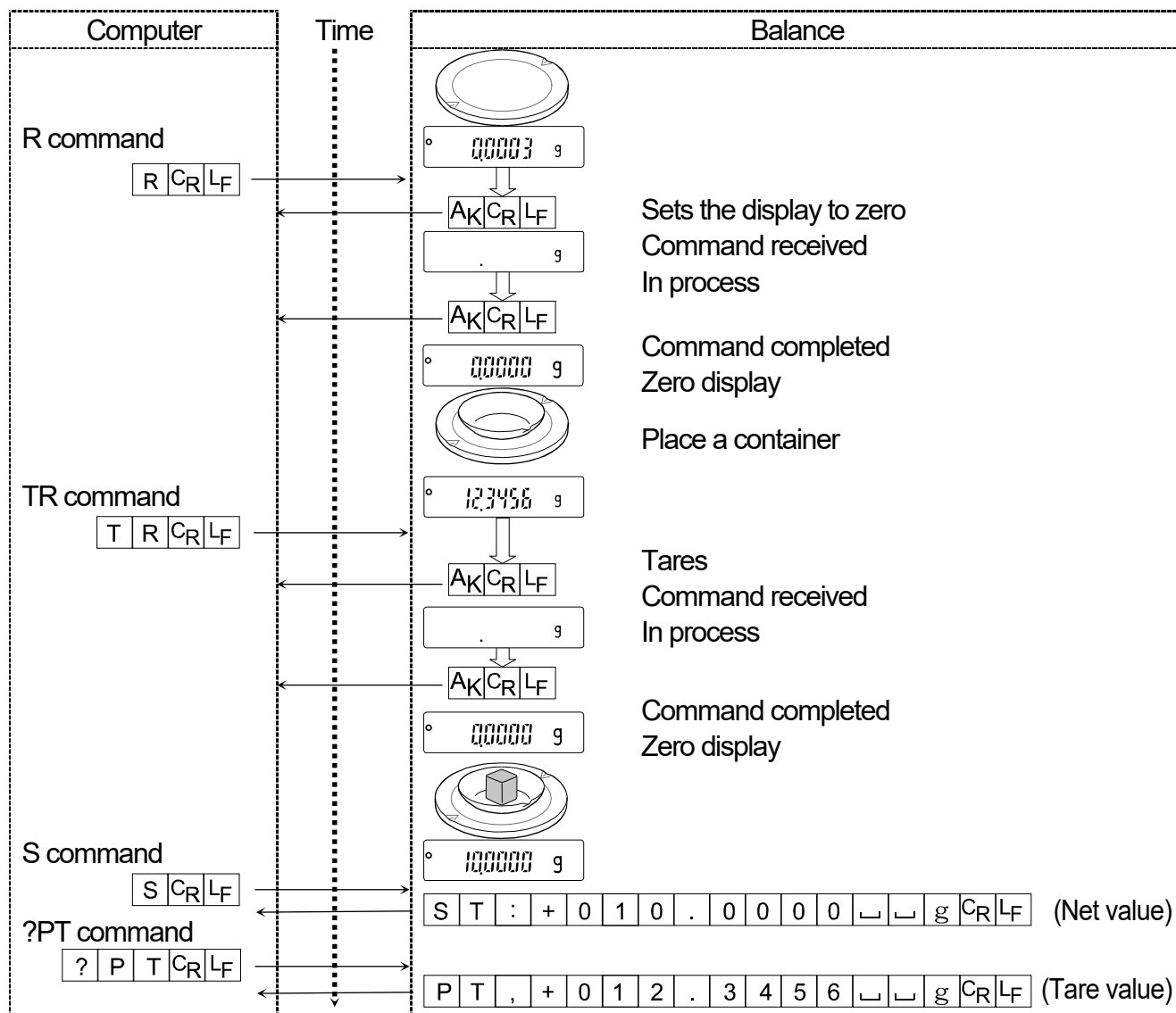

Setting a negative target value and filling with a sample until the display becomes zero

This example uses "ErCd on" of "S iF" so that the <AK> (06h) code is output.

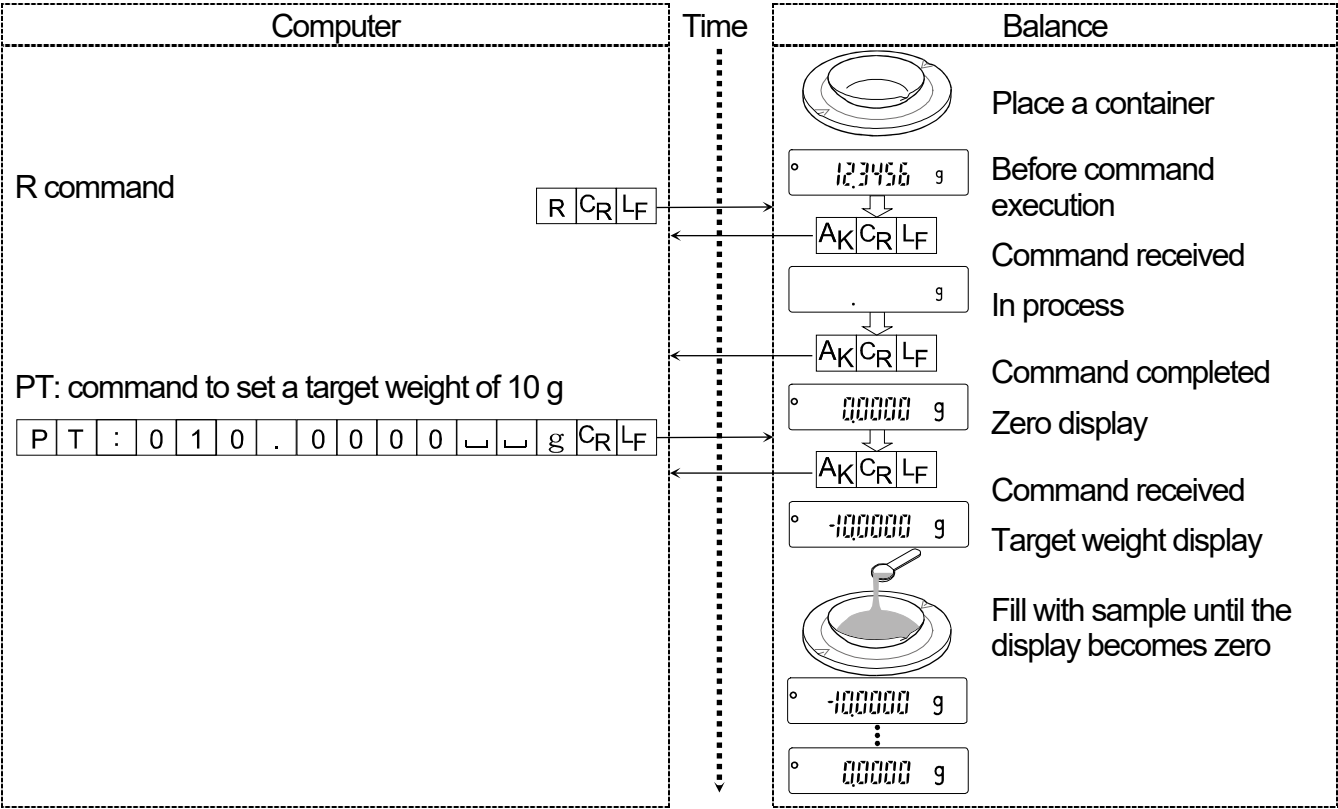

## 17. Key Lock Function

Sending specified commands to the balance will lock the keys on the balance to disable the key functions. This is useful for controlling the balance only by an external device such as a personal computer.

- Even if the keys are locked, key operations using commands are available. For details on commands to operate the keys, refer to “[16.3. Commands](#)”.
- The key lock state can be checked using the command to request the state of the balance keys.
- The key lock state is maintained until the balance receives the key unlocking command or the balance is turned off by removing the AC adapter.
- The key lock function is available for balances with software version 1.660 or later.

### 17.1. Locking All the Keys

Using the KL command, all the keys on the balance can be disabled.

| Command string                               | Description  |
|--|--|
| KL:***<br>(Enter 000 or 001 in place of ***) | KL:000    Unlocks all the keys.<br>KL:001    Locks all the keys  |
| ?KL  | Requests the state of the all key lock function.<br>KL, 000    All keys unlocked.<br>KL, 001    All keys locked. |



## 17.2. Locking the Specified Keys

Entering numerical values \*\*\*\*\* to the LK command will lock specified keys on the balance to disable their key functions. Set the numerical values \*\*\*\*\* by adding the decimal numbers converted from bits assigned to each key as shown below.

| Bits | Decimal | Keys    |
|------|---------|---------|
| 0    | 1       | ON:OFF  |
| 1    | 2       | CAL     |
| 2    | 4       | MODE    |
| 3    | 8       | RANGE   |
| 4    | 16      | PRINT   |
| 5    | 32      | RE-ZERO |
| 6    | 64      | SELECT  |
| 7    | 128     | ION     |

### Example 1: Locking all the keys excluding the **PRINT** key

- 1 Add the decimal numbers corresponding to the keys to be locked.  
 $1 \text{ (ON:OFF)} + 2 \text{ (CAL)} + 4 \text{ (MODE)} + 8 \text{ (RANGE)} + 32 \text{ (RE-ZERO)} + 64 \text{ (SELECT)} + 128 \text{ (ION)}$   
 $= 239$
- 2 Send the LK command to the balance with the sum that is obtained in Step 1. LK:00239

### Example 2: Unlocking all the keys


- 1 Send the LK command to the balance with 0, which indicates no keys to be locked. LK:00000

| Command string   | Description   |
|--|---|
| LK:*****<br>(Enter the numerical values 00000 to 00255 in place of *****.) | Locks the specified keys.<br>e.g.: LK:00239 Locks all the keys excluding the <b>PRINT</b> key.  |
| ?LK  | Requests the state of the specified key lock function.<br>e.g.1: LK, 00239 Specified keys locked.<br>e.g. 2: LK, 00000 All keys unlocked. |

## 18. Checking the Balance Software Version

The balance specifications vary depending on the balance software version. Follow the steps below to check the balance software version.

- 1 Unplug the AC adapter from the balance and plug it back in.
- 2 **P- \*\*\*\*\*** is displayed with the software version in place of \*\*\*\*\*.

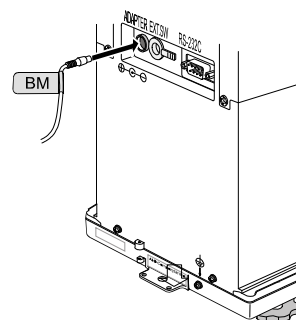




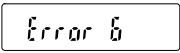
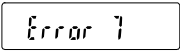
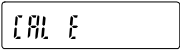
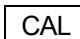
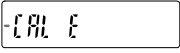

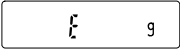
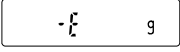

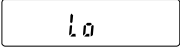
P- 1663











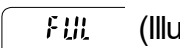
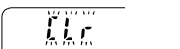
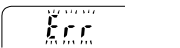
# 19. Maintenance
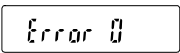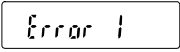
## 19.1. Cleaning and Handling the Balance

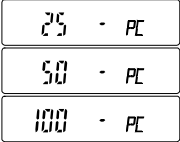- Clean the balance with a lint free cloth that is moistened with warm water and a mild detergent.
- Do not use organic solvents to clean the balance.
- Do not disassemble the balance. Contact the local A&D dealer if the balance needs service or repair.
- Use the original packing material for transportation.
- Consider "[3. Installing the Balance and Precautions](#)" when handling the balance.
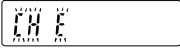
## 19.2. Error Codes

| Display   | Error code | Description   |
|---|------------|---|
|    | EC,E00     | <b>Internal error</b><br>If this error is displayed continuously, contact your local A&D dealer to repair the balance.  |
|    | EC,E11     | <b>Stability error</b><br>The balance cannot stabilize due to an environmental problem. Check around the pan. Prevent vibration, drafts, temperature changes, static electricity and magnetic fields, from influencing the balance.<br>Press the  key to return to the weighing mode. |
|  | EC,E16     | <b>Internal mass error</b><br>Applying the internal mass does not yield a change in the mass value as specified. Confirm that there is nothing on the pan and perform the weighing operation from the beginning again.  |
|  | EC,E17     | <b>Internal mass error</b><br>The internal mass application mechanism does not function properly. Perform the weighing operation from the beginning again.  |
|  | EC,E20     | <b>Sensitivity adjustment weight error</b><br>The sensitivity adjustment weight is too heavy. Confirm the sensitivity adjustment weight value. Press the  key to return to the weighing mode.   |
|  | EC,E21     | <b>Sensitivity adjustment weight error</b><br>The sensitivity adjustment weight is too light. Confirm the sensitivity adjustment weight value. Press the  key to return to the weighing mode.   |
|  |            | <b>Overload error</b><br>A sample beyond the balance weighing capacity has been placed on the pan. Remove the sample from the pan.  |
|  |            | <b>Weighing pan error</b><br>The mass value is too light. Confirm that the weighing pan and pan support are properly installed. Press the  key two times to return to the weighing mode. If the error still persists, adjust the balance.   |
|  |            | <b>Sample mass error</b><br>The balance cannot store the sample for the counting mode or for the percent mode because it is too light. Use a larger sample.   |
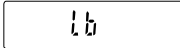
| Display  | Error code | Description   |
|--|------------|---|
|                     |            | <b>Unit mass error</b><br>The sample unit mass for the counting mode is too light. Storing and using it for counting will cause a counting error.<br>Add samples to reach the specified number and press the  key.<br>Pressing the  key without adding samples will shift the balance to the counting mode. But, for accurate counting, be sure to add samples. |
|                     |            | <b>Automatic response adjustment zero error</b><br>The automatic response adjustment cannot be performed because there is something on the pan. Clear the pan. Press the  key to return to the weighing mode.  |
| <br>(Check NG)      |            | <b>Automatic response adjustment unstable error</b><br>The automatic response adjustment cannot be performed because the mass value is unstable.<br>Check the ambient conditions such as breeze, vibration and magnetic fields, also check the weighing pan.<br>Press the  key to return to the weighing mode.   |
|                     |            | <b>Clock battery error</b><br>The clock backup battery has been depleted.<br>Press any key and set the time and date. The clock and calendar function works normally as long as the AC adapter is connected to the balance. If this error appears frequently, contact the local A&D dealer.   |
|                   |            | <b>Low battery error</b><br>Voltage of the power supply drops.<br>Confirm whether the correct AC adapter is used.   |
| <br>(Blinking)    |            | <b>Memory full</b><br>The amount of weighing data in memory has reached the maximum capacity. Delete data in memory to store new data. For details, refer to " <a href="#">12. Data Memory</a> ".   |
| <br>(Illuminated) |            | <b>Memory full</b><br>The amount of sensitivity adjustment or calibration test data in memory has reached the maximum capacity (50 sets).<br>The data in memory will be deleted automatically to store new data. For details, refer to " <a href="#">12. Data Memory</a> ".   |
|                   |            | <b>Memory type error</b><br>Type of memory set in the function table and type of data stored are different. For details, refer to " <a href="#">12. Data Memory</a> ".  |
|                   |            | <b>Memory data error</b><br>A fatal error occurs in the stored data. To use the memory function, delete all the old data and resolve the error. For details, refer to " <a href="#">12.1 Notes on Using Data Memory</a> ".  |
|  | EC,E00     | <b>Communications error</b><br>A protocol error occurred in communications.<br>Confirm the format, baud rate and parity.  |
|  | EC,E01     | <b>Undefined command error</b><br>An undefined command was received.<br>Confirm the command.  |


| Display          | Error code | Description   |
|------------------|------------|---|
|                  | EC,E02     | <b>Not ready</b><br>A received command cannot be processed.<br>Examples: <ul style="list-style-type: none"> <li>□ The balance received a "Q" command, but not in the weighing mode.</li> <li>□ The balance received a "Q" command while processing a RE-ZERO command.</li> </ul> Adjust the delay time to transmit a command. |
|                  | EC,E03     | <b>Timeout error</b><br>When the timeout parameter is set to "t -UP on", the balance did not receive the next character of a command within the time limit of one second. Confirm the communication.  |
|                  | EC,E04     | <b>Excess characters error</b><br>The balance received excessive characters in a command. Confirm the command.  |
|                  | EC,E06     | <b>Format error</b><br>A command includes incorrect data.<br>Example: <ul style="list-style-type: none"> <li>□ The data is numerically incorrect.</li> </ul> Confirm the command.   |
|                  | EC,E07     | <b>Parameter setting error</b><br>The received data exceeds the range that the balance can accept. Confirm the parameter range of the command.  |
| Other error code |            | If an error described above cannot be resolved or other errors are displayed, contact the local A&D dealer.   |

### 19.3. Other Display



When this indicator blinks, automatic sensitivity adjustment is required. The indicator blinks when the balance detects changes in ambient temperature. If the balance is not used for several minutes with this indicator blinking, the balance performs automatic sensitivity adjustment. The blinking duration depends on the operating environment.

**Note** The balance can be used while this indicator is blinking. We recommend that you perform automatic sensitivity adjustment for precision weighing.

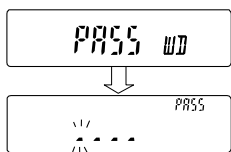

This is the display to prompt you to enter a password. When the password function is used, the correct password must be entered for the balance to enter the weighing mode. For details, refer to "[15. Password Function](#)"

## 19.4. Checking the Balance Performance and Environment

The balance is a precision instrument. When the operating environment or the operating method is inadequate, correct weighing cannot be performed. Place a sample on the pan and remove it, and repeat this several times. If the balance seems to have a problem with repeatability or to perform improperly, check as described below. If improper performance persists after checking, contact the local A&D dealer for repair.

### Checking that the operating environment or weighing method is proper

#### Operating environment

- Is the weighing table solid enough?
- Is the balance level?
- Is the operating environment free from vibration and drafts?
- Is there a strong electrical or magnetic noise source such as a motor near the balance?

#### Weighing method

- Does the weighing pan edge touch anything? Is the weighing pan assembly installed correctly?
- Is the **RE-ZERO** key pressed before placing a sample on the weighing pan?
- Is the sample placed in the center of the weighing pan?
- Is the fine range ring installed for weighing with a readability of 0.01 mg for the **BM-252**?
- Is the fine range ring and small fine range ring installed for weighing with a readability of 0.001 mg for the **BM-5**, **BM-5D**, **BM-20** and **BM-22**?

Is the tabletop breeze break (M) installed for weighing with a readability of 0.001 mg for the **BM-20** and **BM-22**?

- Has the balance been adjusted using the internal mass?
- Has the balance been warmed up for one hour before weighing?

#### Sample and container

- Has the sample absorbed or lost moisture due to the ambient conditions such as temperature and humidity?
- Has the temperature of the container been allowed to equalize to the ambient temperature?
- Is the sample charged with static electricity? Is the built-in ionizer (page 32) used to discharge the sample charged with static electricity.
- Is the sample a magnetic material such as iron? Cautions are required when weighing magnetic materials.

### Checking that the balance performs properly

- Check the balance repeatability using an external weight. Be sure to place the weight in the center of the weighing pan.
- Check the balance repeatability, linearity and calibration test using an external weight with a known value.
- Check whether the sensitivity adjustment is performed correctly. If an error occurs and the sensitivity adjustment cannot be performed quickly, the balance may not measure correctly.

## 19.5. Asking for Repair

If the balance needs service or repair, contact your local A&D dealer.

The balance is a precision instrument. Use much care when handling the balance and observe the following when transporting the balance.

- Use the original packing material for transportation.
- Remove the weighing pan, pan support, rings and dust plate from the balance.

## 20. Specifications

|  |                             | BM-5  | BM-5D              | BM-20                 | BM-22              |
|--|-----------------------------|---|--------------------|-----------------------|--------------------|
| Weighing capacity  |                             | 5.2 g   | 5.2 g              | 22 g                  | 22 g               |
|  |                             |   | 2.1 g              |                       | 5.1 g              |
| Maximum display  |                             | 5.200084 g  | 5.20008 g          | 22.000084 g           | 22.00008 g         |
|  |                             |   | *1 2.100009 g      |                       | *1 5.100009 g      |
| Readability  |                             | 0.001 mg  | 0.01 mg            | 0.001 mg              | 0.01 mg            |
|  |                             |   | 0.001 mg           |                       | 0.001 mg           |
| Repeatability<br>(Standard deviation)                            |                             | *2 0.0012 mg  | 0.01 mg            | *2 0.0025 mg          | 0.01 mg            |
|  |                             |   | *2 0.004 mg        |                       | *2 0.004 mg        |
| Linearity  |                             | ±0.010 mg   | ±0.02 mg           | ±0.010 mg             | ±0.02 mg           |
|  |                             |   | ±0.010 mg          |                       | ±0.010 mg          |
| Stabilization time<br>(Typical at <b>FAST</b> )                  |                             | Approx.<br>10 seconds   | Approx. 8 seconds  | Approx.<br>10 seconds | Approx. 8 seconds  |
|  |                             |   | Approx. 10 seconds |                       | Approx. 10 seconds |
| Sensitivity drift  |                             | ±2 ppm/°C (10 °C to 30 °C, Automatic sensitivity adjustment: OFF)   |                    |                       |                    |
| Operating environment  |                             | 5 °C to 40 °C, 85 % RH or less (No condensation)  |                    |                       |                    |
| Display refresh rate   |                             | 5 times/second or 10 times/second   |                    |                       |                    |
| Environmental sensors  |                             | Resolution: Temperature ±1.5 °C, Humidity ±10 %, Barometric pressure ±10 hPa<br>Applicable range: 5 °C to 40 °C   |                    |                       |                    |
| Counting mode  | Minimum unit mass           | 0.1 mg  |                    |                       |                    |
|  | Number of samples           | 10, 25, 50 or 100 pieces  |                    |                       |                    |
| Percent mode   | Minimum 100% reference mass | 10 mg   |                    |                       |                    |
|  | Readability                 | 0.01 %, 0.1 %, 1 % (Depends on the reference mass stored.)  |                    |                       |                    |
| Interface  |                             | USB and RS-232C   |                    |                       |                    |
| Applicable weights for calibration test / sensitivity adjustment |                             | 5 g   | 5 g                | 20 g                  | 20 g               |
|  |                             | 2 g   | 2 g                | 10 g                  | 10 g               |
|  |                             | 1 g   | 1 g                | 5 g                   | 5 g                |
|  |                             |   |                    | 2 g                   | 2 g                |
|  |                             |   |                    | 1 g                   | 1 g                |
| Size of loading area for weighing mm                             |                             | Weighing pan: φ25<br>Filter weighing: φ50, φ95<br>Micro tube holder: φ12  |                    |                       |                    |
| External dimensions  |                             | 259(W) x 466(D) x 326(H) mm   |                    |                       |                    |
| Power supply & AC adapter type                                   |                             | Power consumption: Approx. 30 VA (supplied to the AC adapter )<br>Confirm that the adapter type is correct for the local voltage and power receptacle type. |                    |                       |                    |
| Mass   |                             | Approx. 10 kg   |                    |                       |                    |

- The operating environment with no excessive changes in ambient temperature and humidity, no vibration and drafts, no influence by magnetic fields and static electricity.
- The internal mass may change due to corrosion or other damage caused by the operating environment, or due to aging. Check the internal mass periodically and correct the internal mass value if necessary.

\*1: Even when a tare weight beyond the precision range is placed, the weighing sample can be weighed in the precision range if using the smart range.


\*2: Repeatability using a 1 gram weight within the weighing capacity (when the tabletop breeze break (M) is used).

|  |                             | BM-252  | BM-200                 | BM-300                          | BM-500                                   |
|--|-----------------------------|---|------------------------|---------------------------------|--|
| Weighing capacity  |                             | 250 g   | 220 g                  | 320 g                           | 520 g                                    |
| Maximum display  |                             | 250.00084 g   | 220.0084 g             | 320.0084 g                      | 520.0084 g                               |
| Readability  |                             | 0.01 mg   | 0.1 mg                 |                                 |  |
| Repeatability<br>(Standard deviation) *3                         |                             | 0.03 mg   | 0.1 mg                 | 0.2 mg                          |  |
| Linearity  |                             | ±0.10 mg  | ±0.2 mg                | ±0.3 mg                         | ±0.5 mg                                  |
| Stabilization time<br>(Typical at <span>FAST</span> )            |                             | Approx.<br>8 seconds  | Approx. 3.5 seconds    |                                 |  |
| Sensitivity drift  |                             | ±2 ppm/°C (10 °C to 30 °C, Automatic sensitivity adjustment: OFF)   |                        |                                 |  |
| Operating environment  |                             | 5 °C to 40 °C, 85 % RH or less (No condensation)  |                        |                                 |  |
| Display refresh rate   |                             | 5 times/second or 10 times/second   |                        |                                 |  |
| Environmental sensors  |                             | Resolution: Temperature ±1.5 °C, Humidity ±10 %, Barometric pressure ±10 hPa<br>Applicable range: 5 °C to 40 °C   |                        |                                 |  |
| Counting mode  | Minimum unit mass           | 0.1 mg  |                        |                                 |  |
|  | Number of samples           | 10, 25, 50 or 100 pieces  |                        |                                 |  |
| Percent mode   | Minimum 100% reference mass | 10 mg   |                        |                                 |  |
|  | Readability                 | 0.01 %, 0.1 %, 1 % (Depends on the reference mass stored.)  |                        |                                 |  |
| Interface  |                             | USB and RS-232C   |                        |                                 |  |
| Applicable weights for calibration test / sensitivity adjustment |                             | 200 g<br>100 g<br>50 g<br>20 g<br>10 g  | 200 g<br>100 g<br>50 g | 300 g<br>200 g<br>100 g<br>50 g | 500 g<br>300 g<br>200 g<br>100 g<br>50 g |
| Size of loading area for weighing mm                             |                             | Weighing pan: φ90<br>Micro tube holder: φ12 (Option)  |                        |                                 |  |
| External dimensions  |                             | 259(W) x 466(D) x 326(H) mm   |                        |                                 |  |
| Power supply & AC adapter type                                   |                             | Power consumption: Approx. 30 VA (supplied to the AC adapter )<br>Confirm that the adapter type is correct for the local voltage and power receptacle type. |                        |                                 |  |
| Mass   |                             | Approx. 10 kg   |                        |                                 |  |

- The operating environment with no excessive changes in ambient temperature and humidity, no vibration and drafts, no influence by magnetic fields and static electricity.
- The internal mass may change due to corrosion or other damage caused by the operating environment, or due to aging. Check the internal mass periodically and correct the internal mass value if necessary.

\*3: Repeatability using a 100 gram weight within the weighing capacity.


## 20.1. External Dimensions

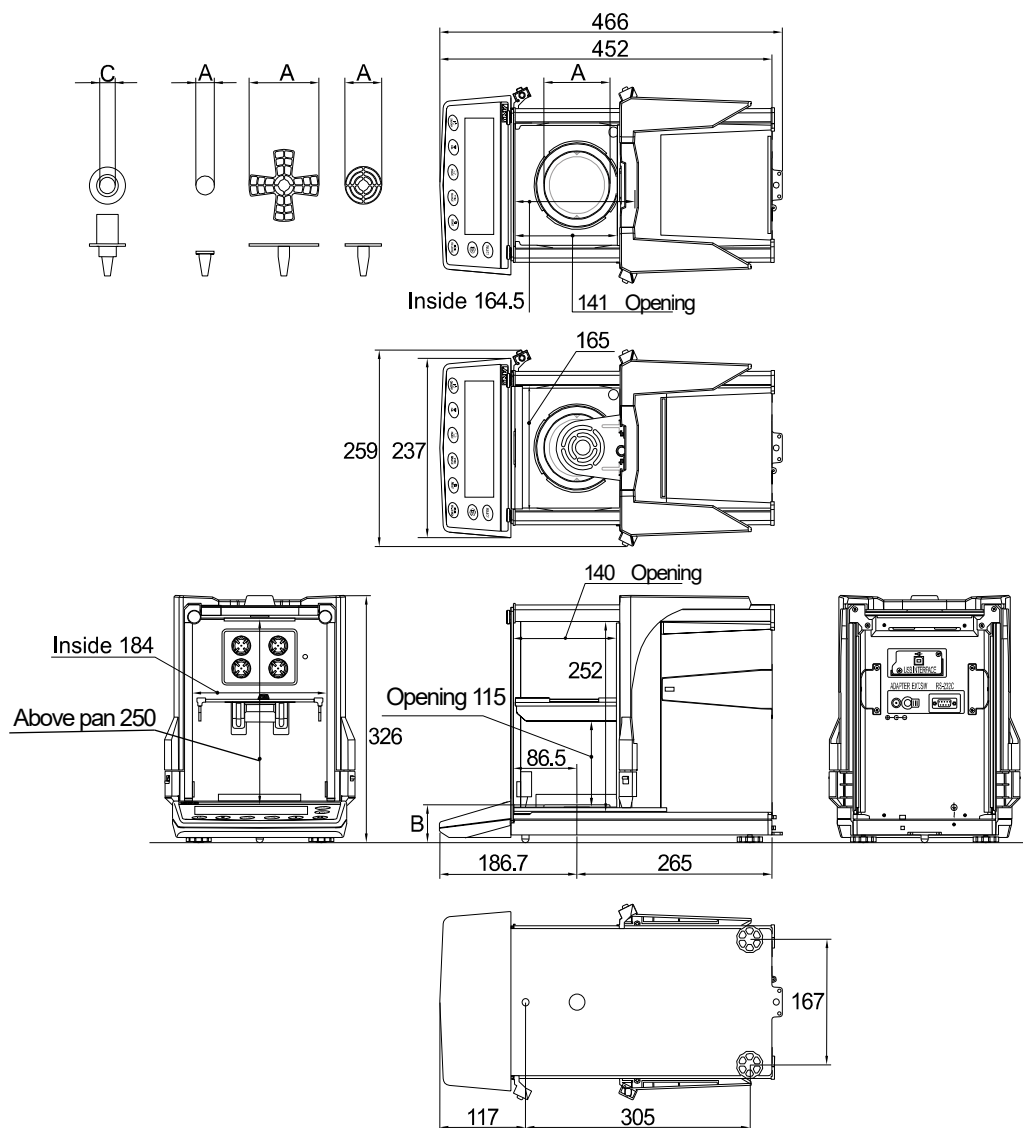

Unit: mm

|                                      | <b>A:</b> Diameter of the weighing pan                                 | <b>B:</b> Height of the weighing pan | <b>C:</b> The inside diameter of the test tube holder |
|--------------------------------------|--|--------------------------------------|---|
| <b>BM-5, BM-5D, BM-20, BM-22</b>     | Weighing pan: $\phi 25$<br>Weighing pan for filter: $\phi 50, \phi 95$ | Weighing pan: $\phi 90$              | $\phi 12$   |
| <b>BM-252, BM-200 BM-300, BM-500</b> | Weighing pan: $\phi 90$  | 50                                   | $\phi 12$ (Option)                                    |

Tabletop breeze break (M)


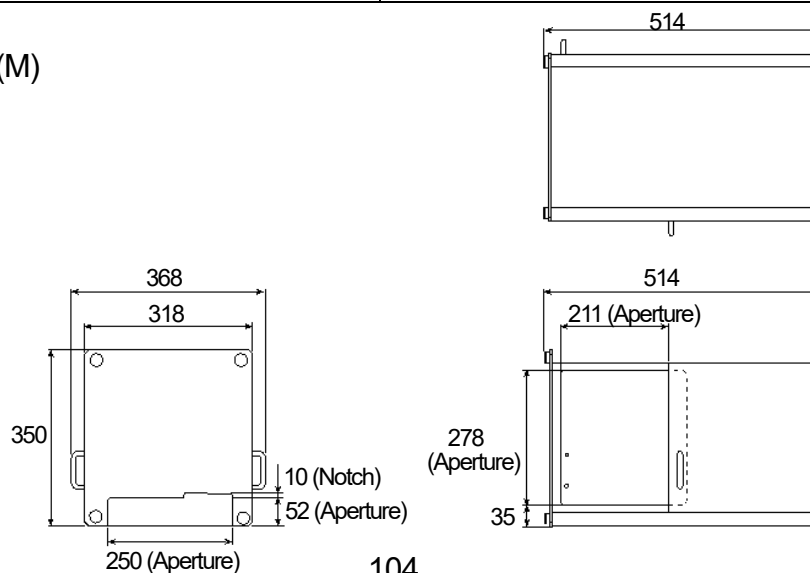




## 20.2. Options and Peripheral Equipment

### Ethernet Interface BM-08


- The interface to connect the balance to a LAN.
- Accessory: "WinCT-Plus" data communication software.
  - The software can acquire data from multiple balances connected to a LAN.
  - The software can control these balances with commands.
  - The software can acquire data transmitted from balances.  
Example: When pressing the **PRINT** key of the balance, data is output and is acquired by the computer.
  - The stored data can be used with Microsoft EXCEL (if installed).

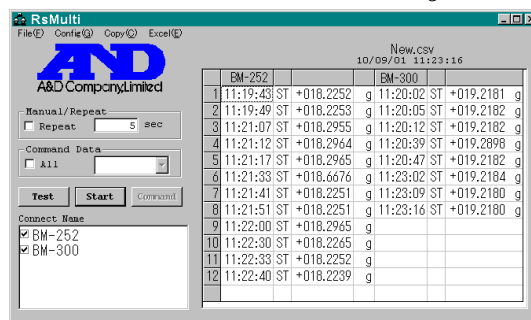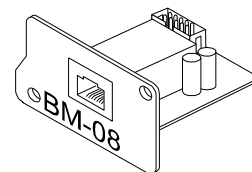

### Pipette Accuracy Tester BM-014


**Note** This option cannot be used for the BM-5 and BM-5D.

- This option can be used to inspect the volume dispensed from the pipette.
- The accessory evaporation trap is used to prevent the evaporation of the test liquid inside it.
- The WinCT-Pipette software converts the test liquid mass into a volumetric value.

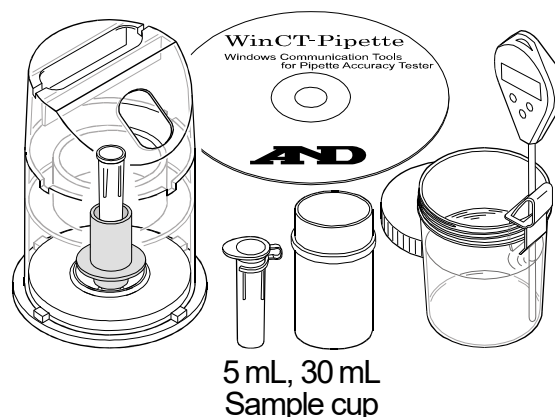

### Electrode Unit of the Ionizer AX-BM-NEEDLESET (4 pieces)

- Use these electrode units for replacement parts of the ionizer built into the balance.
- Replace all four units with new ones at the same time.
- Refer to "6.5. Ionizer (Neutralization Device)" for the replacement.

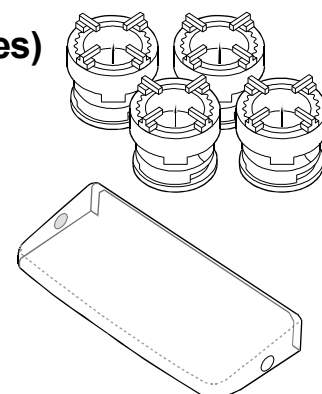

### Display Cover AX-BM-031 (5 pieces)


- This option is the same as the accessory display cover.

### Micro Tube Holder AX-BM-032 (for the BM-5/5D/20/22, standard accessory) AX-BM-034 (for the BM-252/200/300/500)

- This option is used when weighing with the general micro tube or test tube under  $\phi 12$  mm.


**Note** When using the micro tube holder for the BM-5 and BM-5D:

The micro tube holder is approximately 2 g heavier than the standard pan (weighing pan and pan support). When a micro tube holder is used for the BM-5 and BM-5D, the available weighing range will be up to approximately 3 g including the tare such as a micro tube and test tube.

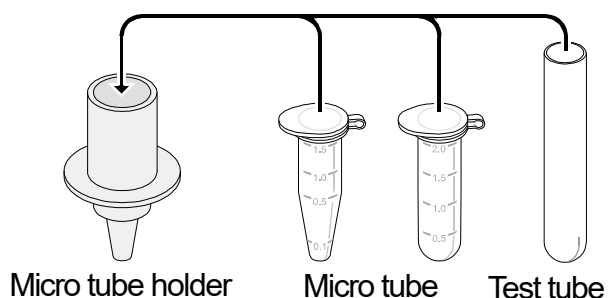


## Body Cover AX-BM-033

- Cover during storage to prevent dust and dirt contamination.
- The cover can be used for the balance with the AC adapter and other cables connected.
- When removing the cover, take care so as not to get its edge caught on the balance.
- Specification
  - Material: Nylon
  - Color : Light gray

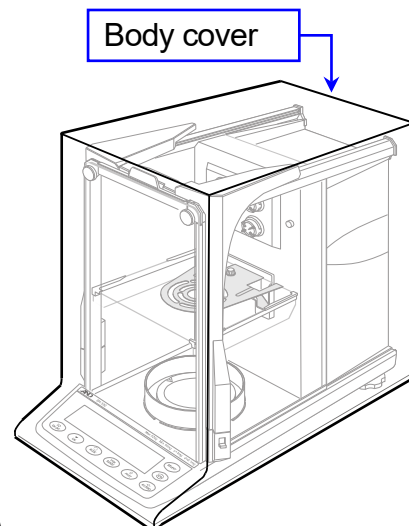

## Slide Breeze Break AX-BM-035 (for the BM-5/5D/20/22)

- This option is attached to the weighing chamber to reduce the effects of the surrounding environment so as to improve weighing accuracy.

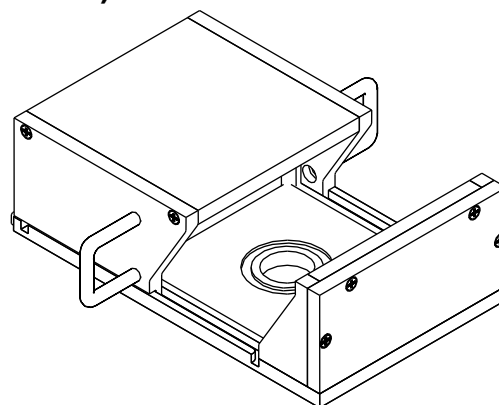

## Aluminum Pans

- This option is the same as the accessory aluminum pans used for weighing a very small amount of sample.

| Name and description              |     |         |            |  | Order code     | View |
|-----------------------------------|-----|---------|------------|--|----------------|------|
| Round aluminum analytical pan (L) | φ15 | 0.8 mL  | 100 pieces |  | AX-ROUND-PAN-L |      |
| Round aluminum analytical pan (M) | φ12 | 0.3 mL  | 100 pieces |  | AX-ROUND-PAN-M |      |
| Round aluminum analytical pan (S) | φ8  | 0.05 mL | 100 pieces |  | AX-ROUND-PAN-S |      |


## AD-1653 Density Determination Kit

**Note** This option cannot be used for the BM-5, BM-5D, BM-20 and BM-22.

$$\frac{\text{Weight in air}}{\text{Weight in air} - \text{Weight in liquid}} \times \text{liquid density} = \text{sample density}$$

Example  $\frac{10.0000 \text{ g}}{10.0000 \text{ g} - 9.5334 \text{ g}} \times 0.9970 \text{ g/cm}^3 = 21.4 \text{ g/cm}^3$

| Temperature | Liquid density            |
|-------------|---------------------------|
| 0 °C        | 0.99984 g/cm <sup>3</sup> |
| 10 °C       | 0.99970 g/cm <sup>3</sup> |
| 20 °C       | 0.99820 g/cm <sup>3</sup> |
| 30 °C       | 0.99565 g/cm <sup>3</sup> |

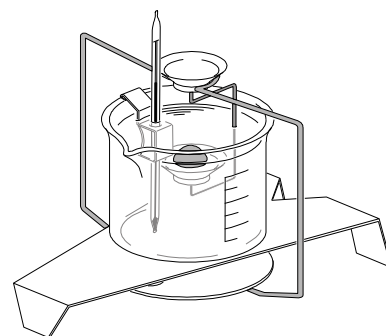

## Option

### **AD-8129TH: Thermal printer**

- Small direct thermal printer that connects to the balance via the RS-232C interface.
- Statistical calculation mode, calendar/clock function, interval printing mode, chart printing mode, dump printing mode

### **AD-1691: Balance environment analyzer**

- During the routine inspection, repeatability and the minimum weighing value can be easily checked.
- Also, this calculates measurement uncertainty and supports environmental evaluation of the balance.
- This analyzer can be easily carried to the balance site, so more than one balance can be controlled using this analyzer.

### **AD-1687: Weighing environment logger**

- A data logger equipped with 4 sensors for temperature, humidity, barometric pressure and vibration that can measure and store environmental data. When connected to the RS-232C interface of the balance, the AD-1687 can store environmental data along with weighing data. Therefore, it is possible to store data in an environment where a computer cannot be used.
- The stored data can be read to a personal computer using USB. As the AD-1687 is recognized as USB memory, special software is not required to read the data.

### **AD-1688: Data logger**

- When connected to the RS-232C interface of the balance, the AD-1688 can store the data in an environment where a personal computer cannot be used.

### **AD-8526: Ethernet converter**

- This option can be used to connect the RS-232C interface of the balance to the Ethernet (LAN) port of a computer. This allows management of the balance weighing data with a computer connected to a network.

### **AD-8527: Quick USB adapter**

- This option transmits the weighing data directly to software applications such as Excel and Word.

### **AD-8920A: Remote display**

- Connected to the balance using the RS-232C interface.

### **AD-8922A: Remote controller**

- This option can be connected to the balance using the RS-232C interface and can control the balance remotely. Various options such as comparator output or analog output are available.

### **AX-USB-9: USB converter**

- Adds a COM port to a personal computer.
- Enables bi-directional communication between the personal computer and the balance when a USB driver is installed.
- Can use serial communication software such as WinCT on a personal computer without COM ports.
- An RS-232C cable is provided to connect the USB converter to the balance.

### **AD-1671: Anti-vibration table**

- Approximately 27 kg, Artificial stone (Terrazzo).
- Use this table when unstable weighing is caused by vibration coming from the floor.
- Use the AD-8922A remote controller to avoid a weighing error by a slight table tilt that key operation may make.

### **AD-1672/AD-1672A Tabletop breeze break (Large)**

- Large sized table breeze break
- Protects the balance from wind from sources such as air conditioning or people passing by, thereby reducing balance weighing errors.
- The transparent panel assembly consists of antistatic plastic material that protects the balance from static electricity.
- Samples can be inserted or removed through the opening on the front.
- AD-1672: 680(W) x 600(D) x 720(H) mm, AD-1672A: 680(W) x 584(D) x 720(H) mm

**AD-1676: Tabletop breeze break (Medium)( for the BM-5/5D/20/22, standard accessory)**

- A tabletop breeze break of a size best suited for the BM series balances.
- Protects the balance from wind from sources such as air conditioning or people passing by, thereby reducing balance weighing errors.
- The transparent panel assembly consists of antistatic plastic material that protects the balance from static electricity.
- Sliding panels on the right and left side enable samples to be inserted or removed from each side.
- 368(W) x 514(D) x 350(H) mm (Including the handle)

**AD-1682: Rechargeable Battery**

- This option allows use of the balance in a place where AC power is not available.

**AD-1683A: Ionizer**

- This ionizer prevents weighing error caused by static electricity on the measured sample.
- DC operation generates a large amount of ions with no wind, making this device suitable for precise weighing of powders and similar materials.
- Can be operated by infrared sensor to eliminate static electricity without making contact. (For details, See “AD-1683A Instruction Manual”)

**AD-1684A: Electrostatic field meter**

- This option measures the amount of the static charge on the sample, tare or peripheral equipment and displays the result.
- If those are found to be charged, discharge them using the AD-1683 DC static eliminator.

**AD-1689: Tweezers for sensitivity adjustment weight**

- A pair of tweezers ideally suited for holding sensitivity adjustment weights of 1 g to 500 g.

**AX-TWEEZERS-25: Tweezers for balance**

- A 250 mm length makes the tweezers suitable for a balance with a large breeze break.
- Used for plate weight (1 mg to 500 mg) / cylindrical weight (1 g to 500 g)
- Material: SUS316 / Front end: PC resin.

**AD-8541-SCALE: Bluetooth converter (for connecting a weighing device)\***

- This option connects a smartphone, a tablet or a PC to the balance wirelessly (via Bluetooth®) up to 10 m.
- The “WinCT-WeiV” dedicated application software for a smartphone and a tablet is available.
- To connect a PC, the AD-8541-PC mentioned below needs to be connected to the PC side.

**AD-8541-PC: Bluetooth dongle (for connecting a PC)\***

- This option connects a PC to the balance wirelessly (via Bluetooth®) up to 10 m.
- Used in a set with the aforementioned AD-8541-SCALE.

**\*Note on AD-8541-SCALE and AD-8541-PC**

The two products may require certification for compliance with local laws pertaining to radio broadcast and equipment for wireless communication. Please refer to the A&D website for countries where the products are already certified for use.

**AX-SW137-PRINT: Foot switch print function with a plug**

- The foot switch has the same function as the PRINT key.


**AX-SW137-REZERO: Foot switch rezero function with a plug**

- The foot switch has the same function as the RE-ZERO key.


### **AX-HOLDER-SET: Sample cup holder**

**Note** This option cannot be used for the BM-5, BM-5D, BM-20 and BM-22.

- A set of holders to secure a sample cup such as a test tube, centrifuge tube or micro tube.

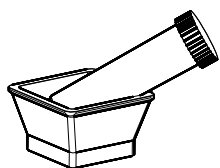


Centrifuge tube  
secured slantingly

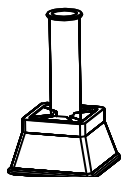

Sample cup holder used  
as test tube stand



Sample cup holder used  
as holder for micro tube

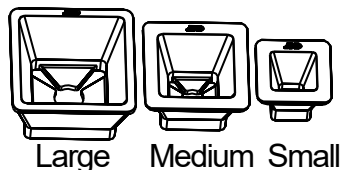

Large

Medium

Small

### **AX-CLEANING-SET: Cleaning set for a balance**

- Used to clean the weighing pan and surroundings.
- The thin and soft fiber material is suitable for cleaning powder material.
- As the fiber material is animal fiber, it does not easily generate static electricity.

\*No function to eliminate static electricity

### **AX-TRAY-10ML: Weighing board for an analytical balance**


- Antistatic treatment prevents static electricity.
- Volume : 10 mL / Mass within 0.49 g / Capacity : 250 sheets

### **AX-SPATULA: Micro spatula for an analytical balance**

- Especially suitable for treating fine powder.
- Length: 182 mm (Mass: 4.3 g / Number included: 5 pcs)

## 21. Terms

|                               |   |
|-------------------------------|---|
| Sensitivity adjustment        | Adjustment of the balance so that it can weigh accurately.  |
| Sensitivity adjustment weight | A weight used for sensitivity adjustment.   |
| Data number                   | Numbers assigned sequentially when weighing data or unit weight is stored.  |
| Digit                         | The minimum weighing value available. Used for the balance, one digit is the smallest mass that can be displayed.   |
| Environment                   | Ambient conditions such as vibration, drafts, temperature, static electricity and magnetic fields which affect the weighing operation.  |
| External weight               | A weight that you have.   |
| GLP                           | Good Laboratory Practice.   |
| GMP                           | Good Manufacturing Practice.  |
| Internal mass                 | Built-in sensitivity adjustment weight.   |
| ISO                           | International Organization for Standardization.   |
| Micro tube                    | 2 mL and 1.5 mL micro tubes in the market can be weighed with the micro tube holder.  |
| Repeatability                 | Variation in measured values obtained when the same mass is placed and removed repetitively. Usually expressed as a standard deviation.<br>Example: "Standard deviation = 1 digit " means that a weighing error may occur within $\pm 1$ digit in the frequency of approx. 68% when weighing the same mass repeatedly.  |
| Re-zero                       | To set the display to zero.   |
| Sensitivity drift             | An affect that a change in temperature causes to the weighing data. Expressed as a temperature coefficient.<br>Example: Temperature coefficient = 2 ppm/°C, If a load is 10 g and the temperature changes by 10°C, the value displayed changes by the following value.<br>$0.000\,002\text{ ppm/}^{\circ}\text{C} \times 10\text{ }^{\circ}\text{C} \times 10\text{ g} = 0.0002\text{ g}$ |
| Stabilization time            | Time required after a sample being placed, until the stabilization indicator illuminates and the weighing data is displayed.  |
| Tare                          | To cancel the mass value of a container which is not to be included in the weighing data. Normally, refers to an operation of placing a container and setting the display to zero.  |
| Target weight                 | An external weight used for calibration test.   |
| Zero point                    | A weighing reference point. Usually refers to the value displayed when nothing is on the weighing pan   |

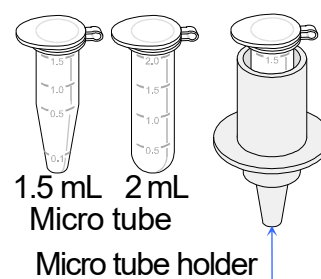

THIS PAGE INTENTIONALLY LEFT BLANK.



### **A&D Company, Limited**

3-23-14 Higashi-Ikebukuro, Toshima-ku, Tokyo 170-0013, JAPAN

Telephone: [81] (3) 5391-6132

Fax: [81] (3) 5391-1566

### **A&D ENGINEERING, INC.**

Headquarters Office: 4622 Runway Boulevard, Ann Arbor, Michigan 48108, U.S.A.

Sales Office: 47747 Warm Springs Boulevard, Fremont, California 94539, U.S.A.

Tel: [1] (800) 726-3364

Weighing Support: [1] (888) 726-5931

Inspection Support: [1] (855) 332-8815

### **A&D INSTRUMENTS LIMITED**

Unit 24/26 Blacklands Way, Abingdon Business Park, Abingdon, Oxfordshire OX14 1DY United Kingdom

Telephone: [44] (1235) 550420

Fax: [44] (1235) 550485

### **A&D AUSTRALASIA PTY LTD**

32 Dew Street, Thebarton, South Australia 5031, AUSTRALIA

Telephone: [61] (8) 8301-8100

Fax: [61] (8) 8352-7409

### **A&D KOREA Limited**

한국에이.엔.디(주)

서울특별시 영등포구 국제금융로6길33 (여의도동) 맨하탄빌딩 817 우편 번호 07331

( 817, Manhattan Bldg., 33. Gukjegeumyung-ro 6-gil, Yeongdeungpo-gu, Seoul, 07331 Korea )

전화: [82] (2) 780-4101

팩스: [82] (2) 782-4264

### **ООО A&D RUS**

ООО "Эй энд Ди Рус"

Почтовый адрес: 121357, Российская Федерация, г. Москва, ул. Вереysкая, дом 17

Юридический адрес: 117545, Российская Федерация, г. Москва, ул. Дорожная, д.3, корп.6, комн. 86

( 121357, Russian Federation, Moscow, Vereyskaya Street 17 )

тел.: [7] (495) 937-33-44

факс: [7] (495) 937-55-66

### **A&D Instruments India Private Limited**

ऐ&डी इन्स्ट्रुमेंट्स इण्डिया प्रा० लिमिटेड

D-48, उद्योग विहार , फेस -5, गुडगांव - 122016, हरियाणा , भारत

( D-48, Udyog Vihar, Phase-V, Gurgaon - 122016, Haryana, India )

फोन : [91] (124) 4715555

फैक्स : [91] (124) 4715599

### **A&D SCIENTECH TAIWAN LIMITED.**

艾安得股份有限公司

台灣台北市中山區南京東路2段206號11樓之2

( 11F-2, No.206, Sec.2, Nanjing E.Rd., Zhongshan Dist, Taipei City 10489, Taiwan, R.O.C. )

Tel : [886](02) 2322-4722

Fax : [886](02) 2392-1794

### **A&D INSTRUMENTS (THAILAND) LIMITED**

บริษัท เอ แอนด์ ดี อินสตรูमेंท์ (ไทยแลนด์) จำกัด

168/16 หมู่ที่ 1 ตำบลรังสิต อำเภอธัญบุรี จังหวัดปทุมธานี 12110 ประเทศไทย

( 168/16 Moo 1, Rangsit, Thanyaburi, Pathumthani 12110 Thailand )

Tel : [66] 20038911