


INSTRUCTION MANUAL

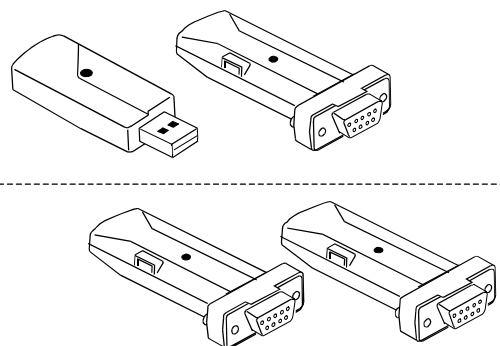

© 2017 A&D Company Ltd. All rights reserved.
No part of this publication may be reproduced, transmitted, transcribed, or translated into any language in any form by any means without the written permission of A&D Company Ltd.

The contents of this manual and the specifications of the instrument covered by this manual are subject to change for improvement without notice.



1WMPD4003537A

For safety use

When this device is operated, mind the following precautions.

CAUTION

Do not disassemble the BT converter. Failure to keep so may result in malfunctions with the BT converter and an elimination of the function, then that is not covered by a warranty.

Precautions on use

CAUTION

- Providing the power from the weighing device, printer or USB to the BT converter, the power can be automatically turned on by connecting the each BT converter to the weighing device, printer or PC.
- This product may require certification for compliance with local laws pertaining to radio broadcast and equipment for wireless communication. Please refer to A&D's website for countries where the product is already certified for use.

1. Introduction

This document is an instruction manual used for AD-8529PC-W Bluetooth converter for PC or AD-8529PR-W Bluetooth converter for printer. Read this instruction manual through before use.


2. Package contents

Either set of the following is packed.

- AD-8529PC-W Bluetooth converter for PC
- AD8529-BA Converter for balance (gray) 1 piece
- AD8529-PC Converter for PC (gray) 1 piece
- Instruction manual (this document) 1 piece
- CD-ROM 1 piece

- AD-8529PR-W Bluetooth converter for printer
- AD8529-BA Converter for balance (gray) 1 piece
- AD8529-PR Converter for PC (blue) 1 piece
- Instruction manual (this document) 1 piece

3. Appearance and each part name

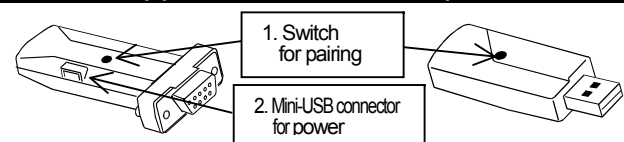

AD8529-BA (gray converter for balance) AD8529-PC (gray converter for PC)
AD8529-PR (blue converter for printer)

- 1. Switch for pairing**
Perform a pairing between the devices.
Use of this switch is normally unnecessary because the pairing has been done beforehand at shipping.
- 2. Mini-USB connector for power**
This is a connector for when using a model that requires the AC adapter sold separately.
The LED for communication confirmation is located on the circuit board, and it blinks in red or green when the power is turned on.

4. Feature


- The BT converter makes the RS-232C cable connection wireless.
- Providing the power from the weighing device to the BT converter, the BT converter can be used by simply inserting to the device.
- The AD-8529PC-W for PC connection is identified as COM port at a PC, so the existing application software such as WinCT to be operated on the Windows can be used.

5. How to use

- Note** - When the converter for PC of set for PC connection is used, it is necessary to install the driver in the PC beforehand. The driver is included in the provided CD-ROM, or download it from the screen for "Software" of "Support" at the A&D's website.
[<http://www.aandd.jp/products/software/software.html>]
Refer to "12. Installing the driver" when confirming installing method for the driver.
Driver's installation is only required when the converter is connected to the PC at the first time.
- When a type of connector for the weighing device is D-Sub25 or DIN7 pin, it is necessary to use a cable described in "9. Accessory".
 - Some weighing devices and any printers other than the AD-8127 separately require an AC adapter, conversion cable for printer connection, etc. Refer to the A&D's website for details.
[A&D website: <http://www.aandd.jp>]

5-1. When using AD-8529PC-W Bluetooth converter for PC


1. Connect the AD8529-BA (gray) to the RS-232C connector for the weighing device to be used.
2. Connect the AD8529-PC (gray) to the USB port on the PC.
3. Confirm that the LED on the AD8529-PC is blinking in green.
If the LED is not blinking in green, refer to "7. When seeming to be malfunctions".
The communication becomes effective by starting up application software such as WinCT on the PC.
Confirm a COM port number indicated at "AND Bluetooth Converter" on the device manager.
Refer to the provided CD-ROM or A&D's website for details.

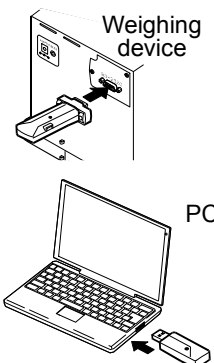

5-2. When using AD-8529PR-W Bluetooth converter for printer

The following procedure is when the AD-8127 printer is used.

1. Connect the AD8529-BA (gray) to the RS-232C connector for the weighing device to be used.
2. Connect the AD8529-PR (blue) to the RS-232C connector for the printer to be used.
3. The communication becomes effective when the LED on the AD8529-PR blinked in green. Data can be sent to the printer by pressing the PRINT key on the weighing device.
* The balances or printers are set so that the BT converter can be connected using factory settings to them.

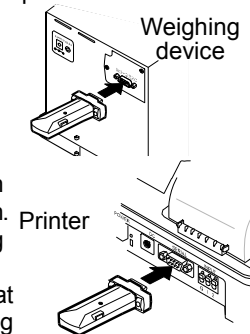


6. Pairing method

The pairing for the BT converter is normally unnecessary because it has been done beforehand.

When using multiple BT converters and having lost track of pairings, it is possible to perform pairing again by the following procedure:

1. Connect the each converters to the device, and press and hold the switch for pairing on the converter for PC or printer until the LED is being lit in orange. (Approx. 3 or 4 seconds)
2. Press and hold the switch for pairing on the converter for balance until the LED on the converter for PC or printer is being blinked in red.
3. The pairing completes when the LED on the converter for PC or printer blinked in green after blinking it in red.

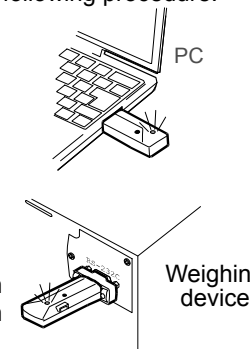

7. When seeming to be malfunctions

Condition	Possible cause	Remedy
When the communication cannot be done	Are the LEDs for each converter blinking in green?	Disconnect the converter and connect it again, or perform the pairing again. Refer to "6. Pairing method".
	Are the settings between the balance and PC matched?	When using AD-8529PC-W... Let the settings between the balance and application software operated on the PC match. When using AD-8529PR-W... Set the balance or printer settings as follows. - Baud rate: 2400 bps - Parity: Even - Bit length: 7bit - Stop bit: 1 - Terminator: CR/LF
	Are the converters connected to the appropriate devices?	The converter for balance is blue and the converter for printer is gray.
	Has the driver been installed in the PC?	When using the AD-8529PC-W, it is necessary to install the driver in the PC. →12. Installing the driver
	Is the proper COM port displayed on the application software that you operate on the PC?	Disconnect the converter and connect it again, and confirm that the proper COM port has been displayed.
When the LEDs do not light up	Is the device a model that requires use of other AC adapter?	Refer to the applicable list for AD-8529 at the A&D's website.

8. Specifications

Communication distance	10m at maximum
Specified power	Not required. (Some balances require the AC adapter sold separately. *1)
Applicable device	A&D's electronic balance with RS-232C interface *1 A&D's printer... AD-8127
Applicable OS	WindowsXP / VISTA / 7 / 8.1 / 10 (compatible with 32bit and 64bit)
Ambient temperature and humidity	5°C to 40°C, Less than 85%R.H. (Do not allow condensation)
Dimensions	Converter for balance or printer... Approx. 19mm (W)×17mm (H)×65mm (D) Converter for PC... Approx. 22mm (W)×12mm (H)×66mm (D)
Weight	Converter for balance or printer... Approx. 18g Converter for PC... Approx. 12g
Wireless communication	CYBLE-022001-00 (Cypress Semiconductor) Bluetooth4.1 (Bluetooth Low Energy)

RS-232C interface	AD8529-BA Converter for balance D-sub9 pin female			
	Pin No.	Signal name	Direction	Description
	2	RxD	Input	Receive data
	3	TxD	Output	Transmit data
	5	GND	-	Ground
	6	DSR	Input	Power
	9	-	Input	Power

RS-232C interface	AD8529-PR Converter for printer D-sub9 pin female			
	Pin No.	Signal name	Direction	Description
	2	TxD	Output	Receive data
	3	RxD	Input	Transmit data
	4	-	Input	Power
	5	GND	-	Ground
	9	-	Input	Power

The converter for balance takes power from the pin 6 or 9.
The converter for printer takes power from the pin 4 or 9.

Settings	AD-8529PC-W	
	Parameter	Value
	Baud rate	: 600, 1200, 2400, 4800, 9600, 14400, 19200, 28800bps
	Bit length / parity	: 7bit/Even, 7bit/Odd, 8bit/None
	Stop bit	: 1, 1.5, 2bit
	AD-8529PR-W	
	Baud rate	: 2400bps
	Bit length / parity	: 7bit/Even
	Stop bit	: 1bit

*1 Refer to those items at the A&D's website.
Windows is registered trademark of the Microsoft Corporation.

9. Accessories

- AX-KO5363-30... Connection cable for a weighing device of D-sub25 pin
- AX-KO5543-100... Connection cable for a weighing device of DIN7 pin

10. About Bluetooth Communication

When using this device for data transfer ensure an unobstructed distance of no more than 10 meters. This device uses Bluetooth for wireless communication and obstacles such as walls can reduce data transmission distance.

Devices such as the following that use the 2.4GHz bandwidth may cause interference when used close to this device.

- Wireless LAN
- Bluetooth capable devices (cellular phone, etc.)
- Home appliances such as microwave ovens

Therefore, this device is not suitable for applications where reliability is the primary concern.

11. Installing the driver

When the converter for PC of the set for PC connection (AD8529-PC) is used, it is necessary to install the driver in the PC. Refer to "Installing the driver" described in the provided CD-ROM for details of installation method for Windows Vista or XP.

11-1. When installing the driver at Windows 10


1. Connect the AD-8529-PC (converter for PC) to the PC.
2. Windows10 leads so as to install the driver.
3. When installation is finished, click the [Close] button.
Device name is "USB Serial Device".

11-2. When installing the driver at Windows7or Windows 8.1


1. Connect the AD-8529-PC (converter for PC) to the PC.
2. In case of Windows8.1... Navigate from "Start" → "PC settings" → "Control panel" to display the "Control Panel" and click the "Hardware and Sound".

In case of Windows7... Navigate from "Start" → "Control panel" to display the "Control Panel" and click the "Hardware and Sound".


3. Click the "Device manager".
4. Confirm the "AND Bluetooth Converter" is displayed.

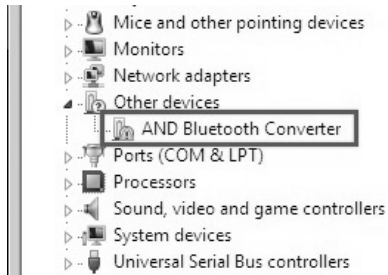


5. Right click "AND Bluetooth Converter", and click "Update Driver Software" on pull-down menu.

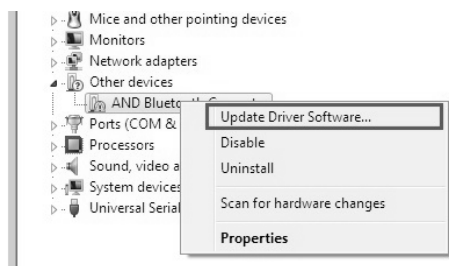

6. Click "Browse my computer for driver software" in the Update Driver Software dialog box.

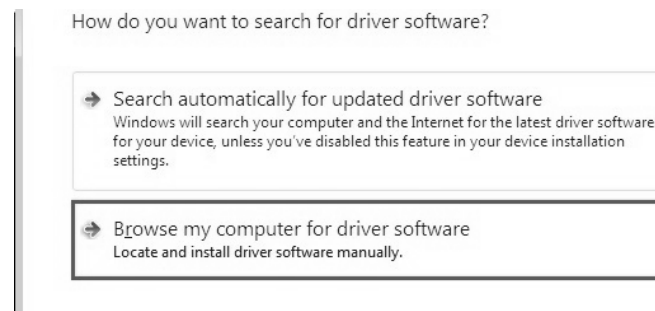


7. Click the "Browse" in the dialog box shown below.

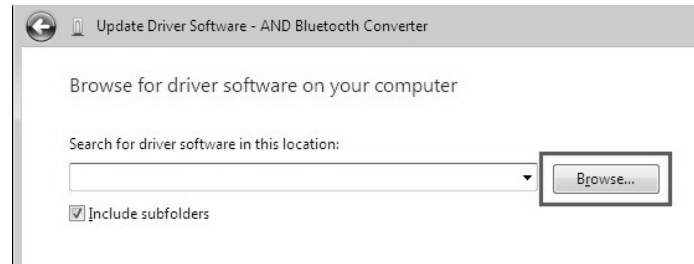


8. Select "AD-8529 DRIVER" in the folder of the provided CD-ROM and click [OK] button.



9. When the dialog box for specifying file location reappears, click the [Next] button.

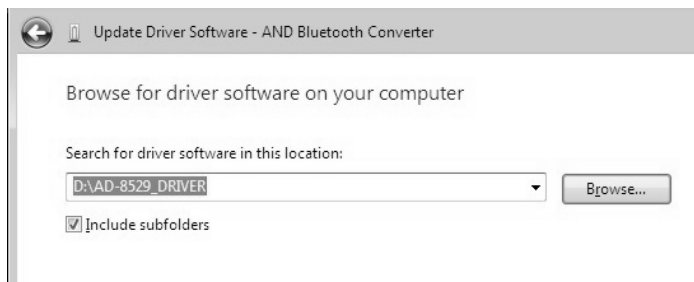

10. The driver will begin to install. If the security popup below appears during installation, click the [Install] button.

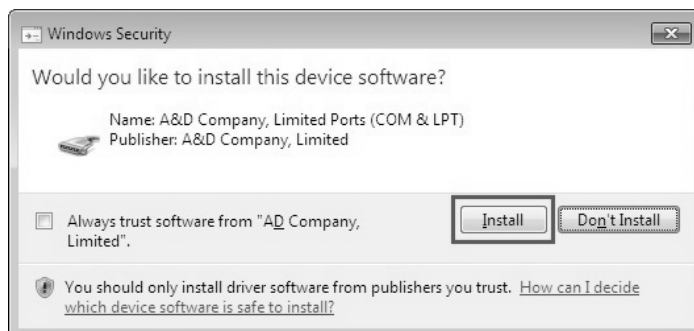

11. When installation is finished, click the [Close] button.

12. Using the WinCT Application Software

The WinCT setup file is included in the CD-ROM.

WinCT (Windows Communication Tools) is a free software for transmitting the data from the A&D electronic balance to a PC (Windows data communication software). WinCT consists of three applications: RsCOM, RsKey and Rsweight. Refer to Readme with WinCT for details of usage with WinCT.

12-1. Installing the WinCT


1. Close all open applications in Windows .
2. Run [WinCT_Setup.msi] in WinCT folder in the provided CD-ROM.
3. Install WinCT following the instructions on the screen

12-2. Communication example using WinCT (RsCOM)

1. Connect the converter for balance to the balance, and connect the converter for PC to the PC. Confirm that the LED on the converter for PC is blinking in yellow.
2. Run "RsCOM" menu. Default menu address is "Start"→"Program"→"A&D WinCT"→"RsCom".

3. Set the computer's RS-232C interface parameters as follows to adjust to the balance's parameters
 - [Port:] Adjust the value to COM port of "AND Bluetooth Converter". (When using Windows 10: "USB Serial Device")
 - [Baud Rate] Baud rate.
 - [Parity] Parity.
 - [Length] Data length.
 - [Stop Bit] Stop bit.
 - [Terminator] Terminator.

* The default settings for RsCom communication parameters are the same as the factory settings for A&D's balances

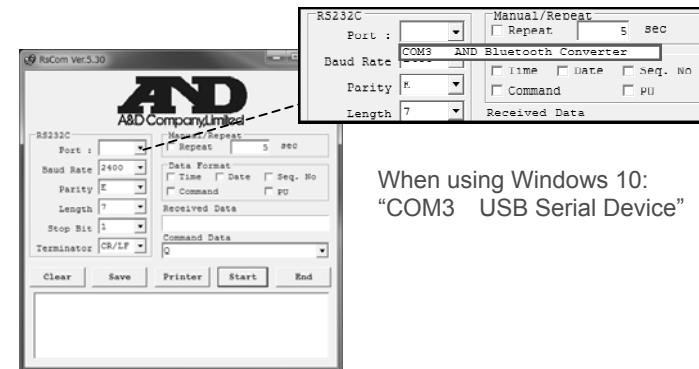

4. By clicking the [Start] button, blinking "o" mark at the top of the display is started, which means that the communication becomes effective.
5. When the stability mark of a balance is displayed and the [PRINT] key of a balance is pressed, the current data is transmitted to a computer.
6. Check the weighing data at [Received Data] in the RsCom window.



7. Select a command on [Command Data] in the RsCom window. Select (click) the [Command] key. Then command is sent to the balance.



8. Select (click) the [Stop] key in the RsCom window. The "o" mark stops blinking on the window.
9. Click the [End] button to close RsCom.

13. Precautions related to radio wave

13-1. FCC

Contains FCC ID:WAP2001

This device complies with Part 15 of the FCC Rules. Operation is subject to the following two conditions: (1) this device may not cause harmful interference, and (2) this device must accept any interference received, including interference that may cause undesired operation

FCC WARNING

Changes or modification not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment.

Note: This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.

13-2. IC

IC RADIATION EXPOSURE STATEMENT FOR CANADA
This device complies with Industry Canada licence-exempt RSS standard(s). Operation is subject to the following two conditions: (1) this device may not cause interference, and (2) this device must accept any interference, including interference that may cause undesired operation of the device.

Le présent appareil est conforme aux CNR d'Industrie Canada applicables aux appareils radio exempts de licence. L'exploitation est autorisée aux deux conditions suivantes : (1) l'appareil ne doit pas produire de brouillage, et (2) l'utilisateur de l'appareil doit accepter tout brouillage radioélectrique subi, même si le brouillage est susceptible d'en compromettre le fonctionnement.

A&D

A&D Company, Limited

3-23-14 Higashi-Ikebukuro, Toshima-ku, Tokyo 170-0013, JAPAN
Telephone: [81] (3) 5391-6132 Fax: [81] (3) 5391-6148