

GXA-12 Animal Weighing Bowl Kit Instruction Manual

Applicable to all GX-A and GF-A series balances except GX-203A and GF-203A.

Total mass of the animal weighing bowl kit is approximately 190 g.

Accessories	Animal weighing bowl	1
	Hook	2
	GXA-12 weighing pan (Use if the minimum division of the balance is 0.01 g or 0.1 g .)	1

Step 1 Select "the usable **weighing pan**".

- If the minimum division is 0.001 g, use "the **weighing pan** of the balance".
- If the minimum division is 0.01 g or 0.1 g, use "the **GXA-12 weighing pan**" and remove "the **weighing pan** of the balance".

Step 2 Place "the usable **weighing pan**" selected on a flat surface such as a desk. Place the animal weighing bowl on it.

Step 3 Hang the upper end of the hook to the animal weighing bowl.

Step 4 Slip the lower end of the hook under one side of the weighing pan at the center.

Step 5 In the same way as described in Step 3 and Step 4, attach the other hook as shown so that the two hooks are placed diagonally.

Step 6 Confirm that the two hooks are attached firmly. Place the assembled unit on the balance.

⚠ Caution

- To remove the animal weighing bowl:
First, remove the assembled unit consisting of the weighing pan, animal weighing bowl and hooks from the balance.
Then, remove the animal weighing bowl.
- If the hooks are loose and come off easily, bend the hook ends to adjust the angles.

